EN - English	1
FR - Français	18
ES - Español	36
IT - Italiano	54
DE - Deutsch	72
PT - Português	90
NL - Nederlands	108
NO - Norsk	113
DA - Dansk	118
SV - Svenska	123
FI - Suomi	128
HU - Magyar	133
PL - Polski	138

European Catalogue Numbers:

Orbital Shakers	Euro Plug	UK Plug	Swiss Plug
3500 Standard Orbital Shakers	444-2903	444-2904	444-2905
3500 Advanced Orbital Shakers	444-2906	444-2907	444-2908
3750 Advanced Reciprocating Shakers	444-2909	444-2910	444-2911
5000 Standard Orbital Shakers	444-2912	444-2913	444-2914
5000 Advanced Orbital Shakers	444-2915	444-2916	444-2917
10000-1 Advanced Orbital Shakers	444-2918	444-2919	444-2920
10000-2 Advanced Orbital Shakers	444-2921	444-2922	444-2923
15000-1 Advanced Orbital Shakers	444-2924	444-2925	444-2926
15000-2 Advanced Orbital Shakers	444-2927	444-2928	444-2929

TABLE OF CONTENTS

Package Contents	1
Warranty	1
Installation	1
Maintenance & Servicing	2
Intended Use	2
Equipment Disposal	2
Environmental Conditions	2
Safety Instructions	3
Warning	3
Standards & Regulations	3
Specifications, Model 3500, 3750	4
Specifications, Models 5000, 10000	5
Specifications, Models 15000	6
Operating Instructions, Standard Orbital Shaker	7
Control Panel, Advanced Orbital Shaker	8
Operating Instructions, Advanced Orbital Shaker	9-11
Technical Service	12
Troubleshooting	12
Accessories	13-17
EC Declaration of Conformity	143-144

PACKAGE CONTENTS

Orbital or Reciprocating Shaker, Non-skid rubber mat,
234cm detachable power cord, Instruction manual

WARRANTY

VWR International warrants that this product will be free from defects in material and workmanship for a period of two (2) years from date of purchase. If a defect is present, VWR will, at its option, repair, replace, or refund the purchase price of this product at no charge to you, provided it is returned during the warranty period. This warranty does not apply if the product has been damaged by accident, abuse, misuse, or misapplication, or from ordinary wear and tear. For your protection, items being returned must be insured against possible damage or loss. This warranty shall be limited to the replacement of defective products. IT IS EXPRESSLY AGREED THAT THIS WARRANTY WILL BE IN LIEU OF ALL WARRANTIES OF FITNESS AND IN LIEU OF THE WARRANTY OF MERCHANTABILITY.

INSTALLATION

Upon receiving the VWR Orbital or Reciprocating Shaker, check to ensure that no damage has occurred during shipment. It is important that any damage that occurred in transport is detected at the time of unpacking. If you do find such damage the carrier must be notified immediately. After unpacking, place the shaker on a level bench or table, away from explosive vapors. Ensure that the surface on which the unit is placed will withstand typical heat produced by the unit and place the unit a minimum of 15cm from vertical surfaces. Always place the unit on a sturdy work surface. The Orbital or Reciprocating Shaker is supplied with a power cord that is inserted into the IEC connector on the back of the unit first, then it can be plugged into a properly grounded outlet. The 230V unit plugs into a 230 volt, 50/60 Hz source. To prevent the rubber mat from sliding, remove backing from adhesive on the underside of mat, reposition in tray, and firmly press.

MAINTENANCE & SERVICING

The Orbital or Reciprocating Shaker is built for long, trouble-free, dependable service. No lubrication or other technical user maintenance is required. It needs no user maintenance beyond keeping the surfaces clean. However at least every three (3) months you should:

- Unplug the unit.
- Remove any accumulated dirt from the base and tray.
- Check all accessible items to make sure they are properly tightened.

The unit should be given the care normally required for any electrical appliance. Avoid wetting or unnecessary exposure to fumes. **DO NOT** use a cleaning agent or solvent on the front panel which is abrasive or harmful to plastics, nor one which is flammable. Always ensure the power is disconnected from the unit prior to any cleaning. If the unit ever requires service, contact your VWR representative.

INTENDED USE

These Shakers are intended for general laboratory use.

EQUIPMENT DISPOSAL

This equipment must not be disposed of with unsorted waste. It is your responsibility to correctly dispose of the equipment at life-cycle-end by handing it over to an authorized facility for separate collection and recycling. It is also your responsibility to decontaminate the equipment in case of biological chemical and/or radiological contamination, so as to protect the persons involved in the disposal and recycling of the equipment from health hazards.

For more information about where you can drop off your waste of equipment, please contact your local dealer from whom you originally purchased this equipment. By doing so, you will help to conserve natural and environmental resources and you will ensure that your equipment is recycled in a manner that protects human health.

ENVIRONMENTAL CONDITIONS

STANDARD MODELS

Operating Conditions: Indoor use only.

* For use in dry CO₂ environments, incubators, or cold rooms. Not for use in CO₂-incubators with high humidity.

Temperature: 0 to 40°C

Humidity: maximum 80% relative humidity, non-condensing

Altitude: 0 to 2000 M above sea level

Non-Operating Storage:

Temperature: -20 to 65°C

Humidity: maximum 80% relative humidity, non-condensing

Installation Category II and Pollution Degree 2 in accordance with IEC 664.

ADVANCED & RECIPROCATING MODELS

Operating Conditions: Indoor use only.

* For use in dry CO₂ environments, incubators, or cold rooms. Not for use in CO₂-incubators with high humidity.

Temperature: -10 to 60°C

Humidity: maximum 80% relative humidity, non-condensing

Altitude: 0 to 2000 M above sea level

Non-Operating Storage:

Temperature: -20 to 65°C

Humidity: maximum 80% relative humidity, non-condensing

Installation Category II and Pollution Degree 2 in accordance with IEC 664.

* **Avoid cold starts:** Unit is not designed to start after being in a cold room environment. Bring unit into cold room from a room temperature environment, operate and remove unit from cold room as soon as operation is complete.

SAFETY INSTRUCTIONS

Please read the entire instruction manual before operating the Orbital Shaker.

WARNING! DO NOT use the Orbital Shaker in a hazardous atmosphere or with hazardous materials for which the unit was not designed. Also, the user should be aware that the protection provided by the equipment may be impaired if used with accessories not provided or recommended by the manufacturer, or used in a manner not specified by the manufacturer.

Always operate unit on a level surface for best performance and maximum safety.

DO NOT lift the unit by the tray.

CAUTION! To avoid electrical shock, completely cut off power to the unit by disconnecting the power cord from the unit or unplug from the wall outlet. Disconnect unit from the power supply prior to maintenance and servicing.

Spills should be removed promptly. **DO NOT** immerse the unit for cleaning.

DO NOT operate the unit if it shows signs of electrical or mechanical damage.

Earth Ground - Protective Conductor Terminal

Alternating Current

Pinch Point - Keep fingers clear during operation.

STANDARDS & REGULATIONS

VWR international hereby declares under its sole responsibility that the products comply with the requirements of the following directives and associated standards:

Associated EU directives:

EMC directive	2014/30/EU
LVD directive	2014/35/EU
ROHS directive	2011/65/EU
WEEE directive	2012/19/EU

Safety standards:

EN 61010-1
EN 61010-2-051
EN 50581

EMC standards:

EN 61326-1

SPECIFICATIONS

MODEL 3500

Overall dimensions LxWxH	41,3 x 35,6 x 14,6cm	
Tray dimensions LxW	33 x 28cm	
Electrical	230 volts, 2,5 amps, 75 watts	
Fuses	5mm x 20mm, 5 amp quick acting	
Speed Range	(Advanced)	15 to 500 rpm
	(Standard)	25 to 500 rpm
Speed Accuracy	(Advanced)	
	Above 100rpm	±1% of set speed
	Below 100rpm	±1rpm
Orbit	19mm	
Capacity	~16kg @ 75 rpm ~2,3kg @ 300 rpm	
Timer	(Advanced)	1 second to 160 hours
	(Standard)	1 to 120 minutes
Controls	(Advanced)	See page 8
	(Standard)	See page 7
Ship Weight	22,2kg	
European Catalog Numbers:	(Standard)	(Advanced)
	Euro plug:	444-2903 444-2906
	UK plug:	444-2904 444-2907
	Swiss plug:	444-2905 444-2908

MODEL 3750

Overall dimensions LxWxH	41,3 x 35,6 x 14,6cm	
Tray dimensions LxW	33 x 28cm	
Electrical	230 volts, 2,5 amps, 40 watts	
Fuses	5mm x 20mm, 5 amp quick acting	
Speed Range	20 to 300 rpm	
Speed Accuracy	(Advanced)	
	Above 100rpm	±1% of set speed
	Below 100rpm	±1rpm
Orbit	19mm Total Travel	
Capacity	6,8kg @ 300 rpm	
Timer	1 second to 160 hours	
Controls	See page 8	
Ship Weight	22,2kg	
European Catalog Numbers:	(Advanced)	
	Euro plug:	444-2909
	UK plug:	444-2910
	Swiss plug:	444-2911

SPECIFICATIONS (CONT'D)

MODEL 5000			MODEL 10000		
Overall dimensions LxWxH	59,9 x 67,8 x 15,2cm		Overall dimensions LxWxH	71,8 x 67,8 x 17,8cm	
Tray dimensions LxW	45,7 x 61cm		Tray dimensions LxW	61 x 61cm	
Electrical	230 volts, 2,5 amps, 75 watts		Electrical	230 volts, 2,5 amps, 80 watts	
Fuses	5mm x 20mm, 5 amp quick acting		Fuses	5mm x 20mm, 5 amp quick acting	
Speed Range	(Advanced)	20 to 500 rpm	Speed Range	10000-1 =15 to 500 rpm	
	(Standard)	25 to 500 rpm		10000-2 =15 to 300 rpm	
Speed Accuracy	(Advanced)		Speed Accuracy	(Advanced)	
	Above 100rpm	±1% of set speed		Above 100rpm	±1% of set speed
	Below 100rpm	±1rpm		Below 100rpm	±1rpm
Orbit	25.4mm		Orbit	10000-1 = 2,54cm	
Capacity	~22,3kg @ 125 rpm			10000-2 = 5,0cm	
	~4,5kg @ 300 rpm		Capacity	~45,3 @ 100 rpm	
Timer	(Advanced)	1 second to 160 hours		~18,1 @ 300 rpm	
	(Standard)	1 to 120 minutes	Timer	1 second to 160 hours	
Controls	(Advanced)	See page 8	Controls	See page 8	
	(Standard)	See page 7	Ship Weight	90,7kg	
Ship Weight	49,5kg		European Catalog Numbers:	(10000-1 Advanced)	(10000-2 Advanced)
European Catalog Numbers:	(Standard)	(Advanced)	Euro plug:	444-2918	444-2921
Euro plug:	444-2912	444-2915	UK plug:	444-2919	444-2922
UK plug:	444-2913	444-2916	Swiss plug:	444-2920	444-2923
Swiss plug:	444-2914	444-2917			


SPECIFICATIONS (CONT'D)

MODEL 15000

Overall dimensions LxWxH	71,8 x 67,8 x 17,8cm	
Tray dimensions LxW	61 x 91,4cm	
Electrical	230 volts, 2,5 amps, 80 watts	
Fuses	5mm x 20mm, 5 amp quick acting	
Speed Range	15000-1 =15 to 500 rpm 15000-2 =15 to 300 rpm	
Speed Accuracy	(Advanced)	
	Above 100rpm	±1% of set speed
	Below 100rpm	±1rpm
Orbit	15000-1 = 2,54cm 15000-2 = 5,0cm	
Capacity	~68,0kg @ 100 rpm ~18,1kg @ 300 rpm	
Timer	1 second to 160 hours	
Controls	See page 8	
Ship Weight	104,3kg	
European Catalog Numbers:	(15000-1 Advanced)	(15000-2 Advanced)
	Euro plug: 444-2924	444-2927
	UK plug: 444-2925	444-2928
	Swiss plug: 444-2926	444-2929

OPERATING INSTRUCTIONS - STANDARD ORBITAL SHAKER

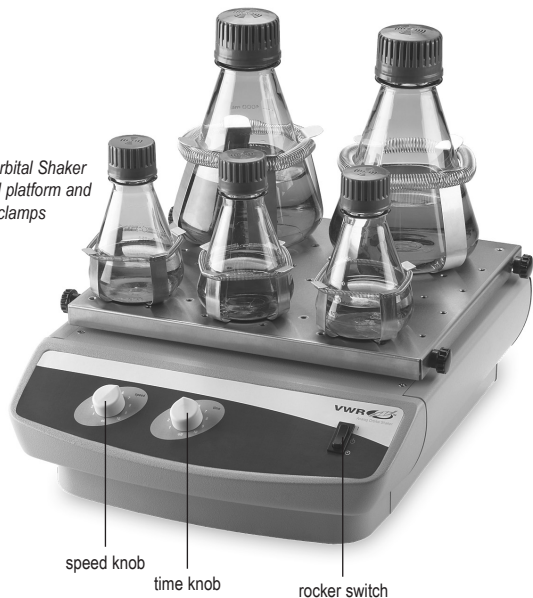Standard Orbital Shaker
with optional platform and
flask clamps

The Standard Orbital Shaker is used for general laboratory shaking needs.

Models 3500 and 5000

1. Getting ready:

- a. The speed and time knobs have a built in on-off switch at their extreme counter-clockwise position. Turn both knobs to the off position.
- b. Make sure the rocker switch is in the standby position.
- c. Plug the cord into a properly grounded outlet.

2. Setting speed:

- a. To run in continuous mode, push rocker to the run position. Set speed knob to desired setting and adjust if necessary. Unit will run until you move the rocker switch to the standby position. The microprocessor speed control slowly ramps to set speed to avoid splashing.

3. Setting time:

- a. To run in timed mode, set timer and the speed knob to the desired settings. Press the rocker switch to the time position and release. The shaker will now run for the set time.
- b. To exit the timed mode turn the time knob to the extreme counter-clockwise position.

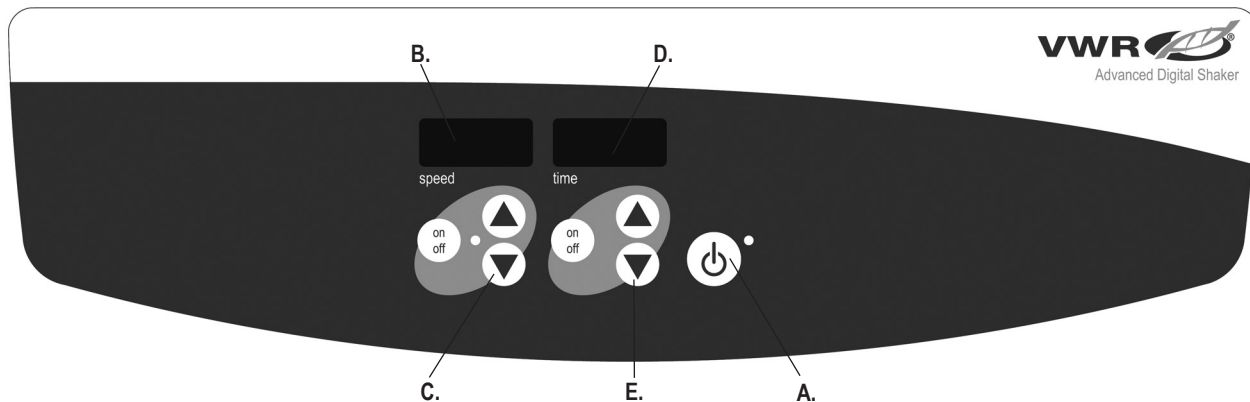
4. Turning unit off:

- a. To stop shaking function, turn the speed knob to the extreme counter-clockwise position and push the rocker switch to the standby position. The shaker should be kept in the standby position when not in use. To completely cut off power to the unit, disconnect the power cord from the unit or unplug from the wall outlet.

OPERATING TIPS

Centering your sample and even weight distribution on the tray helps with balance and stability. As a safety feature, a built-in program will shut power off to the motor if the tray is prevented from rotating, or the unit is overloaded beyond its recommended weight capacity. To reset the unit, press the rocker switch to standby and then press the rocker switch to the on position. The shaker will automatically restart after a power interruption. Built-in memory maintains the last used speed and time settings during a power interruption.

To prevent the rubber mat from sliding, remove backing from adhesive on the underside of mat, reposition in tray, and firmly press.

CONTROL PANEL - ADVANCED ORBITAL SHAKER

The front panel of the Advanced Orbital Shaker contains all the controls and displays needed to operate the unit.

- A. Standby button/standby indicator light:** The standby indicator light will illuminate when the unit is plugged in. The unit will be in standby mode. Press the standby button to start the speed and time functions. The standby indicator light will shut off. Press the standby button again and the unit will once again be in standby mode.
- B. Speed display:** Displays the speed of the shaker.
- C. Up/down arrows:** For set-point control. On/off button starts/stop shaking function.

- D. Time display:** Displays accumulated time (continuous mode) or how much time is remaining (timed mode). The display range is from 0 to 9999 minutes in one (1) second increments. The display will indicate minutes and seconds until the timer reaches 99 minutes and 59 seconds (99:59), then the display will automatically display minutes up to 9999.
- E. Up/down arrows for setpoint control.** On/off button starts/stops the timer function.

Advanced Orbital Shaker
with optional platform

The Advanced Orbital Shaker is used for general laboratory shaking needs.

Advanced Models

The Advanced Orbital Shaker has been designed for the speed and time functions to work independently of one another. The speed can be reset without resetting the timer and the timer can be stopped and started without interrupting the shaking function.

1. Getting ready:

- a. Plug the cord into a properly grounded outlet. The standby indicator light will illuminate, verifying power to the shaker.
- b. Press the standby button to move the unit from standby mode. The standby indicator light will turn off and the speed and time displays will illuminate, displaying the previously used settings.

2. Setting speed:

- a. Press the up/down arrow below the speed display until you reach the desired speed. When you release the button, the display will blink off and then on indicating the new set speed has been accepted.
- b. Press the on/off button to start the shaking function. The indicator light below the speed display will illuminate and blink until the setpoint is reached. Once the set-point is reached the light will stop blinking and remain lit until shaking has ceased. The microprocessor controlled ramping feature slowly increases speed until the set-point is reached which helps to avoid splashing, and provides excellent low end control.
- c. Speed adjustments can be made without interrupting shaking by using the up/down arrows below the speed display. After the change has been made and you release the button, the display will blink off and then on indicating the new set speed has been accepted.
- d. To stop the shaking function press the on/off button below the speed display. The speed indicator light will turn off.

3. Setting time to zero (0:00) and continuous mode: Accumulated time.

- a. Press and hold the on/off button below the time display. After three (3) seconds, the display will indicate the previous set time.

OPERATING INSTRUCTIONS - ADVANCED ORBITAL SHAKER (CONT'D)

- b. Simultaneously press both the up and the down arrows, the display will indicate zero (0:00). The unit time is now set to zero (0:00) minutes. Alternately, you can use the up/down arrows to get to zero (0:00).
- c. Press the on/off button below the time display. The display will indicate accumulated time. The up/down arrows will become inactive. To stop timer, press the on/off button again. **IMPORTANT:** This will **NOT** interrupt the shaking function. Press the on/off button below the speed display to interrupt the shaking function.
- d. To reset, press and hold the on/off button below the time display. After three (3) seconds the display will indicate the previous set time, which was zero (0:00).

4. Setting timed mode: Programmed time.

- a. Press the up/down arrows below the time display until you reach the desired time.
- b. Start this function by pressing the on/off button below the time display. The unit will run for the selected time, the up/down arrows will become inactive while the timer is running. The unit will stop shaking when time display reaches zero (0:00). Four (4) audible beeps will indicate the count down function is complete. The time display will default back to the set time. To repeat for the same time, simply depress the on/off button again.
- c. To interrupt an automatic timing cycle before it is completed, press the on/off button below the time display. The display will flash off and on to indicate the time function is on "hold".

IMPORTANT: This will **NOT** interrupt the shaking function. Press the on/off button below the speed display to interrupt the shaking function. Restart the timer by pressing the on/off button below the time display. Unit will continue counting down to zero (0:00). When the display reaches zero (0:00), you will hear the four (4) audible beeps that indicate the count down function is complete and the shaking function will cease.

5. Turning unit off:

- a. To turn the unit off, press the standby button. The speed and time displays will be blank, the standby indicator light will illuminate. The Orbital Shaker should be kept in standby mode when not in use. To completely cut off power to the unit, disconnect the power cord from the unit or unplug from the wall outlet.

LOAD SENSING FUNCTION

The Advanced Orbital Shaker Models 3750 and up is equipped with a load sensing function that can be activated by the user. This function provides protection against improper positioning of load and maximum load being exceeded. When activated, the unit will automatically sense improper load conditions and slow to a safe running speed, then display that speed followed by E04 on the speed display. The unit will also beep three (3) times every 60 seconds until the error is reset by pressing the speed on/off button. To activate the load sensing function use the following steps:

1. Place the unit in standby mode.
2. Press and hold the speed on/off button and press the standby button. The unit will beep two (2) times, confirming the function is enabled.
3. To restore normal operation, remove AC power to the unit for ten (10) seconds and then restore. If the E04 error occurs be sure the load is within the maximum specification and properly balanced (centered on tray) and/or reduce sample size/speed before restarting the unit. If the E04 occurs due to acceptable sample vibration or another vibration source, the vibration sensing function can be disabled as described above.

OPERATING INSTRUCTIONS - ADVANCED ORBITAL SHAKER (CONT'D)

ADDITIONAL LOAD SENSING FEATURE

The Advanced Orbital Shaker Models 3750 and up, are equipped with an additional overload protection feature that lets the user know when the unit's set speed is higher than the achievable speed of the unit. The unit speed display will show an E7. The unit will also beep three (3) times every 60 seconds until the error is reset by pressing the speed on/off button.

To disable this feature perform the following steps:

1. Place the unit in standby mode by pressing the standby button.
2. Press and hold the speed up button and press on/off button. The unit will beep two (2) times, confirming the function is disabled.
3. To restore unit to normal operation, remove AC power to unit (unplug power cord from wall outlet) for ten (10) seconds and restart.

BEEPER PREFERENCE

To silence beeper operation (except for error codes), with the unit in standby mode, press and hold the time on/off button and press the standby button. To restore normal beeper operation, remove AC power to unit for ten (10) seconds and then restore.

CALIBRATION PROCEDURE (Advanced Orbital Shaker Models 3500 and up)

This procedure is used to self calibrate the Advanced Orbital Shaker. The tray should be free of any samples, vessels, and accessories prior to calibrating.

1. Turn unit on. Speed and time displays will be illuminated.
2. Press and hold the standby button and momentarily press the speed on/off button. The speed display should read "CAL".
3. The unit will run for approximately one (1) minute and automatically calibrate.

RS-232 SERIAL PORT (Advanced Orbital Shaker Models 3500 and up)

RS-232 serial port provides two-way communications for data logging and unit control. If you need additional details, please contact your local VWR After Sales Service Department.

OPERATING TIPS

Centering your sample and even weight distribution on the tray helps with balance and stability. As a safety feature, a built-in program will shut power off to the motor if the tray is prevented from rotating, or the unit is overloaded beyond its recommended weight capacity. The shaker will automatically restart after a power interruption. Built-in memory maintains the last used speed and time settings during a power interruption.

TECHNICAL SERVICE

For more information or technical assistance contact your local VWR representative or visit www.vwr.com.

Web resources: Visit VWR's website at www.vwr.com for:


- Complete technical service contact information
- Access to VWR's Online Catalogue, and information about accessories and related products
- Additional product information and special offers

TROUBLESHOOTING


Problem	Cause	Solution
Unit will not run	Mechanical obstruction Motor obstruction	Add or replace fuse as necessary. If problem persists, please contact your VWR representative for repair.
Unit is excessively noisy	Sensor fan misaligned Motor misaligned	Ensure that tray is secured tightly. If problem persists, please contact your VWR representative for repair.
Unit not shaking at proper speed	-	Perform speed calibration test on page 11 If problem persists, please contact your VWR representative for repair.
E3	Mechanical obstruction Drive system failure Ceased bearing Drive belt broken	Remove mechanical obstruction. If problem persists, the reason may be the drive system and should not be addressed by the end user. Contact your VWR representative for repair.
E4	Improper positioning of load Maximum load exceeded	Ensure the load is evenly distributed and does not exceed the maximum load capacity for the unit. See "Load Sensing Function" on page 10. If problem persists, please contact your VWR representative for repair.
E7	Unable to reach set speed	Reduce the speed setting or weight. This function can be disabled by following the instructions listed under "Additional Load Sensing Feature" on page 11. (NOTE: This function is only available on models 3750 and up and is not available on Model 3500 shakers.)
E8	Electronics error	This error cannot be fixed by the end user. Please contact your VWR representative for repair.

ACCESSORIES

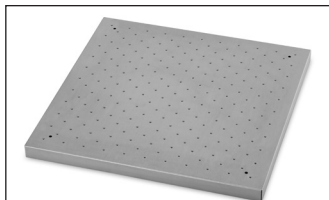Platform Usage Chart

Universal Platform

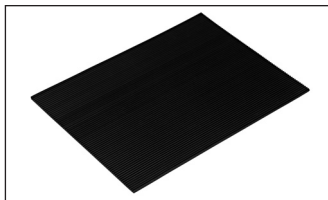Rubber Mat

Dedicated Platform

Tier Braces with Universal Platform

Universal Platform Size	Description	Part Number	Used On Shaker Model Numbers
28 x 33cm	Universal platform with holes for mounting accessories	444-7087	Model 3500, Model 3750
33 x 33cm	Universal platform with holes for mounting accessories	444-2930	Model 3500, Model 3750
46 x 46cm*	Universal platform with holes for mounting accessories	444-2931	Model 3500, Model 3750**
46 x 61cm*	Universal platform with holes for mounting accessories	444-2932	Model 3500, Model 3750**, Model 5000**
46 x 76cm*	Universal platform with holes for mounting accessories	444-2934	Model 5000
61 x 61cm	Universal platform with holes for mounting accessories	444-2933	Model 10000
61 x 91cm	Universal platform with holes for mounting accessories	444-2935	Model 15000
28 x 33cm	Rubber Replacement Mat	444-2966	Model 3500, Model 3750
46 x 61cm	Rubber Replacement Mat	444-2967	Model 5000
61 x 61cm	Rubber Replacement Mat	444-2968	Model 10000
61 x 91cm	Rubber Replacement Mat	444-2969	Model 15000

DEDICATED Platform Size	Dedicated Platforms with PVC Flask Clamp Capacities	Part Number	Used On Shaker Model Numbers
33 x 33cm (125mL)	16 x 125mL	444-2945	Model 3500, Model 3750
33 x 33cm (250mL)	12 x 250mL	444-2946	Model 3500, Model 3750
33 x 33cm (500mL)	8 x 500mL	444-2947	Model 3500, Model 3750
33 x 33cm (1L)	4 x 1L	444-2948	Model 3500, Model 3750
46 x 46cm (125mL)*	27 x 125mL	444-2949	Model 3500, Model 3750 **
46 x 46cm (250mL)*	20 x 250mL	444-2950	Model 3500, Model 3750 **
46 x 46cm (500mL)*	13 x 500mL	444-2951	Model 3500, Model 3750 **
46 x 46cm (1L)*	9 x 1L	444-2952	Model 3500, Model 3750 **
Tier Braces	Provides 25.4cm of clearance between stacking platforms	444-2965	

* Two Tier Ready

** PLATFORM STACKING IS NOT AVAILABLE ON MODEL 3750 AND MODEL 5000

ACCESSORIES (CONT'D)

Platform Usage Chart

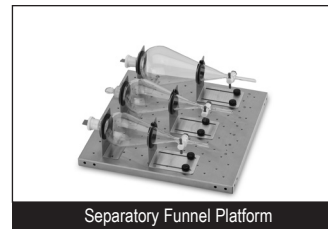
Culture Platform

Adjustable Platform

Large Vessel Carrier

Separatory Funnel Platform

Platform Size	Platform Type	Part Number	Used On Shaker Model Numbers
46 x 46cm*	Culture Platform with anti slip mat for petri dishes	444-2936	Model 3500, Model 3750**
46 x 61cm*	Culture Platform with anti slip mat for petri dishes	444-2937	Model 3500, Model 3750**, Model 5000**
46 x 46cm	Adjustable Platform 3 Roller Bars	444-2938	Model 3500
46 x 61cm	Adjustable Platform 4 Roller Bars	444-2939	Model 5000
46 x 76cm	Large Vessel Carrier with 4 Roller Bars	444-2940	Model 5000
61 x 61cm	Large Vessel Carrier with 4 Roller Bars	444-2941	Model 10000
61 x 91cm	Large Vessel Carrier with 5 Roller Bars	444-2942	Model 15000
46 x 46cm	Separatory Funnel Platform accomodates 3 funnels 500mL to 2L	444-2953	Model 3500, Model 3750
Tier Braces	Provides 25.4cm of clearance between stacking platforms	444-2965	
46cm	Replacement Rollerbar	444-2943	444-2938, 444-2939, 444-2940
61cm	Replacement Rollerbar	444-2944	444-2941, 444-2942

* Two Tier Ready

** PLATFORM STACKING IS NOT AVAILABLE ON MODEL 3750 AND MODEL 5000

ACCESSORIES (CONT'D)

Test Tube Rack Platform Capacity

Full Size Stationary

Rack Size: 9.5 x 41.9 x 10.2cm Platform	10 to 14mm Test Tube Rack capacity = 48 444-7037	16 to 20mm Test Tube Rack capacity = 33 444-7038	21 to 25mm Test Tube Rack capacity = 21 444-7039
Platform Size	Part Number		
28 x 33cm	444-7087	N/A	N/A
33 x 33cm	444-2930	N/A	N/A
46 x 46cm	444-2931	3	3
46 x 61cm	444-2932	5	5
46 x 76cm	444-2934	6	6
61 x 61cm	444-2933	7	7
61 x 91cm	444-2935	5	5

Full Size Pivoting

Internal Rack Size: 12.9 x 27.4 x 9.9cm External Rack Size: 12.7 x 27.6 x 12.7cm Platform	13mm Test Tube Rack capacity = 90 444-2955	16mm Test Tube Rack capacity = 60 444-2956	20mm Test Tube Rack capacity = 40 444-2957	25mm Test Tube Rack capacity = 24 444-2958	30mm Test Tube Rack capacity = 12 444-2959
Platform Size	Part Number				
28 x 33cm	444-7087	1	1	1	1
33 x 33cm	444-2930	2	2	2	2
46 x 46cm	444-2931	2	2	2	2
46 x 61cm	444-2932	3	3	3	3
46 x 76cm	444-2934	4	4	4	4
61 x 61cm	444-2933	4	4	4	4
61 x 91cm	444-2935	6	6	6	6

Half Size Stationary


Rack Size: 12.7 x 17.8 x 10.2cm Microtube Rack: 4.3cm Platform	1.5 to 2mL Microtube Rack capacity = 70 444-0276	10 to 13mm Test Tube Rack capacity = 63 444-0278	14 to 16 mm Test Tube Rack capacity = 48 444-0279	18 to 20 mm Test Tube Rack capacity = 35 444-0280	22 to 25 mm Test Tube Rack capacity = 24 444-0281	15mL Centrifuge Test Tube Rack capacity = 35 444-0283	50mL Centrifuge Test Tube Rack capacity = 12 444-0284
Platform Size	Part Number						
28 x 33cm	444-7087	2	2	2	2	2	2
33 x 33cm	444-2930	2	2	2	2	2	2
46 x 46cm	444-2931	4	4	4	4	4	4
46 x 61cm	444-2932	6	6	6	6	6	6
46 x 76cm	444-2934	8	8	8	8	8	8
61 x 61cm	444-2933	8	8	8	8	8	8
61 x 91cm	444-2935	7	7	7	7	7	7

Full Size Stationary

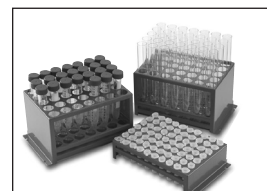Full Size Pivoting

Half Size Stationary

ACCESSORIES (CONT'D)

Flask Clamp Platform Capacity

PVC Erlenmeyer Flask Clamps

Stainless Steel Erlenmeyer Flask Clamps

Stainless Steel Media Bottle Clamps

PVC Erlenmeyer Flask Clamps

Platform Size	Platform Part Number	125mL 444-2960	250mL 444-2961	500mL 444-2962	1L 444-2963	2L 444-2964
28 x 33cm	444-7087	10	8	5	2	N/A
33 x 33cm	444-2930	12	10	6	4	3
46 x 46cm	444-2931	20	18	12	8	4
46 x 61cm	444-2932	28	25	16	10	6
46 x 76cm	444-2934	36	33	20	14	8
61 x 61cm	444-2933	41	35	24	13	9
61 x 91cm	444-2935	61	55	38	22	13

Stainless Steel Media Bottle Clamps

Platform Size	Platform Part Number	500mL 444-7035	1L 444-7036
28 x 33cm	444-7087	5	2
33 x 33cm	444-2930	6	5
46 x 46cm	444-2931	16	10
46 x 61cm	444-2932	20	13
46 x 76cm	444-2934	28	18
61 x 61cm	444-2933	25	18
61 x 91cm	444-2935	40	30

Stainless Steel Erlenmeyer Flask Clamps

Platform Size	Platform Part Number	10mL 444-7040	25mL 444-7041	50mL 444-7042	125mL 444-7043	250mL 444-7044	500mL 444-7045	1L 444-7030	2L 444-7031	2.8L 444-7032	4L 444-7033	6L 444-7034
28 x 33cm	444-7087	60	25	13	10	9	7	4	N/A	N/A	N/A	N/A
33 x 33cm	444-2930	60	30	15	12	12	8	4	3	1	1	1
46 x 46cm	444-2931	113	64	32	20	20	13	8	5	2	4	2
46 x 61cm	444-2932	158	88	44	28	28	20	12	6	3	4	3
46 x 76cm	444-2934	203	112	56	36	36	26	15	8	3	6	4
61 x 61cm	444-2933	221	121	61	41	41	25	16	9	5	5	5
61 x 91cm	444-2935	336	160	94	61	64	40	24	14	7	9	7

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The undersigned states that this declaration is issued under the sole responsibility of the manufacturer, VWR International bvba, and that the object(s) of the declaration described in annex 1 are in conformity with the relevant Union harmonization legislation.
In the event of unauthorized modification of any of the products listed in the annex 1, this declaration becomes invalid.

Object Name: VWR Shakers

Relevant EC Directives:

Low Voltage Directive 2014/35/EU
EMC Directive 2014/30/EU
RoHS 2011/65/EU


Harmonized Standards:

EN 61326-1:2013
EN 61010-1:2001
EN 61010-2-051:2003
EN 50581:2012

Document Number: 14871.2017.01

Place & date of issue: Leuven, 3/03/2017

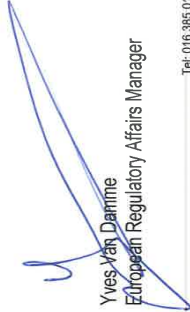Signed for and on behalf of VWR International bvba:

Yves Van Damme
European Regulatory Affairs Manager

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India

VWR Lab Products Private Limited
No.139, BDA Industrial Suburb,
6th Main, Tumkur Road, Peenya Post,
Bangalore, India – 560058
Tel.: +91-80-28078400
Fax +91-80-4117120
Email: vwr_india@vwr.com

Ireland / Northern Ireland

VWR International Ltd / VWR International
(Northern Ireland) Ltd
Orion Business Campus
Northwest Business Park
Ballycoolin
Dublin 15
Tel.: 01 88 22 222
Fax: 01 88 22 333
Email sales.ie@vwr.com

Italy

VWR International S.r.l.
Via San Giusto 85
20153 Milano (MI)
Tel.: 02-3320311/02-487791
Fax: 02-332031307/02-40090010
Email: info.it@vwr.com

The Netherlands

VWR International B.V.
Postbus 8198
1005 AD Amsterdam
Tel.: 020 4808 400
Fax: 020 4808 480
Email: info.nl@vwr.com

Norway

VWR International AS
Haavard Martinsens vei 30
0978 Oslo
Tel.: 22 90 00 00
Fax: 815 00 940
Email: info.no@vwr.com

Poland

VWR International Sp. z o.o.
Limbowa 5
80-175 Gdansk
Tel.: 058 32 38 210
Fax. 058 32 38 205
Email: info.pl@vwr.com

Portugal

VWR International - Material de Laboratório, Lda
Centro Empresarial de Alfragide
Rua da Indústria, nº 6
2610-088 Alfragide
Tel.: 21 3600 770
Fax: 21 3600 798/9
Email: info.pt@vwr.com

Singapore

VWR Singapore Pte Ltd
18 Gul Drive
Singapore 629468
Tel: +65 6505 0760
Fax: +65 6264 3780
Email: sales.sg@vwr.com

Spain

VWR International Eurolab S.L.
C/ Tecnología 5-17
A-7 Llinars Park
08450 - Llinars del Vallès
Barcelona
Tel.: 902 222 897
Fax: 902 430 657
Email: info.es@vwr.com

Sweden

VWR International AB
Fagerstagatan 18a
163 94 Stockholm
Tel.: 08 621 34 00
Fax: 08 621 34 66
Email: info.se@vwr.com

Switzerland

VWR International GmbH
Lerzenstrasse 16/18
8953 Dietikon
Tel.: 044 745 13 13
Fax: 044 745 13 10
Email: info.ch@vwr.com

Turkey

VWR International Laboratuvar
Teknolojileri Ltd.Şti.
Orta Mah. Cemal Gürsel Caddesi
Ördekcioglu İşmerkezi No.32/1
34896 Pendik - Istanbul
Tel.: +90216 598 2900
Fax: +90216 598 2907
Email: info.tr@vwr.com

UK

VWR International Ltd
Customer Service Centre
Hunter Boulevard
Magna Park
Lutterworth
Leicestershire
LE17 4XN
Tel.: 0800 22 33 44
Fax: 01455 55 85 86
Email: uksales@vwr.com

Legal Manufacturer:

VWR International bvba • Researchpark Haasrode 2020 • Geldenaaksebaan 464 • B-3001 Leuven • + 32 16 385011 • <http://be.vwr.com>
715089-00 (Rev 11)