# TABLE OF CONTENTS

SECTION 1 INTRODUCTION1
1.1 The Coulter Principle 1
1.2 About the Multisizer <sup>™</sup> 3 COULTER COUNTER <sup>®</sup>
1.3 About this Manual 4
1.3.1 Conventions
1.3.2 Warnings and Cautions5
1.4 Warnings 5
1.4.1 Electrical
1.4.2 Chemical
1.4.3 Mechanical5
1.4.4 Fire
1.5 Cautions
1.5.1 Electrical
1.5.2 Chemical
1.6 Warming up
1.7 Sources of Error
1.7.1 Calibration constant
1.7.2 Damaged apertures7
1.7.3 Deposits on and around aperture wafers7
1.7.4 External sound and mechanical vibration7
1.7.5 Electrical interference
1.7.6 Particles behind the aperture7
1.7.7 Sample dispersion
1.7.8 Excessive aperture current

SEC1	ΓΙΟΝ 2 INSTALLATION	8
2.1	Location	8
2.2	Unpacking	8
2.3	Connections	9
SECT	FION 3 BASIC OPERATIONS	14
3.1	Starting the Instrument and its Control Program	14
3.2	Instrument Controls	15
3.3	Shutting down the Multisizer 3	17
SECT	FION 4 GETTING STARTED	18
4.1	Locating and Identifying Controls and Parts You Need to Know	18
4.1.	1 The Front Panel of the Multisizer 3	18
4.1.2	2 The Sample Compartment	18
4.2	Fitting the Electrode and Stirrer	20
4.3	Fitting the Aperture Tube	21
4.4	Setting Up the Aperture Tube	22
1.	Removing the currently installed aperture tube.	23
2.	Installing a New Aperture Tube	
3.	Placing a beaker of clean electrolyte on the platform	25
4.	Closing the Sample Compartment door	25
5.	Selecting the electrolyte	26
6.	Selecting the aperture tube	26
7.	Filling the System	27
8.	Adjusting the Metering Pump	27
9.	Setting Current and Gain.	28
10	D. Measuring Noise Level.	
11	1. Calibration and Verfication of the Calibration	29
4.5	Aligning the Optics	30

SECI	TION 5 CALIBRATION AND VERIFICATION OF CALIBRATION	31
5.1	Choosing the Calibrator	
5.2	Preparing the Calibrator	
5.3	Running the Calibrator	33
5.4	Obtaining the mean Kd	35
5.5	Entering the mean value for the Kd in the aperture list	35
5.6	Verifying the calibration	
SECT	TION 6 SAMPLE ANALYSIS	40
6.1	Introduction	40
6.2	Changing Standard Operating Method (SOM).	41
6.3	Convert Pulses to Size Settings	
6.4	Saving Standard Operating Method	50
6.5	Entering Sample Information	51
6.6	Running a Background.	55
6.7	Subtracting the Background	56
6.8	Running the Sample	57
6.10	Cleaning the System	58
SECT	TION 7 REVIEWING AND SAVING THE RESULTS	60
7.1	Introduction	60
7.2	Opening a File	60
7.3	Overlaying Files	63
7.4	Averaging Files	67
7.5	Creating a Size Trend	72
7.6	Overlapping Files from Different Aperture Sizes.	76
SECT	TION 8 PRINTING A REPORT	80
8.1	Introduction	80

8.2 S	electing the Content of the Printed Report	80
8.2.1	Sample Info	81
8.2.2	Size Section of the Printed Report	82
8.2.3	Size Statistic Section of the Printed Report	87
8.2.4	Pulses Section of the Printed Report	89
8.2.5	Trend Section of the Printed Report	
8.2.6	Printer Section	
8.2.7	Closing the Printing Report Dialog Box	93
8.3 C	Customizing the Preferences	
8.3.1	Size	
8.3.2	Trend	101
8.3.3	Pulses	102
8.3.4	Other Options for the Preference File	103
8.3.5	Averaging and Trend	104
8.3.6	Number and Date Formats	105
8.3.7	Data Export	106
8.3.8	Graph Options	106
8.3.9	Page Setup	107
8.3.10	Pick Fonts	107
8.3.11	Windows Colors	108
8.3.12	Customize Toolbar	108
8.3.13	Selecting the Language	109
8.3.14	Saving Preferences	109
SECTI	ON 9 CREATING A STANDARD OPERATING	

	PROCEDURE (SOP)1	11
9.1	Introduction	111

9.2	Creating and Saving an SOP	111
9.2.1	Selecting the SOM file.	112

9.2.2 Selecting the Preferences File	114
9.2.3 Saving the SOP	115
9.3 Using An SOP	116
9.4 Removing Active SOP	117
SECTION 10 REGULATORY COMPLIANCE	118
10.1 21 CFR Part 11	118
10.2 Electronic records	118
10.3 FDA Requirements	118
10.4 Implementation of electronic records and electronic signatures	118
10.5 Controls for electronic records	119
10.6 Establishing an electronic record control with the Multisizer™ 3	119
10.6.1 Setting up security	119
10.6.2 Adding a user account	120
10.6.3 Selecting user names and passwords	121
10.6.4 File History and Audit Trail	122
10.6.5 Setting Audit Trail	124
10.6.6 Selecting user privileges	125
10.6.7 Electronic Signature	126
10.6.8 Generating electronic signatures	126
10.6.9 Applying electronic signatures	127
10.6.10 Additional security features	128
10.7 Starting the Multisizer 3 software with enabled security	128
SECTION 11 SPECIFICATIONS	129
11.1 Dimensions and weight	129
11.2 Environmental Conditions	129
11.3 Operational Parameters	129
11.4 Regulatory Compliance	126

# **SECTION 1 INTRODUCTION**

### 1.1 <u>The Coulter Principle</u>

In the late 1940's a technique that allowed particles homogeneously suspended in a conducting liquid to be simultaneously counted and sized was developed and patented by Wallace Coulter. Instruments employing this technique were originally designed to facilitate blood cell analyses but over the years have become invaluable tools in industrial applications as well. The Coulter Principle, also known as the Electrical Sensing Zone (or ESZ) method, is referenced in many national and international standards, and analyses as well as analytical instruments based on the Coulter Principle often serve as references in the evaluation of both other particle size analyzers and sizing techniques.

In the ESZ method a suspension is made to flow through a small cylindrical opening (the aperture) separating two electrodes between which an electric current flows. Although the magnitude of this current may be small, (typically about 1 mA), the resistance, or "pinch", created by the restriction separating the electrodes produces a considerable current density within the aperture. As each particle passes through the aperture (or "sensing zone") it displaces its own volume of conducting liquid, momentarily increasing the impedance of the aperture.

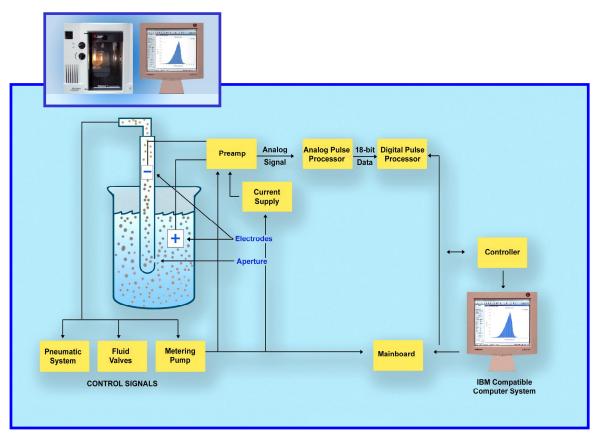


Figure 1.1-1 Diagrammatic illustration of the Coulter Principle applied in the Multisizer 3

This change in impedance produces a tiny but proportional current flow into an amplifier that converts the current fluctuation into a voltage pulse large enough to measure accurately. The Coulter Principle states that the amplitude of this pulse is directly proportional to the volume of the particle that produced it. Scaling these pulse heights in

volume units enables a size spectrum to be acquired and displayed. In addition, if a metering apparatus is used to draw a known volume of the suspension through the aperture, a count of the number of pulses will yield the concentration of particles per unit volume in the suspension.

In industrial applications historically it has been common practice to report particle 'size' in linear terms, and so the volume of a particle is converted into an equivalent spherical diameter (ESD). In this case, the "particle size" is the diameter of a sphere whose volume is equal to that of the particle. The biological and biomedical sciences traditionally (when reporting the results of cellular analyses) are more accustomed to using actual volumetric metric units, and in these applications "particle size" is usually presented in terms such as femtoliters or cubic microns (fL, or  $\mu$ L<sup>3</sup>).

Particle surface areas can be calculated, with the assumption that the surface area is that derived from (and equates to) the surface area of a perfect sphere whose ESD is the same as that of the volume of the particle.

The International Standard ISO 13319 Determination of Particle Size Distributions-Electrical Sensing Zone Methods describes and gives guidance on the measurement of the size distribution of particles using the Coulter Principle.

#### 1.2 <u>About the Multisizer™ 3 COULTER COUNTER®</u>

The BECKMAN COULTER Multisizer 3 is a flexible, multi-channel analyzer employing the Coulter Principle (or Coulter electrical impedance method) together with state-of-theart Digital Pulse Processing (DPP) technology to provide both particle sizing and counting within an overall size range of 0.4 µm to 1200 µm. Data is acquired and processed using proprietary DSP circuitry and may be archived, displayed and reanalyzed on any PC equipped with a Pentium<sup>®</sup> II or similar microprocessor running Beckman Coulter's Multisizer<sup>™</sup>3 software on Microsoft's Windows<sup>®</sup> operating systems versions 95, 98, 2000 or NT 4.0.

The Multisizer 3 uses a unique, mercury-free system to draw suspensions through the aperture at a steady rate and to meter these liquid volumes precisely enough to yield accurate particle concentration measurements. The instrument is also equipped with a continuous flow alternative for applications where accuracy of liquid metering is unimportant or less critical, as well as options that allow the user to define analysis parameters based on elapsed time or particle count.

By using the very latest in digital pulse measurement techniques, an entire analysis (which may consist of many thousands of particle pulses) can be stored either in memory, or on disk, for recall and reprocessing. This technology allows, for the first time, the resolution of an analysis to be increased up to a hundred-fold by narrowing the size range of interest in real time, eliminating the need to re-run the sample. Using the "Convert Pulses to Size" feature in the software, measurements from the lowest to the highest possible resolution can be made on just one pass of sample through the aperture.

Storing the individual values of pulse heights means that it is not necessary to predetermine the number of size classes (also called "channels" or "size bins") into which any size distribution can be divided. This choice is at the discretion of the operator and can be chosen either before or after the analysis or even changed during analysis. Any number of channels from 4 up to 300 may represent a partial or full range of an aperture. A smaller number may improve the appearance of a size distribution by reducing statistical scatter in the results, while details in the results may be enhanced by

increasing the number of size classes in regions of particular interest. A higher number of channels provides more detail about the size of the particles The advanced digital electronics used in the Multisizer 3 allow the number of channels to be changed in real time without losing any of the information acquired with other channel spacing.

Results can also be stored in a fully processed (abridged) form that conserves disk space. Although in this form the raw pulse data is lost and so cannot be re-analyzed at a different resolution, the processed results still can be enlarged, or a region of interest selected, using the "Zoom" facility in the software.

Both number data and size distributions are corrected by proprietary algorithms for the occurrence of multiple particle passage (coincidence) in the aperture. An indicator provides real time particle concentration in the analysis suspension. The Multisizer 3 software includes a pulse monitor window that offers a guide to the electronic behavior of the aperture. The time base of this window can be adjusted to show either pulse concentration or the shapes of individual pulses. Electronic pulse editing facilities ensure that very narrow size distributions can be highly resolved by minimizing the effect of line width broadening in the sensing aperture.

Data can be displayed graphically as number, volume, mass and surface area size distribution plots, or as tables of values versus size. Both display types can show data as differential or cumulative frequencies. Full data reduction facilities are included for statistical and trend analysis, with export and import capabilities to text and spreadsheet programs via disk files or the Windows Clipboard. All results can be printed/plotted on any printer that has an installed driver compatible with the PC's Windows platform.

A report facility allows results to be combined with sample information in a user selectable way to produce a systematic reporting configuration. Report formats can be individually named for ease of archiving and retrieval. Number and date formats are supported in both conventional as well as in ISO 9276-1 forms.

The Multisizer 3 system, illustrated in figure 1.2-1, comprises an integrated fluid handling and pulse processing unit (the Analyzer Module), and a personal computer with an XVGA monitor. (A keyboard, mouse and an ink-jet printer are also part of the system but are not shown in the figure.)

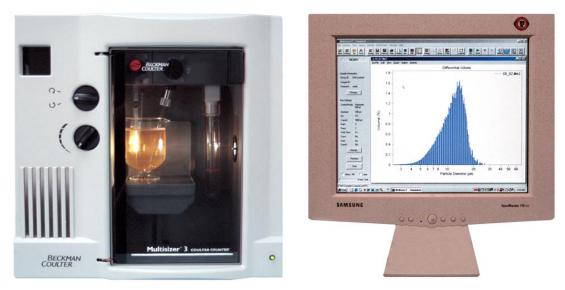


Figure 1.2-1 The Multisizer 3 system

With the pre-amplifier design used by Beckman Coulter in the Multisizer 3, there is no voltage rise. A sensitive oscilloscope attached to the aperture electrodes will detect no pulses at all. Instead, the pre-amplifier offers a momentary parallel path for the current that cannot pass through the aperture. The sum of these two currents, one through the aperture, the other through the amplifier, exactly equals the constant current that flows when no particle is present. The electrical current flowing in the aperture electrode circuit therefore always remains fixed.

A current pulse is converted into a series of voltage pulses of different heights by main amplifiers of different but exactly known gain. By scanning and digitizing all of these pulses at extremely high speed and storing the parameters that are required, it is possible to describe the original pulse. It creates for each pulse a unique set of numerical integers that represent the pulse characteristics to the required precision.

In the Multisizer 3, these numbers are large enough to describe any particle pulse within the aperture measurement range to better than 1% accuracy. All pulse parameters are recorded as a file of data in the order in which they arrive (list mode data). If for any reason, pulse data arrive quicker than they can be stored, the values are held temporarily in a buffer that continues to be emptied to the main file after an analysis has been completed. This updating process ensures that no data is lost except for the statistical losses due to coincident particle arrival within the aperture's sensing zone.

Because pulses are digitized to maximum resolution, the resolution of the Multisizer 3 can be increased or decreased before, during or after an analysis. The number of "channels" or "size classes" into which the data is divided (as well as how these intervals are spaced over the aperture range) represents the extent of the instrument's resolution. The same size distribution can be shown as 100 channels of results, 256 channels, or any other number from as few as 3 to as many as 300 size classes over any aperture's working range as required by the operator. All measurements, however presented, actually contain the maximum resolution and can be saved with this high degree of precision, though this will increase the size of the data file significantly. Using a setting of 256 channels (as opposed to the maximum the software will allow) reduces the statistical scatter (white noise) and improves the appearance of the results. If the number of particles measured is not sufficient to provide a size distribution without significant scatter the Multisizer 3 allows the analysis to be continued without losing previous data, even when the continuation is at a different level of resolution than the initial result.

# 1.3 About this Manual

This manual is intended to provide the user with all the information needed to operate and maintain the Multisizer 3 system safely and effectively within the stated working range and conditions given in <u>Section 11</u>.

### 1.3.1 Conventions

 The fluid handling and pulse processing unit of the Multisizer 3 system is referred to as the "Analyzer Module". The PC, keyboard, monitor and printer are all referred to collectively as "the computer".

- C means "click the mouse button". Unless the mouse has been inverted for lefthanded users, this will be the left mouse button.
- ①① means "double click the (left) mouse button".
- 🖾, 🗹 and ⊙ mean "check the box or radio button", and and O mean "uncheck the box or button" or leave unchecked.

#### 1.3.2 Warnings and Cautions

Indicates a situation or condition that might result in personal injury if not corrected.

Indicates a situation or condition that, if ignored or proceeded with, may result in damage to the instrument or serious error in the results produced by it.

(Help note) contains hints and useful information to help optimize results quality and ease of use.

#### 1.4 <u>Warnings</u>

(STOP)

#### 1.4.1 Electrical

- (1) High voltages are present inside the instrument. Always disconnect the instrument from the power supply before removing the cover.
- (2) The instrument must be grounded correctly.

#### 1.4.2 Chemical

- (1) Do not use any electrolyte solutions or diluents that are not compatible with the materials of construction. Consult Beckman Coulter Inc. or its local representative before using any concentrated acids or alkalis in the instrument.
- (2) Toxicity, safety, and proper handling procedures for diluents and reagents used in particle and cell analysis should be adhered to at all times. Consult appropriate safety manuals and Material Safety Data Sheets for the items.
- (3) Care should be taken when mixing or exchanging diluents. Reactions can occur between incompatible solvents and electrolyte salts, and may be violent.
- (4) Do not use azide and cyanide salts in acid solutions.
- (5) Flammable electrolyte solutions should be prepared for use in an appropriate environment and brought to the instrument only when required for an analysis.

#### 1.4.3 Mechanical

- (1) Do not interfere with or attempt to disable the interlocks incorporated in the door and beaker platform assembly.
- (2) Do not use force to remove or replace any glass item. In the event of difficulty, consult Beckman Coulter Inc. or its local representative.
- (3) Replace an aperture tube empty of diluent. Do not attempt to pre-fill a tube before locating it in the aperture tube connector.

## 1.4.4 Fire

- (1) Many non-aqueous electrolyte solutions are flammable and contain organic salts that can support or enhance fire. Where possible choose less flammable alternatives.
- (2) Prepare and filter flammable solvents and diluents away from the Multisizer 3 itself, preferably under a suitable fume hood. Extinguish all open flames in the vicinity and observe normal precautions against ignition by hot surfaces or electrical sparks.

## 1.5 <u>Cautions</u>

### 1.5.1 Electrical

- (1) Use the Windows<sup>®</sup> Shut Down command to turn the computer off before switching off power. Switching off power first will generate a Windows error message at next computer start up.
- (2) Power must be largely free from interference. The Multisizer 3 is CE, UL and CSA compliant with regards to electrical interference but some circumstances may benefit from a constant voltage transformer or regulator.

## 1.5.2 Chemical

- (1) If Beckman Coulter ISOTON<sup>®</sup> II diluent is used, the Multisizer 3 must be drained and flushed through with distilled water before carrying out a bleach disinfection procedure.
- (2) If the Multisizer 3 has been subjected to a bleach (sodium hypochlorite solution) disinfection procedure, the instrument must be drained and flushed through with distilled water before refilling with ISOTON® II.
- (3) Take care disconnecting diluent lines -- open-ended tubing may allow liquid to siphon out of containers.
- (4) Never stand containers of liquids on top of the Multisizer 3. Repair of instruments damaged or affected by spilt liquids will not be covered by any warranty or guarantee.

### 1.6 Warming up

As with all sensitive electronic instruments, the Multisizer 3 components offer best performance once they have reached a steady working temperature. This may typically take 15 minutes from power on.

### 1.7 <u>Sources of Error</u>

### 1.7.1 Calibration constant

Size distributions are only correct if the calibration constant (Kd) used to construct the size scale is correct. Instructions on how to calibrate an aperture are contained in <u>Section 5</u>. If a different constant is entered manually, the software will use the manually entered value despite the fact that this may result in an incorrect calculation of the size range appropriate for that particular aperture.

#### 1.7.2 Damaged apertures

Aperture damage, such as a cracked wafer, can exhibit a variety of symptoms. Usually, it results in sporadic bursts of high counts at the lower aperture size limit. It is often the cause of deterioration in blank count in a tube previously giving quiet and reliable blank counts.

#### 1.7.3 Deposits on and around aperture wafers

Protein and others biological materials may create a deposit on apertures that could alter their calibration and background count characteristics. Apertures used for these applications should be subjected to appropriate cleaning procedures at frequent intervals. It is advisable to perform a cleaning procedure to avoid any material deposit or microbial growth inside the system (see <u>Section 6.10</u>).

#### 1.7.4 External sound and mechanical vibration

Loud sounds and vibrations may register in the aperture as measurable pulses.

#### **1.7.5** Electrical interference

Electrical interference both from the power line and radiated may, if outside the scope of EMC compliance requirements, produce erroneous measurements.

#### **1.7.6** Particles behind the aperture

In the case of either very large or unusually dense materials, particles may settle inside the aperture tube when the instrument is standing idle. Sediments of such particles have been known to swirl up behind the aperture when the instrument is then used and cause interference with subsequent counts and size distributions.

#### 1.7.7 Sample dispersion

Measurements of powder dispersions are only as good as the dispersed state of the sample and the sampling procedure allows. Incorrectly dispersed or unstable dispersions will give rise to incorrect results.

#### 1.7.8 Excessive aperture current

Excessive current causes the electrolyte solution in the aperture to heat, or even to boil, causing noise (see <u>Section 4.4.9</u>).

**Back to TOC** 

# SECTION 2 INSTALLATION

# 2.1 Location

Ideally the instrument should be located in a dust free room. Dusty environments should be avoided to minimize sample contamination. The instrument should be placed on a bench that is not subject to:

- (1) Excessive airborne dust
- (2) Strong vibrations
- (3) Spillage or splashing
- (4) Sounds of high intensity (e.g., ball mills, ultrasonic baths or probes); electrical noise (e.g., electrical switching gear, water baths or gas chromatographs)

## 2.2 Unpacking

The Multisizer 3 will normally be installed only by Beckman Coulter approved and trained installation engineers. Unless otherwise agreed to by Beckman Coulter, **DO NOT UNPACK** the Multisizer 3.

The Multisizer 3 is a heavy instrument. Follow the instructions below. Failure to do so may result in personal injury and/or damage to the instrument. Such damage may not be covered by warranty.

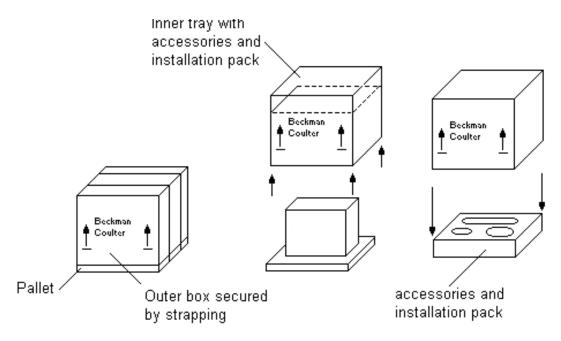


Figure 2.2-1 Unpacking the Multisizer 3 Analyzer Module

- (1) Lift the Analyzer Module (requires two people) off the pallet and onto the bench taking care not to exert any force on or lift by the plastic front compartment molding (see Figure 2.2-1).
- (2) Turn the outer box over to gain access to the Accessories and Installation Pack fastened to the inside top of the outer box. Slide out the Pack and inspect all installation items for damage. Compare the contents with the Packing List.

(3) Report any missing items or damage to any part of the instrument to Beckman Coulter or their local distributor, and to the carrier responsible for delivery. Do not send back any items before receiving instructions to do so from the manufacturer.


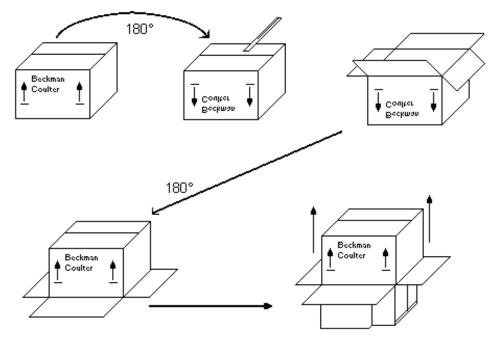


Figure 2.2-2 Unpacking the computer monitor

- (4) Open the carton containing the computer monitor and remove the monitor as shown in figure 2.2-2 by inverting the carton and slitting the tape along the box bottom flaps. Turn the carton back upright and slide it upward from the monitor.
- (5) Unpack the computer and printer by placing upright on the bench and slitting the tape around all box flaps. Lift the items out from the boxes. Compare the computer and printer manufacturers' packing lists and report any deficiencies or damage as in (3) above.

### 2.3 <u>Connections</u>

- 1. Remove the tubing and containers kit from its package.
- 2. Prepare the lids for each container. The waste lid uses one curve tube to connect the container to the waste outlet of the instrument and has only one connection cable. The electrolyte lid uses two tubes and it has a connector cable and ground cable. The straight tube, the intake tube, goes from the inner side of the lid to the bottom of the electrolyte container. The curve tube connects the container to the electrolyte intake on the side of the instrument (See Figure 2.3.1 and 2.3.2).
- 3. Leave the waste container empty and screw the lid with the sensor and tube. Fill the electrolyte container with the required electrolyte and screw the lid down.
- 4. Attach the appropriate tube from each container to the fittings marked "Diluentl" (Electrolyte) and "Waste" at the top of the left side panel (See Figure 2.3.2).

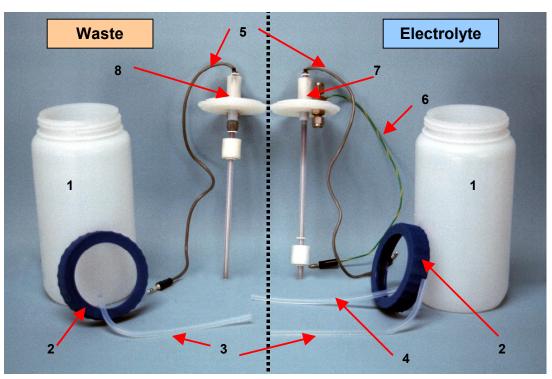

- 1. Containers
- 2. Lids
- 3. Connecting Tubes
- Electrolyte Intake Tubing
   Connecting Cables
- 6. Ground Cable
- 7. Electrolyte Sensor
   8. Waste Sensor



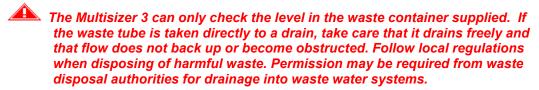


Figure 2.3.1 Assembly of the Electrolyte/Waste System



Figure 2.3-2 Left side panel showing waste and diluent connections

5. Plug the connector cables from the Waste and Diluent sensors into the dual sockets immediately behind the pipe fittings on the left-hand side of the instrument. Either cable may go into either socket. Plug the ground cable from the electrolyte sensor into the ground socket (See Figure 2.3.2).



- 6. Use the EtherNet cable supplied with the instrument to connect the Multisizer 3 to the PC's network card. (See Figure 2.3.3). The parallel and serial connectors are intended to be used by Beckman Coulter authorized service personnel.
- 7. Check that the voltage selection indicator located on the rear side of the instrument is set for the voltage (line voltage) to which the instrument is to be connected. (see Figure 2.3-4).


All items comprising the Multisizer 3 system must be set for the correct voltage. Connecting the instrument to an incorrect voltage may cause damage to the instrument and invalidate any warranty. Different voltages require appropriate fuses.

- 8. If the selector does not indicate the correct voltage setting, change it as follows: Remove the fuse holder and cover plate by inserting a small bladed screwdriver (~3 mm blade width) into the recess at the top (figure 2.3-4). Lever upwards to "pop out" the holder. Using needle nosed pliers, pull out the voltage selector board. Do not pull on the plastic voltage indicator pin. Voltage values are marked around the four corners of the plate (Figure 2.3-4). Holding the board as shown in Figure 2.3-4B, pull out the plastic pin as far as it will go (Figure 2.3-4C) and rotate it so that it points outwards towards the cover plate. Figure 2.3-4D shows the orientation of the pin for 120 V.
  - (a) Push the pin back in towards the board (the pin will locate in a recess unique to the chosen voltage). Turn the board until the voltage required faces inwards (towards the instrument). Replace the selector board and push firmly into place.
  - (b) If fuses need replacing use only 250 V, 3.15 A, "T" (time delay), IEC standard 20 mm fuses.

(c) Fit the cover plate so that the appropriate small hole at its base fits over the white plastic indicator pin and push firmly at the top until it clicks into its seating. The white pin should now indicate the correct voltage for the supply. The positions left to right are: (See Figure 2.3-3).



- 9. Make sure all the power switches of the Multisizer 3, the computer, and the printer are set to O (=OFF) then connect these items to power with the cables supplied.
- 10. First switch the Multisizer 3 analyzer unit power switch to | (=ON), followed by the computer and printer. To find out how to use the instrument's control program turn to <u>Section 3.3.1</u>

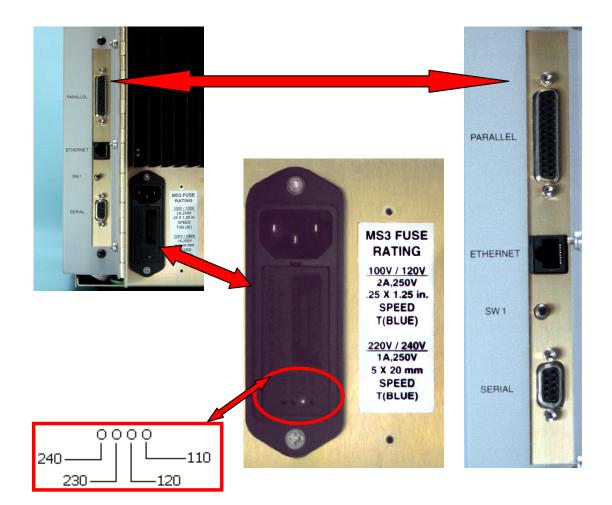

**Figure 2.3-3** Detail from the bottom left side of the rear panel showing the connections to the PC and the power socket with the fuse holder

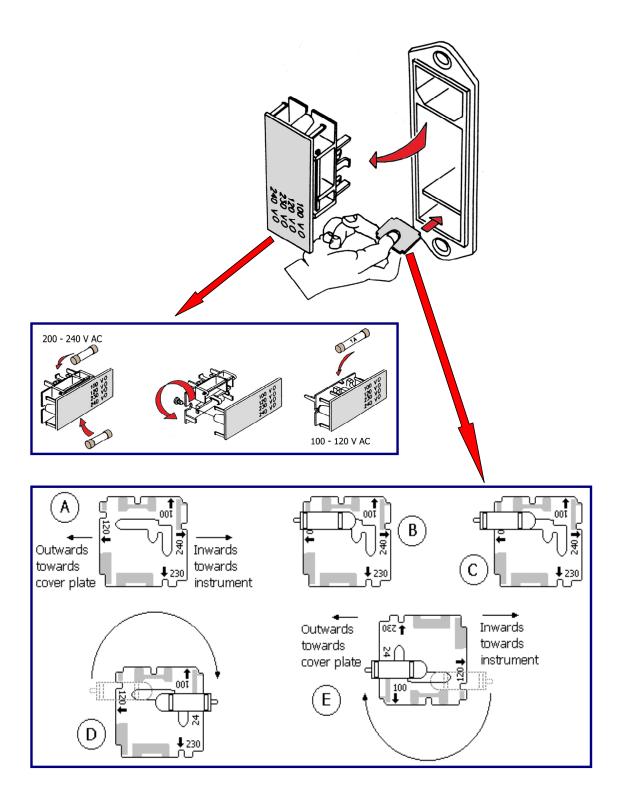


 Figure 2.3-4 The voltage selector board: B shows the orientation for 240 V; C & D show

 the conversion from 240 V to 120 V, and E the final orientation for 120 V

 before re-insertion.

 Back to TOC

# **SECTION 3 BASIC OPERATIONS**

The software controls all the operations on the Multisizer 3. Only the stirrer speed and the sample platform need to be adjusted manually on the analyzer unit.

#### 3.1 Starting the Instrument and its Control Program

1. If the analytical system has been started before the program has been launched, the first view presented by the program is the copyright acknowledgement box (Figure 3.1-1). If the checkbox is unchecked, the program will behave purely as a data processing program and the Multisizer 3 control menu options will be either "greyed out" or not displayed.

Beckman Cou	lter Multisizer 3	×
•: <b>U</b> I	Version 3.33 23 Jul 2001	
Сору	vright© 1990-2001 Beckman Co	oulter, Inc.
	BECK	ER
J PA	RTICLE CHARACTERIZ	ATION
19:20:41	Eastern Daylight Time	31 Jul 2001
Connect I	to Multisizer 3	OK
	Configured for no security	

Figure 3.1-1 Copyright Acknowledgement Box

2. Check the box **Connect to Multisizer 3**, *A*→**OK** to launch the control program.

#### ✓ ☺ If the analyzer unit has not already been switched on, do not check the "Connect to Multisizer 3" box. Connection can be made later under "Run" in the Main Menu Bar.

- 3. Once connection has been made, a series of status messages illustrate the initializing process of the Analyzer Module. These normally start with "Waiting" and/or "Initializing valves" but the precise sequence will depend on the actual start condition of the Analytical Module. After initializing the complete system, the status message will change to "Ready" (see Figure 3.1-2).
- 4. The Security Set Up dialog box is shown the first time that the program is started. At this moment the security level can be selected. These security options can also be selected at a later time (see <u>Section 10.7.1</u>). It is advisable to read <u>Section 10</u> before selecting the security level.
- 5. To set up the security options at a later time, *<sup>•</sup>*⊕→Cancel in the Security Set Up dialog box (see Figure 3.1-3). Refer to Section 10 to set up security at a later time.
- To select the security level at this time choose the required level option (see Figure 3.1-3). <sup>A</sup>→OK and go to <u>Section 10</u>.

7. Disconnection from the Multisizer 3 can be made at any time by *A*→**Run** then *A*→**Disconnect from Multisizer 3**. (When disconnection has been made this way the fact is "remembered" when the program is closed down. The checkbox in figure above will be then be unchecked next time the program is launched.)

Ele Window Run System Settings Preferences Security Help
Image: Construction         Image: Construction
READY         Sample Information         Gioup ID: 22         Sample ID:         Comment:         Dhange.         SOM         Control Mode: Volumentic         SoupLine         SoupLine         Aperture:         100 µA         Gain:         2         Ruma:         100 µA         Gain:         2         Preview         Statt         Preview         Statt         Vest Tank

Figure 3.1-2 Screen view when is "Ready".

Figure 3.1-3 Security Set Up dialog box

## 3.2 Instrument Controls

1. Control of the Multisizer 3 is essentially accomplished via the drop-down menus enabled when a connection to the Analyzer Module has been established. Without a connection, the Main Menu bar defaults to the standard menu set for data review,

editing, and printing. In this case the control menus are no longer visible. (The Multisizer 3 software, because of the digitization of pulse data, has the ability to reprocess previously acquired results, as long as the data was stored with the option to include the pulse data selected.)

Multisizer 3	- Simulated			
<u>F</u> ile <u>W</u> indow	<u>R</u> un S <u>y</u> stem	<u>S</u> ettings	Preferences	Sec <u>u</u> rity <u>H</u> elp
Open Overlay	L A Up Fonts	Overlag	Save Save as	Print Preview Info
RE	ADY	1		
, ,				
Sample Inform	ation	-		
Group ID:	D54 Control2			
Sample ID:				
	water			
	ange			
Run Settings-		-		
Control Mode	: Volumetric			
	500 µL			
Aperture:	100 µm			
Kd:	111			
Current:	1600 μA			
Gain:	2			
Runs:	1			
Wait Time:	0 s			
Save:	No			
Print:	No No			
Export:	NO			
Cha	ange			
Pro	eview			
	Start			
<u></u>				
🗖 Stirrer: O	ff 🗌 Lamj	5 E		
	Waste Ta	nk		

Figure 3.2-1 The Status Panel

- 2. Provided connection to the Analyzer Module has already been established, stopping and starting an analysis is controlled by "Buttons" on the Status Panel shown on the left side of the Control Page. These buttons change their function and annotation according to the status of the analysis.
- **3.** The instrument can also be directly controlled via the drop-down menus, for example:

😈 Multisizer 3 - Simulated							
<u>F</u> ile <u>W</u> indow	<u>R</u> un System	<u>S</u> ettings	<u>P</u> references	Sec <u>u</u> rity <u>H</u>			
Open Overlay	<u>S</u> tart <u>P</u> ause Stop		F2 F4 F3	Preview			
R			F1 F5	0			
Sample Inforr	Reset		F8 F9				
Group ID:	Change Ap	erture Tub	e <u>W</u> izard				
Sample ID: Comment:	<u>C</u> alibrate Ap ⊻erify Apert		tion				
Cł	Disconnect	from Multi	sizer <u>3</u>				
- Run Settings Control Mode	e: Volumetric 500 μL						

Figure F3.2-2 The Run Menu

#### 3.3 Shutting down the Multisizer 3

At any time, the software program can be closed down by any of the standard Windows applications closing actions such as:

**A**→**File** on the Main Menu Bar then **A**→**Close**,

or  $\neg \ominus \rightarrow \boxtimes$  on the Window top right-hand corner.

- (1) When the program has been closed, the Analyzer Module power switch can be switched off.
- (2) Closing down the program leaves the Multisizer 3 connection in its last state. If connection has been enabled, the copyright acknowledgement box will show "Connect to Multisizer 3" checked next time the program is started. If the connection has already been disabled through "Run", the copyright acknowledgement box will show the "Connect to Multisizer 3" box unchecked.


#### **Back to TOC**

# **SECTION 4 GETTING STARTED**

# 4.1 Locating and Identifying Controls and Parts You Need to Know



# 4.1.1 The Frontal Panel of the Multisizer 3

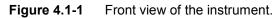

On the front panel of the instrument you will find:

- Aperture Viewer: This screen allows visual monitoring of the aperture.
- Stirrer Rotation Control: This knob changes the direction the stirrer rotates.
- Stirrer Speed Control: This knob is used to set the rotation speed of the stirrer and to turn it on and off. The stirrer can also be turned on or off from the Multisizer 3 software.

# 4.1.2 The Sample Compartment

The sample compartment is enclosed in a Faraday Cage to avoid electric interference. Opening the glass door of the sample compartment you can find the Sample Platform and the places where the electrode and stirrer have to be attached (See Figures 4.1.2 and 4.1-3).


Inside the Sample Compartment you will find:

- Electrode
- Stirrer

•

- Aperture Tube
- Sample Platform

- **Trap:** Its function is to prevent large and dense particles to go into the system. When the sample sediment reaches the red mark it is necessary to clean the trap.
- Aperture Tube Knob: This knob is used to hold or release the Aperture Tube.
- Stirrer Knob: This knob is used to fix the position of the stirrer.

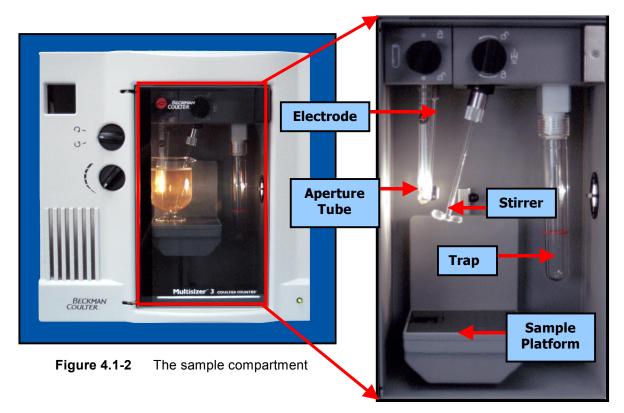

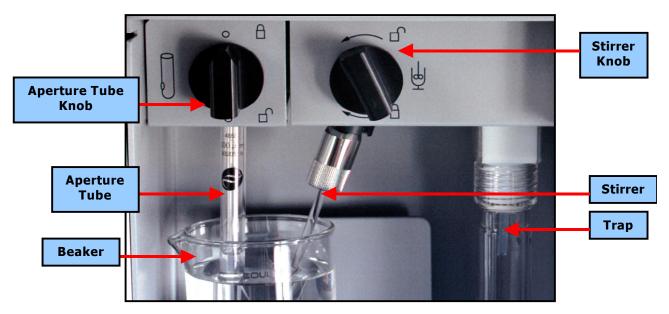


Figure 4.1-3 Inside the sample compartment

# 4.2 Fitting the Electrode and Stirrer

- 1. Carefully remove the electrode and stirrer from their packages.
- 2. Hold the electrode only by its metal connector and insert it into the electrode socket located on the upper side of the sample compartment. Rotate it until the platinum spade is positioned roughly parallel with the lamp lens axis behind the aperture tube then push up to complete the connection. (See Figure 4.2-1).

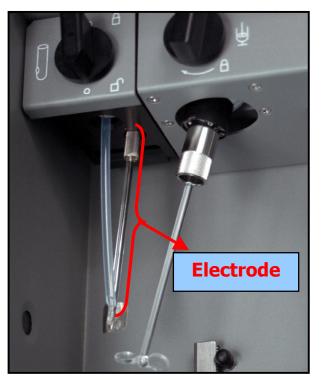


Figure 4.2-1 Electrode.

The electrode stem is made of glass and is therefore fragile. If the connector does not enter its socket easily, do no use undue force. Do not twist, push or pull the glass stem in any way. Hold the electrode only by the connector.

The wire in a broken electrode which may appear to have electrical continuity but the break may still cause erroneous results.

3. Hold the stirrer, paddle downwards, about 2 cm from the tapered top, and push firmly into the collet of the stirrer motor as far as it will go. It helps to support the collet with the finger and thumb of the other hand. The black elastomer band around the collet will provide enough securing tension to hold the stirrer in place (see <u>Figure 4.2-2</u>).



If the stem does not enter the collet easily, do not use undue force. Do not twist or flex the glass stem in any way. Hold only as described.

4. Adjust the vertical position control of the stirrer motor until the stirrer paddles sweep at or just below the bottom of the aperture tube. Test the clearance



available to the paddles by rotating the collet between finger and thumb. The stirrer must not encounter the electrode or the aperture tube when it rotates.

Figure 4.2-2 Stirrer.



Do not hold the stirrer by its paddles or push or pull on them to provide additional leverage.

Glass stirrers are fragile and knocking them against beaker walls or subjecting them to ultrasound from probes and baths easily damages the paddles. Multisizer 3 glass stirrers cannot be used with very small beakers (e.g., less than 150 mL). A small helical agitator can often be used with beakers down to 50 mL squat form. Neither can be used with ACCUVETTE® Il vials.

#### 4.3 <u>Fitting the Aperture Tube</u>

The first time you use the instrument familiarize yourself with the operations for fitting and changing an aperture tube. You should practice several times until you get used to performing this routine.

1. Wash all items in filtered liquid.

For all aqueous diluents, well-filtered distilled quality (DQ) water will be found suitable for washing and rinsing items. For most non-aqueous electrolyte solutions, filtered propan-2-ol (iso-propanol, I.P.A.) is a useful rinsing liquid. When using electrolyte solutions that are incompatible with solvents containing -OH groups, filtered butan-2-one (also known as methyl ethyl ketone} or a low boiling point petroleum can be used instead of I.P.A.

- 2. <u>Figure 4.1-1</u> shows a frontal view of the instrument. Open the sample compartment door. The trigger release for the beaker platform is located under the front lip of the platform itself. This trigger is designed to be squeezed while supporting the platform with fingers and thumb. Move the platform down until it is at its lowest position.
- 3. Rotate the aperture tube clamp counter-clockwise to the "open seal" position. Insert the aperture tube in the connection head as far as it will go with the logo facing the front towards you. **DO NOT GREASE** the aperture tube. The Multisizer 3 aperture tube connection is sealed by O-rings and forms a greaseless seal.
- Tighten the Aperture Tube Knob clockwise to the closed (tighten seal) position (see Figure 4.3-1).

There is no need to over tighten the knob to obtain a leak proof tight seal. A correct seal is obtained when the aperture tube cannot easily be rotated or pulled down when gripped by fingertips alone. Using just enough lock on the aperture tube will prevent excessive wear and tear on the O-ring seal and extend its life.

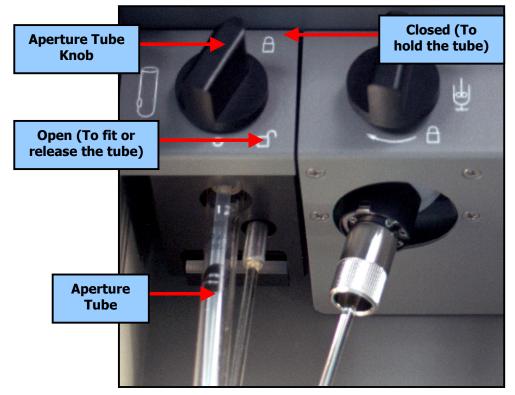


Figure 4.3-1 Aperture Tube

### 4.4 <u>Setting Up the Aperture Tube</u>

Before setting up the new aperture tube, the instrument must be in the following condition:

- (1) The electrolyte container is sufficiently filled (enough to raise the level sensor).
- (2) The Waste container is empty (enough to lower the level sensor).
- (3) The Analyzer Module and the computer have been powered on with the control program running.

📴 Multisizer 🕻	3 - Advanced Operato	or mode – Simulato	ed	
File Window	Run System Settings	Preferences Secu	urity Help	
Open Overlay	Start Pause Stop Cancel Continue Resume Reset	F2 F4 F3 F10 F5	Preview Info	Edit Graph Listing
Sample Infor	Preview	F9		
Group ID	Change Aperture Tub	e Wizard	$\triangleright$	
Sample ID: Comment:	Calibrate Aperture Verify Aperture Calibr	ation		
Cł	Disconnect from Multi	sizer 3		


On the Main Menu Bar, ⊕→Run, then, ⊕→Change Aperture Wizard.

The wizard will guide you through the process.

# 1. Removing the currently installed aperture tube.

Rotate the aperture tube clamp counter-clockwise to the "open seal" position and remove the aperture tube (see Figure 4.2-2)  $\bigcirc \rightarrow$  Next

Change Aperture Tube Wiza < Back Next >				
Remove the currently installed aperture tube				
C Install a new aperture tub	е			
C Place a beaker of clean e	electrolyte o	n the p	olatform	
C Close the door				
${f C}$ Select the electrolyte	Select the electrolyte		lectrolyte	
C Select the new aperture t	Select the new aperture tube		erture Tube	
${f C}$ Fill the system	C Fill the system		ill System	
${f C}$ Adjust the metering pump		Met	ering Pump	
C Set current and gain		Cur	rent & Gain	
C Measure noise level		Mea	asure Noise	
C Calibration	Verify		Calibrate	
C All done			Done	
C	lose			

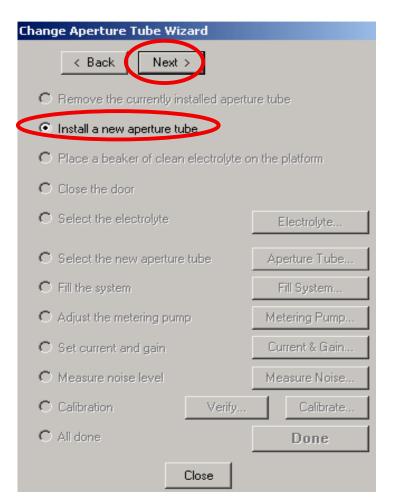

# 2. Installing a New Aperture Tube.

From the table below, select an aperture tube appropriate to the sample to be analyzed.

Aperture Size (µm)	Analysis Range (µm)
20	0.4 - 12.0
30	0.6 - 18.0
50	1.0 - 30.0
70	1.4 - 42.0
100	2.0 - 60.0
140	2.8 - 84.0
200	4.0 - 120.0
280	5.6 – 168.0
400	8.0 - 240.0
560	11.2 – 336.0
1000	20.0 - 600.0 •
2000	40.0 - 1200.0♦

• Range depends on sample density

Install the selected tube following the procedure described in Section 4.2.  $\textcircled{} \rightarrow \mathbf{Next}$ .

	Change Aperture Tube Wizard	
<ol><li>Placing a beaker of clean electrolyte or the platform.</li></ol>	< Back Next >	
After placing the beaker with clean electrolyte on the platform, fill the electrolyte container	<ul> <li>Hemove the currently installed apert.</li> </ul>	ire tube
with the same kind of electrolyte being used in the beaker.	<ul> <li>Install e cerv operture tube</li> <li>Place a beaker of clean electrolyte o</li> </ul>	n the platform
	C Close the door	In the platform
	C Select the electrolyte ISOTON II	Electrolyte
	C Select the new aperture tube	Aperture Tube
	C Fill the system	Fill System
	C Adjust the metering pump	Metering Pump
	${f C}$ Set current and gain	Current & Giain
	C Measure noise level	Measure Noise
	C Calibration Verify	Calibrate
	C All done	Done
	Close	
ange Aperture fube Wizard		
Back Next >		

Change Aperture Tube Wizard			
< Back Next >			
C Remove the currently installed apertu	ire tube		
C Install a new aperture tube			
C Place a beaker of clean electrolyte o	n the platform		
Close the door			
C Select the electrolyte ISOTON II	Electrolyte		
C Select the new aperture tube	Aperture Tube		
C Select the new operate table	Apertare Tabe		
C Fill the system	Fill System		
C Adjust the metering pump	Metering Pump		
C Set current and gain	Current & Gain		
C Measure noise level	Measure Noise		
C Calibration Verify	Calibrate		
C All done	Done		
Close			


## 4. Closing the Sample Compartment door.

Raise the platform with the beaker to the appropriate height and close the door.  ${}^{\circ} \rightarrow$ Next.

# 5. Selecting the electrolyte.

 $\mathcal{A} \rightarrow \mathsf{Electrolyte}$ . From the drop down list select the electrolyte to be used. If the electrolyte is not on the list, type the name in the box. You can also select the dispersant to be used, or type its name if it is not on the list. After selecting the electrolyte and dispersant,  $\mathcal{D} \rightarrow OK$ . Be Sure to have the beaker and electrolyte container filled with the same electrolyte selected


Change Aperture Tube Wizard		
< Back Next >		
C Remove the currently installed apertu	re tube	Select Electrolyte and Dispersant
${f C}$ install a new aperture tube		Electrolyte: SOTON II
C Place a beaker of clean electrolyte or	n the platform	
O Physicae and a start of the		Dispersant: ISOTON II - Glycerol 6:4 0.9% Sodium Chloride in Water
		4% Sodium Chloride in Water 2% Ammonium Thiocyanate in Isopropyl Alcohol
<ul> <li>Select the electrolyte</li> <li>ISOTON II</li> </ul>	Electrolyte	5% Lithium Chloride in Methyl Alcohol 4% Zinc Chloride in Acetone
C Select the new aperture tube	Aperture Tube	5% Lithium Chloride in Acetone 4% Lithium Iodide in Acetone
${f C}$ Fill the system	Fill System	4% Ammonium Thiocyanate in Ethanol
${f C}$ Adjust the metering pump	Metering Pump	
C Set current and gain	Current & Gain	
C Measure noise level	Measure Noise	
C Calibration Verify	Calibrate	
C All done	Done	
Close		

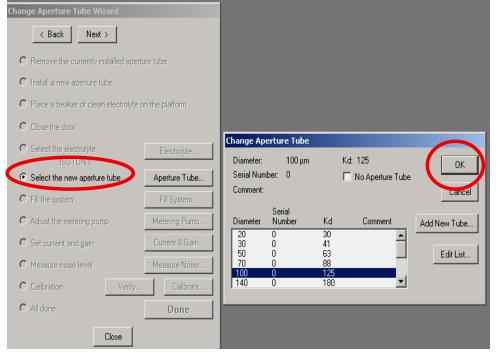

# 6. Selecting the aperture tube.

7. Filling the System.

 <sup>^</sup>⊕→ Fill System

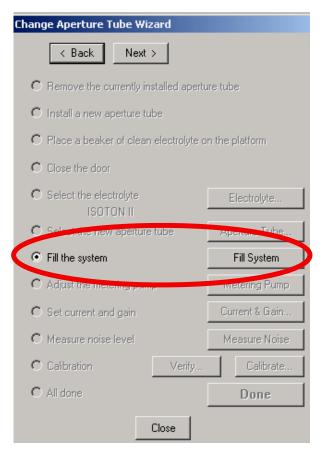

ange Aperture Tube Wizard		
< Back Next >		
C Remove the currently installed aperture tube		
C Install a new aperture tube		
C Place a beaker of clean electrolyte or	n the platform	
C Close the door		
C Select the electrolyte ISOTON II	Electrolyte	
C Select the new aperture tube	Aperture Tube	
C Fill the system	Fill System	
<ul> <li>Adjust the metering pump</li> </ul>	Metering Pump	
C Set current and gain	Current & Gain	
C Measure noise level	Measure Noise	
C Calibration Verify	Calibrate	
C All done	Done	

Adjusting the Metering Pump.
 Ad→ Metering Pump

## 9. Setting Current and Gain.

 $^{+}$ → Current & Gain,  $^{+}$ → Auto Set. When the instrument is done setting the current and gain,  $^{+}$ → OK. If you are analyzing cells,  $^{+}$ → Cells, and  $^{+}$ → OK.

Change Aperture Tube Wizard	
< Back Next >	
C Remove the currently installed aperture tube	
C Install a new aperture tube	
${f C}$ Place a beaker of clean electrolyte on the platform	Current and Gain Settings
C Close the door	Aperture Current: 1600 µA. OK
C Select the electrolyte Electrolyte	Preamp Gain
C Select the new aperture tube Aperture Tube	
C Fill the system Fill System	C 4 Auto-Set.
O Admittle metering pump metoring Pump	
Set current and gain     Current & Gain	C 32
C Measure noise level Measure Noise	
C Calibration Verify Calibrate	
C All done Done	
Close	
	Automatic Current and Gain 🛛 🔀
	Flow Rate: 38.4 µL/s
	Resistance: 14.6 k ohms
	Current Gain
	Particles: 1600 μA 2
	Maximum Cell⊻olume: 0 μm³
	Cancel OK

If cells are to be analyzed, it is very important to select cells in this step. Not selecting cells could cause their damage due to a higher current setting.

### 10. Measuring Noise Level.


A→ Measure Noise.

Change Aperture Tube Wizard				
< Back Next >				
C Remove the currently installed apertu	C Remove the currently installed aperture tube			
C Install a new aperture tube				
C Place a beaker of clean electrolyte o	C Place a beaker of clean electrolyte on the platform			
C Close the door				
C Select the electrolyte ISOTON II	Electrolyte			
${f C}$ Select the new aperture tube	Aperture Tube			
$igodoldsymbol{C}$ Fill the system	Fill System			
${f C}$ Adjust the metering pump	Metering Pump			
C Secourrent and gain	const.º. Gain			
Measure noise level	Measure Nois			
C Calibration Verify	Calibrate			
C All done	Done			
Close				

# 11. Calibration and Verification of the Calibration

If the installed aperture is going to be used for first time, it needs to be calibrated. Now is the time to do this. For a previously calibrated aperture, the verification of the calibration can be done at this time. For details on the calibration and verification procedures, refer to <u>Section 5</u>. If calibration is not necessary,  $\textcircled{} \rightarrow Next, \textcircled{} \rightarrow Done$ 

hange Aperture Tube Wiza	rd		
< Back Next >			
C Remove the currently installed aperture tube			
C Install a new aperture tub	8		
C Place a beaker of clean e	lectrolyte o	n the platform	
C Close the door			
C Select the electrolyte		Electrolyte	
C Select the new aperture t	ube	Aperture Tube	
C Fill the system		Fill System	
C Adjust the metering pump		Metering Pump	
C Set current and gain		Current & Gain	
C Measure noise level		Measure Noise	
C Calibration	Verify	Calibration	
All done		Done	
	ose	$\overline{}$	

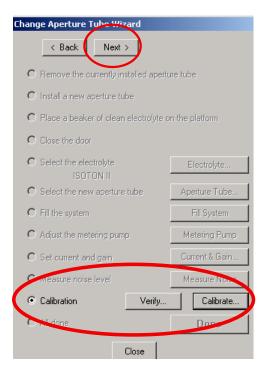

#### 4.5 Aligning the Optics

The three knobs on the side of the instrument are used to align and focus the aperture image in the display window. The two bottom knobs are labeled with a horizontal and a vertical arrow. The horizontal adjusting knob causes the aperture image to move left and right on the display window. The vertical adjusting knob causes the aperture image to move up and down on the display window. The top knob labeled with an "eye", focuses the image on the display window. Adjust the three knobs until the aperture image is centered on the Aperture Viewer (see Figure 4.5.1).

Re-adjust the focus to obtain a well contrasted aperture image. If good contrast is impossible to obtain, it may be necessary to reposition the beaker slightly, particularly if the beaker has a flat bottom and tall form.

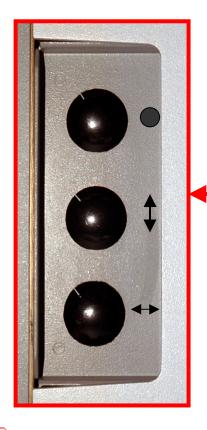


Figure 4.5.1Optics Alignment Knobs



Remember that the image in a microscope is always upside down and horizontally reversed - it always moves the opposite way to the adjustment.

Back to TOC

## SECTION 5 CALIBRATION AND VERIFICATION OF CALIBRATION

### 5.1 <u>Choosing the Calibrator</u>

After installing the aperture as described in <u>Section 4</u>, choose the appropriate calibrator size according to the installed aperture. The calibrator must to be standardize for the **modal value and provide its value**. Beckman Coulter offers a variety of calibrator suitable for the Multisizer 3.

Aperture Size (μm)	Calibrator Nominal Modal Size (μm)	Recommended Beckman Coulter CC Calibrator Nominal Modal Size (µm)
20	2 – 4	2
30	3 – 6	3
50	5 – 10	5
70	7 – 14	10
100	10 – 20	10
140	14 – 28	20
200	20 – 40	30
280	28 – 56	30
400	40 - 80	43
560	56 –128	65
1000	60 – 150	90
2000	90 – 260	90

**Table 5.1-1** Calibrators and size range for the calibration run


# Beckman Coulter is not responsible for erroneous results if the proper calibrator and calibration procedure is not used.

#### 5.2 <u>Preparing the Calibrator</u>


1 Place an ACCUVETTE<sup>®</sup> or a beaker<sup>\*</sup> with the electrolyte and a few drops of the calibrator into the analyzer unit.

\*An ACCUVETTE<sup>®</sup> (20 mL) is appropriate for apertures up to  $100\mu$ m, for larger apertures a larger beaker should be used.

2 On the Status Panel
 <sup>A</sup>→Preview, or in the Main Menu, <sup>A</sup>→Run,<sup>A</sup>→Preview.

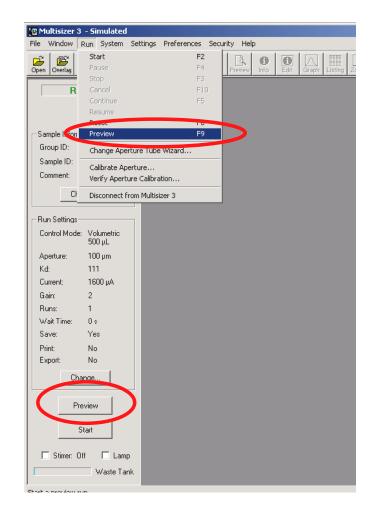

3. On the concentration index bar, verify that the concentration indicates OK; try to be as close as possible to 10%. If the concentration is high it can lead to an incorrect calibration. If the concentration is too low, the calibration will take a longer time to be completed. Adjust the concentration if it is necessary. When done,  $\textcircled{} \rightarrow Cancel$ .

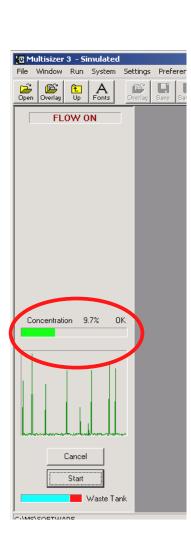

## 5.3 Running the Calibrator

U Multisizer	3 - Simulated											-	
File Window	Run System Sett	ings Preferences	Security Help										
Open Overlay	Start	F2		a a [	A Listing		Color Cursor Le		nfo Start		→!     🗖		
Open Overlay	Pause	F4	Preview	0 0 Info	iraph Listing 2	R X Zoom Stats	Color Cursor Le	gend	nfo Start	Stop Cancel	Junblock Emp	Calib.	Verify
	Stop Cancel	F3 F1											
R	Continue	F1 F5										$\sim$	
	Resume												
	Reset	F8											
Sample Infor	Preview	F9											
Group ID:	Change Aperture	Tube Wizard											
Sample ID.													
Comment	Calibrate Apertur Verify Aperture C												
	Disconnect from I	Multisizer 3											
- Run Settings													
-	e: Volumetric												
Control mod	500 µL												
Aperture:	100 µm												
Kd:	111												
Current:	1600 μA												
Gain:	2												
Runs:	1												
Wait Time:	0 s												
Save:	Yes												
Print	No												
Export:	No												
	hange												
	nange												
	Preview												
	TEVIEW												
[	Start												
🔲 Stirrer:	Off 🗖 Lamp												
	Waste Tank												
, Calibrata tha an	orturo Kel usina o colil	heation standard											

2. On the Calibrator Size box enter the modal value of the calibrator. This value appears on the assay sheet of the calibrator.

🔽 Multisizer 3 - Simulated	
	ettings Preferences Security Help
Open Overlay Up Fonts	Image: Save as     Image: Save a
READY	
	Calibrator Size
Sample Information Group ID: Sample ID: Comment: Change	Aperture:       100 μm       Serial Number:       0         Modal diameter of calificator:       μm         Recommended calibrator size:       10 μm         Concentration should be 5% to 10%         Cancel       Start
Control Mode: Volumetric 500 μL Aperture: 100 μm Kd: 111 Current: 1600 μA Gain: 2	

## 3. **∕∂→Start.**

_														
	😈 Multisizer 3	- Simulated												
	File Window R	Run System	Settings	Prefe	erences	Securi	ty Help	)						
		L A Up Fonts	Overlag	Save	Save as	Print	Preview	<b>B</b> Info	<b>B</b> Edit	Giraph	Listing	R	X Stats	
Ī	RE	ADY												
				Cali	brator !	5ize								
	Sample Informa Group ID: Sample ID: Comment: Cha	nge				·lodal d ommen	um liameter ded calib entration	of calibi prator si	ze: 10 μ	0.06 Im to 20 o 10%	μm μm Start			
	Control Mode: Aperture: Kd: Current:	Volumetric 500 μL 100 μm 111 1600 μA	I											

4. When the calibration run is completed, A→Accept New Kd. The instrument is calibrated

It is convenient as a future reference to clearly identify the calibration files. These files can be saved in the folder Calibration located in C:\MS.

File       Edit       View       Favorites       Tools       Help         Back       Forward       Up       Search       F         Address       MS       MS         Folders       X         Desktop       My Documents         My Documents       My Computer         My Computer       3% Floppy (A)         HP_PAVILION (C:)       P. RESTORE         ADDBEAPP       ATI         Cdrom       dosboot         HP       HP         HP_Calc       HP/Calc
Back Forward Up Search F Address MS Folders X Desktop My Documents My Computer PAVILION (C:) PADBEAPP Carom Carom Carom Carom PUSBDRV PAVILION PAVILION (C:) PADIBEAPP PAVILION C:) PADIBEAPP PAVILION C:) PADIBEAPP PAVILION PATI PADIBEAPP PAVILION PATI PADIBEAPP PATI PATI PADIBEAPP PATI PATI PADIBEAPP PATI PATI PADIBEAPP PATI PATI PADIBEAPP PATI PATI PATI PADIBEAPP PATI PATI PADIBEAPP PATI PATI PATI PADIBEAPP PATI PATI PATI PADIBEAPP PATI PATI PATI PATI PATI PATI PADIBEAPP PATI PA
Back Forward Up Search F Address MS Folders X Desktop My Documents My Computer PAVILION (C:) PAVILION (C:) PAVILION (C:) PAVILION (C:) PAUSBEAPP Carom Carom Carom PUSBDRV PHP AUDIO
Folders         ×           Image: Desktop         Image: Desktop
My Documents     My Computer     My Computer     All State Stress     My Computer     All State Stress     ADOBEAPP
Wy Computer     Wy Comput
3% Floppy (A:)     4%
HP_PAVILION (C:)     RESTORE     ADOBEAPP     ATI     cdrom     dosboot     EUSBDRV     HP     HP      AUDIO
Contraction Contr
dosboot December Dec
⊕- <u>C</u> HP ⊕- <u>C</u> HP AUDIO
HP AUDIO
- Comouse
Calibration
Demodeta
- 🗀 Languages
Samples
Preferences
SOM
HSOffice     My Documents
My Documents      My Photos

#### 5.4 Obtaining the mean Kd


For more consistent results we recommend the first time an aperture is calibrated to perform the following procedure:

- (1) Complete steps <u>5.1</u> through <u>5.3</u> ten times.
- (2) Record the Kd value for each calibration. Calculate the mean for the 10 values.
- (3) The mean for the Kd is only valid for the set aperture tube/electrolyte/instrument used for the calibration. In future calibrations the Kd should always be within  $\pm 4\%$  of the mean value from Step 5.4.2. The next table shows an example using a 10  $\mu$ m calibrator and a 100 $\mu$ m aperture tube.

Run	Kd
1	124.15
2	125.82
3	125.13
4	125.23
5	125.43
6	125.13
7	124.25
8	125.62
9	124.25
10	125.03
MEAN	125.00


#### 5.5 Entering the mean value for the Kd in the aperture list

1. On the Main Menu Bar,  $\mathcal{D} \rightarrow \mathbf{Run}$ , then  $\mathcal{D} \rightarrow \mathbf{Change Aperture Wizard}$ .

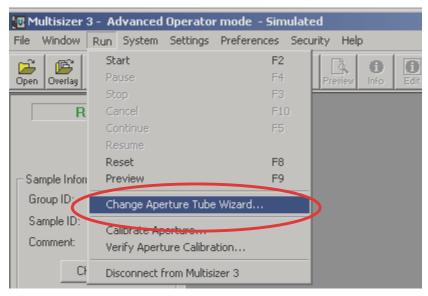


On the Change Aperture Wizard, <sup>⊕</sup>→Next to skip the steps up to Aperture Tube.
 <sup>⊕</sup>→Aperture Tube, in the new box, <sup>⊕</sup>→Edit List. On the Edit List Box, <sup>⊕</sup>→on the list the aperture size you just calibrated. Enter the serial number for the aperture tube, mean Kd value and <sup>⊕</sup>→Update.

Multisizer 3 - Advanced Op File Window Run System Se	erator mode - Simulated
🖻 🖻 🖪	Image: Save Save as       Print       Preview       Info       Edit       Image: Save Save as       Info       State       Image: Save Save as       Image: Save as
READY	Change Aperture Tule Tube Change Aperture Tube Diameter: 100 µm Kd: 111 OK
Sample Information Group ID: Sample ID: Comment:	C Remove the currently installed aperture tube C Install a new aperture tube C Place a beaker of clean electrolyte on the platform Diameter Number Kd Comment 20 0 30 Add New Tube
Change Run Settings Control Mode: Volumetric 500 µL	C Close the door       2       20       0       30       ▲         C Close the door       2       30       0       41       ▲         Select the electrolyte       Electrolyte       Electrolyte       Electrolyte       Electrolyte       Edit List         ISOTON II       ISOTON II       Electrolyte       Electrolyte       Electrolyte       ■         ISOTON II       Electrolyte       Electrolyte       Electrolyte       ■       ■         ISOTON II       Electrolyte       Electrolyte       ■       ■       ■         IO       0       100.856       ■       ■
Aperture: 100 µm Kd: 111 Current: 1600 µA Gain: 2 Runs: 1 Wait Time: 0 s Save: Yes Print: No Export: No Change Preview	<ul> <li>C Fill the system</li> <li>Adjust the metering pump</li> <li>Metering Pump</li> <li>Set current and gain</li> <li>Current &amp; Gain</li> <li>C Measure noise level</li> <li>Measure Noise</li> <li>C Calibration</li> <li>Verify</li> <li>C Calibrate</li> <li>C All done</li> <li>D one</li> <li>Close</li> <li>C Comment</li> <li>C C Comment</li> <li>C C Comment</li> <li>C C C C Comment</li> <li>C C C C C C C C C C C C C C C C C C C</li></ul>
Stirrer: Off Lamp	

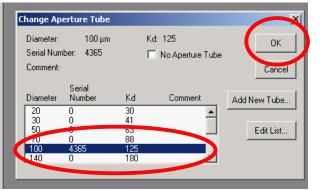

3. ∕**⊕→OK** 

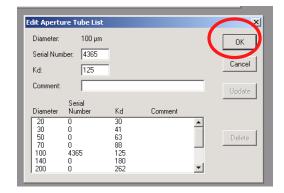

Change Aperture Tube Wizard	×		
< Back Next >			
C Remove the currently installed apertu	re tube	5.	<b>∕ी</b> →Close
C Install a new aperture tube			
C Place a beaker of clean electrolyte or	n the platform		
C Close the door			
C Select the electrolyte ISOTON II	Electrolyte		
C Select the new aperture tube	Aperture Tube		
<ul> <li>Fill the system</li> </ul>	Fill System		
C Adjust the metering pump	Metering Pump		
C Set current and gain	Current & Gain		
C Measure noise level	Measure Noise		
C Calibration Verify	Calibrate		
C All done	Done		
Close			


#### 5.6 Verifying the calibration


As the Multisizer 3 uses constant current, the instrument holds the calibration practically for indefinite time provided that the same aperture tube/electrolyte combination is used and the working conditions are stable. The calibration can be verified using the following procedure:

1. Place an ACCUVETTE® or a beaker\* with the electrolyte and a few drops of the calibrator into the analyzer unit.


\*An ACCUVETTE<sup>®</sup> (20 mL) is appropriate for apertures up to 100 $\mu$ m, for larger apertures a larger beaker should be used.

 On the Status Panel <sup>A</sup>→Preview, or in the Main Menu, <sup>A</sup>→Run,<sup>A</sup>→Preview.

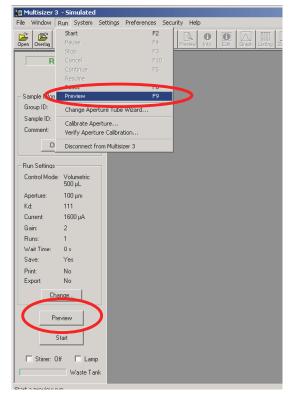

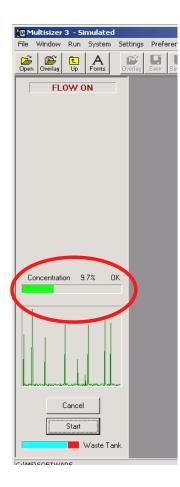

3. On the concentration index bar, verify that the concentration indicates OK. Try to be as close as possible to 10%. If the concentration is high it can lead to an incorrect result. If the concentration is too low, the verification will take a longer time to be completed. Adjust the concentration if it is necessary. When done,  $\neg \oplus \rightarrow Cancel$ .

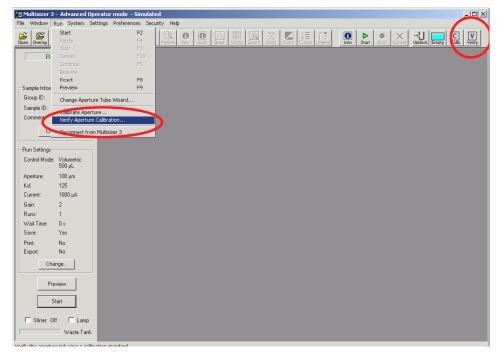

6.

5. On the Calibrator Size box enter the modal value of the calibrator. This value appears on the assay sheet of the calibrator. To choose the appropriate calibrator, refer to <u>Table 5.1.1</u>

Calibrator Size	×
Aperture: 100 μm	Serial Manuber: 4365
Modal diameter o	f calibrator:
Recommended calib	rator size: 10 μm το zo μm
Concentration s	hould be 5% to 10%
	Cancel Start
∕ð-→Start.	Calibrator Size
	Aperture: 100 μm Serial Number: 4365
	Modal diameter of calibrator: 10.02 μm
	Recommended calibrator size: 10 µm to 20 µm
	Concentration should be 5% to 10%
	Cancel Start

When the run is completed, the dialog box will indicate the status of the calibration. If it indicates verification is OK, <sup>A</sup>→Close. If the instrument is out of calibration, go to Step <u>5.2</u> and <u>5.3</u>.

It is convenient as a future reference to clearly identify the verification files. These files can be saved in the folder Verification located in C:\MS.

S MS						
<u> </u>	⊻iew	F <u>a</u> vo	rites	Tool	s <u>H</u> elp	)
. ⇔	$\rightarrow$	_	t.		Q	
Back		rd	Up	5	Search	F
Address 🗋	MS					
Folders						×
🚮 Desktop						
🗄 🗠 My Doc	uments					Π
🗄 🖳 My Con	nputer					
🕀 🚽 3½ F	loppy (/	A:)				
B- HP_						
	RESTO					
	DOBE4	\PP				
📄 🗎 🖻 🗖 A						
- <u></u>						
	losboot PUSBE					
		JHV				
		NO				
	IPCalc	10				
		à				
🗄 🧰 k	mouse					
🗄 🧰 К	PCMS					
- 🗀 n	nouse					
🖨 🔂 N	1S					-
	Calib					
	Demo					
	🛾 Langi					
	] Samp	oles				
ė- <b>G</b>	SOP	,				
		eferen	ces			
6						
	Verific					
📙 🗄 🔂 N	ly Photo	os				-

## **SECTION 6 SAMPLE ANALYSIS**


#### 6.1 Introduction

At the start of each analysis, it is good practice to check each item in the following list even though no change may be required from previous instrument settings.

Before running an analysis, the instrument must be in the following condition:

- (4) The system has been filled with electrolyte.
- (5) The correct aperture has been fitted.
- (6) A beaker with clean electrolyte has been placed on the platform.
- (7) The aperture is immersed into the electrolyte contained in the beaker
- (8) The electrolyte container is sufficiently filled (enough to raise the level sensor).
- (9) The Waste container is empty (enough to lower the level sensor).
- (10) The Multisizer 3 and the computer have been powered on with the control program running.
- (11) The appropriate analysis settings have been selected.

These settings are the essential parameters that must be established before the analysis to ensure that the particle pulses measured by the Multisizer 3 are optimally positioned within the digital converter range and have the best signal to noise ratio. Other parameters that in previous instrument designs have had to be preset, such as the number of channels of data, can be changed after, or even during an analysis. The **Analysis Settings** menu can be found on the main menu bar.

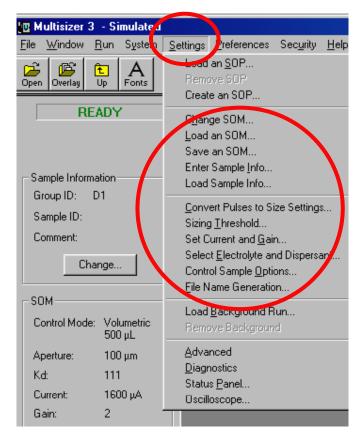

## 6.2 <u>Changing Standard Operating</u> <u>Method (SOM).</u>

	L Up Fonts	Load an <u>S</u> OP Remove SOP Create an SOP
Sample Informa		Change SOM Load on COM Save an SOM Enter Sample Info Load Sample Info
Sample ID: Comment: Chai	nge	<u>C</u> onvert Pulses to Size Settings Sizing <u>I</u> hreshold Set Current and <u>G</u> ain Select <u>E</u> lectrolyte and Dispersant Control Sample <u>O</u> ptions <u>F</u> ile Name Generation
SOM Control Mode:	Volumetric 500 μL	Load <u>B</u> ackground Run Remove Background
Aperture: Kd: Current:	100 μm 111 1600 μA	Advanced  Diagnostics Status Panel Uscilloscope
Gain: Runs:	2 · 1	
Wait Time:	0 s	
Save:	Yes	
Print:	No	

File Window Run System Settings Preferences Security Help

🖪 Multisizer 3 - Simulated

Standard Operating Me	thod (SOM)	×
Control Mode	i0 💌 econds	ОК
C Volumetric:	i00 🔻 µL	Cancel
🔿 Total Count: 🔤	0000 particles	<u>S</u> ave
🔿 Modal Count: 🔤 🗍	000 perícles	
C Manual		
Number of Flame: 2	Wait Between Rur	
Save Average of A	All <u>R</u> uns 🔽 Print A	verage Report
Aperture: 100 μm	Kd: 111	
Sizing Threshold: 1.63 Noise Level: 1.63		<u>T</u> hreshold
Current: 1600 µA	Gain: 2	Current & <u>G</u> ain
ISOTO	NII	Electrolyte
After Each Run		
Flush Aperture Tube		
🔽 Sa <u>v</u> e File 🔽	Include Pulse Data	
<(	G10>_ <s10>_<u2>.<x:< td=""><td>File Name</td></x:<></u2></s10>	File Name
✓ Export Data	CVS	6 Export
Print Report	Acrobat Distille	r Report
	(not specified)	Directory

## 6.2.1 Selecting the Control Mode

Select the appropriate control radio button on the Control sub-panel (refer to <u>Table 6.2-1</u>). Enter the variable associated with the method in the field.

• **Time.** Any time up to 99999 seconds can be entered. The analysis will stop when the selected time has elapsed.

• **Volumetric.** Any volume up to  $2000\mu$ L can be entered. The analysis will stop when the selected volume has been analyzed. Use this mode only with apertures not larger than  $280\mu$ m. (See Section 6.8).

• **Total Count.** The number for total count can be selected up to 525,000. The analysis will stop when the number corrected for coincidence corresponding to the selected raw count number has been accumulated.

• **Modal Count.** This is the number of particles counted in the modal channel. The modal count number can be selected up to 525,000 counts.

• Manual. The analysis has to be manually stopped.

Analysis for:	Control Mode				
	Manual Modal Count Total Count Time Volumetri				
Size Distribution	$\checkmark$		$\checkmark$		$\checkmark$
Particle Concentration	No	No	No	•	

♦ Refer to <u>Section 6.8</u>

Table 6.2.1 Selecting the control mode for the analysis

#### 6.2.2 Selecting the number of runs

You can select to run single or multiple runs. Selecting multiple runs allows you to insert a delay time between runs. When multiple runs are selected, the average of the runs can be automatically saved and printed by checking the appropriate boxes.

Standard Operating Method (SOM)	×
Control Mode	ок
C Volumetric: 500 y µL Ca	ancel
C Total Count: 10000 particles	ave
C Modal Count: 1000 particles	
C Manual	
Mumber of Runs: 3 Wait Between Runs: 60 s ▼ Save Average of All <u>R</u> uns ▼ Print Average Rep	econds ort
Aperture: 100 µm Kd: 111	
Sizing Threshold: 1.63 µm Noise Level: 1.63 µm	i
Current: 1600 µA Gain: 2 Current & G	ain
ISOTON II <u>E</u> lectrolyte	e
After Each Run	
🔽 <u>F</u> lush Aperture Tube	
☑ Save File ☑ Include Pulse Data	
<g10>_<s10>_<u2>.&lt;&gt;&gt; File Nan</u2></s10></g10>	ne
Print Report Acrobat Distiller Report	t
(not specified)	ry

Standard Operating Method (SOM)	×
Control Mode Time: 60 seconds	ОК
C Volumetric: 500 💌 pL	Cancel
C Total Count: 10000 particles	Save
C Modal Count: 1000 particles	
C Manual	
Number of Runs: 3 Wait Between Runs:	
Save Average of All <u>R</u> uns 🔽 Print Av	erage Report
Aperture: 100 μm Kd: 111	
Sizing Threshold: 1.63 μm Noise Level: 1.63 μm	Ihreshold
Current: 1600 µA Gain: 2 C	Current & <u>G</u> ain
ISOTON II	Electrolyte
After Each Run	
✓ Elush Aperture Tube	
Save File 🔽 Include Pulse Data	
<g10>_<s10>_<u2>.&lt;&gt;&gt;</u2></s10></g10>	File Name
☑ Export Data CVS	Export
Print Report Acrobat Distiller	Report
(not specified)	Directory

#### 6.2.3 Selecting the threshold

- The threshold is the size above which particles are counted. It is also the lower limit of the size distribution. This setting CAN NOT be changed during or after the analysis.

### 2. ∕Ĉ→Measuring Noise Level.

The instrument determines the lowest optimum threshold based on the measured noise level. Any number, however, can be manually entered for the threshold. If a number smaller than the size selected by the instrument is chosen, probability of noise interference will increase.

Standard Operating	Method (SOM)			×	
Control Mode			OK		
Time:	60 💌 se	Sizing Th	reshold		×
C Volumetric:	500 💌 μί	Apertu	re: 100 μm		
C Total Count:	10000 pa	<u>T</u> hresho	ld: 1.63	μm	Cancel
C Modal Count:	1000 pa	_			
C Manual		Noise Lev	el: 1.63 μm 1.6% of aper	ture diame	<u>D</u> efaults
Number of Runs: 3	 	6	Measure <u>N</u> oise I	Level	
🔽 Save Average	of All <u>R</u> uns	(Use clea	an electrolyte wri	en measu	ring noise level)
Aperture: 100 μm	Kd: 111				
Sizing Threshold: Noise Level:		Γ	<u>T</u> hreshold		
Current: 1600 μA	. Gain: 2	(	Current & <u>G</u> ain		
ISO	ITON II	Γ	Electrolyte		
After Each Run					
Elush Aperture T	ube				
🔽 Sa <u>v</u> e File	🔽 Include Pulse	Data			
	<g10>_<s10>_</s10></g10>	<u2>.<x></x></u2>	File Name		
🔽 Export Data		CVS	Export		
Print Report	Acro	bat Distiller	Report		
	(not	specified)	Directory		

## 6.2.4 Adjusting current and gain. 1. ∽ີ → Current & Gain

Standard Operating Method (SOM)
Control Mode OK OK
C Volumetric: 500 y pt. Cancel
C Total Count: 10000 particles Save
C Modal Count: 1000 particles
C Manual
Number of Runs: 3 Wait Between Runs: 60 seconds
Save Average of All Runs  Print Average Report
Aperture: 100 μm Kd: 111
Sizing Threshold: 1.63 μm Noise Level: 1.63 μm
Current: 1600 μA Gain: 2 Current & <u>G</u> ain
ISOTON II <u>Electrolyte</u>
After Each Run
✓ Elush Aperture Tube
🔽 Save File 🔽 Include Pulse Data
<g10>_<s10>_<u2>.<x> File Name</x></u2></s10></g10>
Export Data CVS Export
Print Report Acrobat Distiller Report
(not specified) Directory

Standard Operating Method (SOM) ⊂ Control Mode 2. ℃→Auto	Sat
	Jei
C Time: 10 records	
€ Volumetric: 500 ▼ μL Cancel	
C Total Count: 10000 particles Save	
C Modal County Innon Settings	
C Manual	
Aperture Current 1600 pA	
Number of Runs: Preamp Gain-	
C 1	
Aperture: 100 © 2	
Auto-Set	
Sizing Threshology 8	
Current: 160 C 16	
C 32	
After Each Run	
Elush Aperture Tube	
🔽 Sa <u>v</u> e File 🔽 Include Pulse Data	
<g10>_<s10>_<r2>.&lt;&lt;&gt;&gt; File Name</r2></s10></g10>	
Export Data	
Erint Report HP DeskJet 840C Series Report	

3. Select Particles or Cells. ∕∂→OK

If cells are to be analyzed, it is very important to select cells in this step. Not selecting cells could cause them to be damaged due to a higher current setting.

Every time the current and gain are adjusted, it is necessary to measure the noise level.

## 6.2.5 Changing the Electrolyte

If you need to use a different electrolyte with the same aperture tube, perform the following procedure:

Automatic Current and Gain

Current

1600 µA

800 pA

Cancel

Maximum Cell Volume: 0

38.4 µL/s

14.6 k ohms

Gain

2

4

Flow Rate:

Resistance:

Particles:

C <u>C</u>ells:

ΟK

X

	Standard Operating Method (SOM)
_	Control Mode OK
1. ∽ <del>0</del> →Electrolyte	Image: Time:     60     ▼     seconds
	C Total Count: 10000 particles Save
	C Modal Count: 1000 particles
	C Manual
	Number of Runs: 3 Wait Between Runs: 60 seconds
	Save Average of All Buns
	Aperture: 100 µm Kd: 111
	Sizing Threshold: 1.63 μm Noise Level: 1.63 μm
	Current: 1600 µA Gain: 2 Current & Gain
	ISOTON II <u>E</u> lectrolyte
	- After Each Run
	✓ Elush Aperture Tube
	I Save File I Include Pulse Data
	<g10>_<s10>_<u2>.<x></x></u2></s10></g10>
	Print Report Acrobat Distiller Report
	(not specified)
Standard Operating Method (SOM)	×
Control Mode	<b>2.</b> Select the electrolyte by
C Time: 10 💌 seconds	double clicking in the drop
O Volumetric: 500   μL	Cancel down list or typing its name
C Total Count: 10000 particles	in the box.
Select Electrolute and Dispersant	
C Ma	X
C Ma Electrolyte: Fout ON II	
C Ma Electrolyte: SUTON II Dispersant: ISOTON II - Glycerol 6:4	OK Cancel
C Ma Electrolyte: SUTON II SOTON II ISOTON II - Glycerol 6:4 0.9% Sodium Chloride in Water 4% Sodium Chloride in Water	T Cancel
C Mo C Ma Electrolyte: Solution II ISOTON II - Glycerol 6:4 0.9% Sodium Chloride in Water 4% Sodium Chloride in Water 2% Anmonium Thiocyanate in 5% Lithium Chloride in Methyl A	r Isopropyl Alcohol
C Ma Electrolyte: SUTON II Discersant: ISOTON II - Glycerol 6:4 0.9% Sodium Chloride in Water 2% Ammonium Thiocyanate in S% Lithium Chloride in Action 5% Lithium Chloride in Acetone 5% Lithium Chloride in Acetone	r Isopropyl Alcohol Alcohol
C Ma Electrolyte: Solution II Dispersant: Aperture: 100 µm Sizing Threshold 1.6 Sizing Threshold 1.6 Sizing Threshold 1.6 Sizing Threshold 1.6 C Ma Sizing Threshold 1.6 Sizing Threshold 1	r Isopropyl Alcohol Alcohol e Ethanol
C Ma Electrolyte: Solution II Dispersant: Aperture: 100 µm Sizing Threshold 1.6 C Ma Electrolyte: Solution II - Glycerol 6.4 0.9% Sodium Chloride in Water 2% Ammonium Thiocyanate in 5% Lithium Chloride in Acetone 5% Lithium Chloride in Acetone 5% Lithium Chloride in Acetone 4% Lithium Iodide in Acetone 4% Monomium Thiocyanate in 4% Remonium Thiocyanate in 4% Remonium Thiocyanate in 100 µm	r Isopropyl Alcohol Alcohol
C       Ma       Electrolyte:       SutTON II         Number       Disrefsant:       ISOTON II - Glycerol 6:4         Number       Disrefsant:       0.9% Sodium Chloride in Water         Aperture:       1L0 μm       Sizing Thresholo 1:6         Sizing Thresholo 1:6       4% Zinc Chloride in Acetone         Sizing Thresholo 1:6       4% Lithium Chloride in Acetone         Gurrent:       160 μA	r Isopropyl Alcohol Alcohol e Ethanol
C       Ma       Electrolyte:       SutTON II         Number       Disrefsant:       ISOTON II - Glycerol 6:4         Number       Disrefsant:       0.9% Sodium Chloride in Water         Aperture:       1L0 μm       Sizing Thresholo 1:6         Sizing Thresholo 1:6       4% Zinc Chloride in Acetone         Sizing Thresholo 1:6       4% Lithium Chloride in Acetone         Gurrent:       160 μA	r Isopropyl Alcohol Alcohol Ethanol It & <u>G</u> ain
C       Ma       Electrolyte:       SutTON II         Number       Disrefsant:       ISOTON II - Glycerol 6:4         Number       Disrefsant:       0.9% Sodium Chloride in Water         Aperture:       1L0 μm       Sizing Thresholo 1:6         Sizing Thresholo 1:6       4% Zinc Chloride in Acetone         Sizing Thresholo 1:6       4% Lithium Chloride in Acetone         Gurrent:       160 μA	r Isopropyl Alcohol Alcohol B Ethanol It & Gain Strolyte
C       Ma       Electrolyte:       Satition II         Number       Dispersant       ISOTON II       ISOTON II         Number       Dispersant       ISOTON II       O.9% Sodium Chloride in Water         Aperture:       110 µm       S% Lithium Chloride in Methyl A         Aperture:       110 µm       S% Lithium Chloride in Acetone         Sizing Thresholo       1.6       4% Lithium Chloride in Acetone         4% Ammonium Thiocyanate in       Current:       1600 µA         ISOTON II       Elector	r Isopropyl Alcohol Alcohol Ethanol It & <u>G</u> ain
C       Ma       Electrolyte:       Faulton II         Number       Dispersant       ISOTON II - Glycerol 6:4         Number       Dispersant       ISOTON II - Glycerol 6:4         Aperture:       140 µm       Sodium Chloride in Water         Aperture:       140 µm       Sizing Thresholo 1:6         Sizing Thresholo 1:6       4% Zinc Chloride in Acetone         Sizing Thresholo 1:6       4% Lithium Chloride in Acetone         Uurrent:       1600 µA         SizoTON II       Elector	r Isopropyl Alcohol Alcohol B Ethanol It & Gain Strolyte Standard Operating Method (SOM) Control Mode Time: Desconds OK Control Mode
<ul> <li>Ma</li> <li>Electrolyte: Solution II</li> <li>Isotron II</li> <li>Sotian Chloride in Water</li> <li>Aperture: 110 μm</li> <li>Sizing Threshols 1.6</li> <li>Lithium Chloride in Acetone</li> <li>Lithium Chloride in Acetone</li> <li>Ammonium Thiocyanate in</li> <li>Current: 1600 μA</li> <li>Gain: 2</li> <li>Current</li> <li>Isotron II</li> <li>Electron II</li> </ul>	r Isopropyl Alcohol Alcohol stebhanol ht & <u>Gain</u> strolyte Standard Operating Method (SOM) Control Mode Control Mod
<ul> <li>Ma</li> <li>Electrolyte: Button:</li> <li>Soton:</li> <li>So</li></ul>	Image: Conceleration of the second seco
<ul> <li>Ma</li> <li>Electrolyte: Button:</li> <li>Sotton:</li> <li>Sotton:&lt;</li></ul>	Isopropyl Alcohol         Alcohol         Settext         Ethanol         tt & Bain         strolyte         Standard Operating Method (SOM)         Control Mode         OK         Control Mode         Time:         10 y         seconds         OK         Cancel         Total Count:         10000 perticles         Mo         Select Electrolyte and Dispersant
<ul> <li>Ma</li> <li>Electrolyte: Solution II</li> <li>ISOTON II</li> <li>O3% Sodium Chloride in Water</li> <li>Aperture: 110 μm</li> <li>Sizing Threshold 1.6</li> <li>Sizing Threshold 1.6<!--</th--><th>GK         r         Isopropyl Alcohol         Alcohol         Standard Operating Method (SOM)         Control Mode         OK         Cancel         Mo         Select Electrolyte and Dispersant         Mo         Electrolyte:         ISCINII</th></li></ul>	GK         r         Isopropyl Alcohol         Alcohol         Standard Operating Method (SOM)         Control Mode         OK         Cancel         Mo         Select Electrolyte and Dispersant         Mo         Electrolyte:         ISCINII
<ul> <li>Ma</li> <li>Electrolyte: Button:</li> <li>Sotton:</li> <li>Sotton:&lt;</li></ul>	Isopropyl Alcohol         Alcohol         Standard Operating Method (SOM)         Ethanol         Standard Operating Method (SOM)         Control Mode         OK         Control Mode         Time:         10         seconds         Cancel         Mo         Select Electrolyte and Dispersant         Ma         Electrolyte:         ISE-control
<ul> <li>Ma</li> <li>Electrolyte: Solution II</li> <li>ISOTON II</li> <li>O3% Sodium Chloride in Water</li> <li>Aperture: 110 μm</li> <li>Sizing Threshold 1.6</li> <li>Sizing Threshold 1.6<!--</th--><th>r Isopropyl Alcohol Alcohol Ethanol It &amp; Gain strolyte Stordard Operating Method (SOM) Control Mode Time: Sol y L Control Mode Time: Sol y L Cancel Total Count: Total Count: To</th></li></ul>	r Isopropyl Alcohol Alcohol Ethanol It & Gain strolyte Stordard Operating Method (SOM) Control Mode Time: Sol y L Control Mode Time: Sol y L Cancel Total Count: Total Count: To
<ul> <li>Ma</li> <li>Electrolyte: Solution II</li> <li>ISOTON II</li> <li>O3% Sodium Chloride in Water</li> <li>Aperture: 110 μm</li> <li>Sizing Threshold 1.6</li> <li>Sizing Threshold 1.6<!--</th--><th>r Isopropyl Alcohol Alcohol Ethanol It &amp; Gain strolyte Stordard Operating Method (SOM) Control Mode Control Mode</th></li></ul>	r Isopropyl Alcohol Alcohol Ethanol It & Gain strolyte Stordard Operating Method (SOM) Control Mode Control Mode
<ul> <li>Ma</li> <li>Electrolyte: Solution II</li> <li>ISOTON II</li> <li>O3% Sodium Chloride in Water</li> <li>Aperture: 110 μm</li> <li>Sizing Threshold 1.6</li> <li>Sizing Threshold 1.6<!--</th--><th>Image: Concelent of the second se</th></li></ul>	Image: Concelent of the second se
<ul> <li>Ma</li> <li>Electrolyte: Solution II</li> <li>ISOTON II</li> <li>O3% Sodium Chloride in Water</li> <li>Aperture: 110 μm</li> <li>Sizing Threshold 1.6</li> <li>Sizing Threshold 1.6<!--</th--><th>Image: Cancel         Isopropul Alcohol         Isopropul Alcohol         Isopropul Alcohol         Image: Cancel         Ethanol         It &amp; Bain         stolyte         Standard Operating Method (SOM)         Image: Control Mode         &lt;</th></li></ul>	Image: Cancel         Isopropul Alcohol         Isopropul Alcohol         Isopropul Alcohol         Image: Cancel         Ethanol         It & Bain         stolyte         Standard Operating Method (SOM)         Image: Control Mode         <
<ul> <li>Ma</li> <li>Electrolyte: Solution II</li> <li>ISOTON II</li> <li>O3% Sodium Chloride in Water</li> <li>Aperture: 110 μm</li> <li>Sizing Threshold 1.6</li> <li>Sizing Threshold 1.6<!--</th--><th>Image: Cancel         Isopropyl Alcohol         Alcohol         Standard Operating Method (SOM)         It &amp; Gain         Strolyte         Standard Operating Method (SOM)         Image: Control Mode         Image: Contret     </th></li></ul>	Image: Cancel         Isopropyl Alcohol         Alcohol         Standard Operating Method (SOM)         It & Gain         Strolyte         Standard Operating Method (SOM)         Image: Control Mode         Image: Contret
<ul> <li>Ma</li> <li>Electrolyte: Solution II</li> <li>ISOTON II</li> <li>O3% Sodium Chloride in Water</li> <li>Aperture: 110 μm</li> <li>Sizing Threshold 1.6</li> <li>Sizing Threshold 1.6<!--</th--><th>Image: Cancel Cancel</th></li></ul>	Image: Cancel

Every time the electrolyte is changed, it is necessary to adjust current and gain and measure the noise level. If this is the first time using the new electrolyte, it is necessary to calibrate the aperture. If the electrolyte has been used before, it is necessary to verify the calibration (See Section 5)..

#### 6.2.6 Flushing the Aperture Tube

By selecting this option, the instrument will automatically flush the aperture every time a run is completed. To activate auto flush, check the appropriate box.

Standard Operating	Method (SOM)	×
© Time:	60 💌 seconds	ОК
C Volumetric:	500 💌 μL	Cancel
C Total Count:	10000 particles	Save
C Modal Count:	1000 particles	
C Manual		
Number of Runs: 3		uns: 60 seconds
🔽 Save Average	of All <u>R</u> uns 🔽 Print	Average Report
Aperture: 100 μm	Kd: 111	
Sizing Threshold: Noise Level:		<u>T</u> hreshold
Current: 1600 μ4	Gain: 2	Current & <u>G</u> ain
ISC	)TON II	Electrolyte
After Each Run		
Elush Aperture T	ube	
Save File	🔽 Include Pulse Data	
	<g10>_<s10>_<u2>.&lt;&gt;</u2></s10></g10>	K> File Name
Export Data	C	/S Export
Print Report	Acrobat Distil	ler Report
	(not specifie	d) <u>D</u> irectory

Standard Operating Method (SOM)	1
© Time: 60 ▼ seconds	
C Volumetric: 500 y µL Cancel	
C Total Count: 10000 particles Save	
C Modal Count: 1000 particles	
C Manual	
Number of Runs: 3 Wait Between Runs: 60 seconds	
Save Average of All <u>B</u> uns  Frint Average Report	
Aperture: 100 μm Kd: 111	
Sizing Threshold: 1.63 µm Noise Level: 1.63 µm	
Current: 1600 µA Gain: 2 Current & <u>G</u> ain	
ISOTON II <u>E</u> lectrolyte	
After Each Run	
Elush Aperture Tube	
🔽 Save File 🛛 🔽 Include Pulse Data	
(C10;_:(616)_:(U2).(X) File Name	
Export Data CVS Export	
Print Report Acrobat Distiller Report	
(not specified) Directory	

#### 6.2.7 Saving the Data.

To automatically save the results after each analysis, check the **Save File** box. By checking this box, the run files are saved without the pulse data. To save all the original information from the pulses, check the **Save Pulse** Data box. Saving the pulse data allows you to change resolution, range and other parameters of the analysis without needing to re-run the sample.

The size of the files containing the pulse data is significantly large. Depending on the use of the instrument, it is advisable to transfer the files regularly to an appropriate storage media. A CD writer or similar storage device can be used.

#### 6.2.8 Generating the File Name

To identify the files from different analyses, a file name should be generated. The operator can customize it by checking the boxes of the information to be part of the file name. On the **Standard Operating Method** dialog box,  $\bigcirc \rightarrow$  **File Name**. In the File Generation Name Dialog box, customize name,  $\bigcirc \rightarrow$  **Use Example** and  $\bigcirc \rightarrow$  **OK**.

Standard Operating Method (SOM)	
Control Mode	
C Volumetric: 500 💌 pL Cancel	File Name Generation
C Total Count: 10000 particles Save	Image: Complex comple
C Modal Count: 1000 particles	Cancel
C Manual	Clear All
Number of Runs: 3 Wait Between Runs: 60 seconds	Up to 0 characters Up to 0 characters Defaults
Save Average of All Buns Frint Average Report	□ <d> <u>D</u>ate □ <t#> <u>I</u>ime □</t#></d>
Aperture: 100 µm Kd: 111	( 3 Jul 2001) 2: 10-33 💌
Sizing Threshold: 1.63 µm Noise Level: 1.63 µmhreshold	✓ <r#> Bun Number     ✓     Unique Number       2     digits     0     digits</r#>
Current: 1600 µA Gain: 2 Current & <u>G</u> ain	
ISOTON II <u>E</u> lectrolyte	Multisizer 3: #M3 Multisizer II: #M2 Z2: ZZ2
After Each Run	
Elush Aperture Tube	Example: <u10>_<s10>_<b2>.&lt;&gt;&gt;</b2></s10></u10>
🔽 Save File 🔽 Include Pulse Data	Use E <u>x</u> ample
<g10>_<s10>_<u2(<>&gt; File Name</u2(<></s10></g10>	Template: <610>_<\$10>_ <r2>.&lt;&gt;&gt;</r2>
Export Data CVS Export	File names will be 8 to 28 characters long
Print Report Acrobat Distiller Report	
(not specified) Directory	

#### 6.2.9 Exporting Data

To automatically export the analysis data after each run is completed, check the Export Data box. To select the export data format and destination,  $\neg \ominus \rightarrow Export$ , select the options and  $\neg \ominus \rightarrow OK$ .

Standard Operating Method (SOM)	
Control Mode	
© Time: 60 ▼ seconds	
C Volumetric: 500 y µL Cancel	Data Export
C Total Count: 10000 particles Save	
C Modal Count: 1000 particles	Export Extension: CVS OK
C Manual	I Sample Info I Short Cancel
Number of Runs: 3 Wait Between Runs: 60 seconds	I Statistics □ All Clear All
Save Average of All Buns     Print Average Report	Averaged Statistics
	Interpolation I
Aperture: 100 μm Kd: 111	Size Listing C 123456.78
Sizing Threshold: 1.63 μm Noise Level: 1.63 μm Ihreshold	Fulse Listing
Current: 1600 µA Gain: 2 Current & <u>G</u> ain	Data Format
ISOTON II <u>E</u> lectrolyte	Comma delimited Border width: 1
	C Tab delimited Cell spacing: U
After Each Run	C Blank delimited
Elush Aperture Tube	
I Save File I Include Pulse Data	Export Directory
<g10>_<s10>_<u2>.<x> File Name</x></u2></s10></g10>	C Current directory C Curtom directory Browse
Export Data	Custom directory Browse
Princheport Acrobat Distiller Heport	
(not specified) Directory	

#### 6.2.10 Printing the Analysis Report

To automatically obtain a printed report of the analysis after completing each run,  $- \partial \rightarrow \mathbf{Report}$ .

Select the content of the printed report and  $\neg \ominus \rightarrow OK$ .

Standard Operating Method (SOM)		
Control Mode C Time: 60 seconds OK		
C Volumetric: 500 y µL Cancel		
C Total Count: 10000 particles Save	Printed Report	×
C Modal Count: 1000 particles	Short	Clear <u>A</u> ll OK
C Manual	Size	Cancel
Number of Runs: 3 Wait Between Runs: 60 seconds	Graphs Select	- Size Statistics
Save Average of All <u>B</u> uns	Averaged Statistics	Mass Select
Aperture: 100 µm Kd: 111	Compare Statistics	□ Number
Sizing Threshold: 1.63 µm Noise Level: 1.63 µmhreshold	Listing Select	└ Volume 🔽 % 🗂 /mL
Current: 1600 µA Gain: 2 Current & Gain	Interpolation Select	🗖 Surface Area 🗖 % 🗖 /mL
ISOTON II <u>E</u> lectrolyte	P <u>u</u> lses	Irend
After Each Run	Graph Select	Graph Select
Elush Aperture Tube	Listing Select	Listing Select
☞ Sa <u>v</u> e File ☞ Include Pulse Data	Printer	
<g10>_<s10>_<u2>.<x> File Name</x></u2></s10></g10>	Use Default Printer	Select
Export Data CVS Export	C Use Specified Printer:	
Print Report Acrobat Distiller Report		
(not specified)		

#### 6.2.11 Changing the Destination Directory.

Standard Operating Method (SOM)	
Control Mode OK OK	
C Volumetric: 500 PL Cancel	
C Total Count: 10000 particles Save	
C Modal Count: 1000 particles	
C Manual	
Number of Runs: 3 Wait Between Runs: 10 seconds	
Save Average of All Buns  Frint Average Report	Select the Run Directory
Aperture: 100 μm Kd: 111	C Use the current directory
Sizing Threshold: 1.63 μm Noise Level: 1.63 μm	© Use a gustom directory
Current: 1600 µA Gain: 2 Current & <u>G</u> ain	Basant
ISOTON II <u>E</u> lectrolyte	Directories: C:\MS\M3x
After Each Run	Browse
Elush Aperture Tube	
I Save File II Include Pulse Data	
<g10>_<s10>_<u2>.<x> File Name</x></u2></s10></g10>	
Export Data CVS Export	
Print Report Acrobat Distiller Report	
(not specifier Directory	


At this moment the Standard Operating Method has been selected. In the Standard Operating Method Panel, ⊸∂→OK.

Standard Operating Me	ethod (SOM)	×
Control Mode		ОК
Time:	50 💌 seconds	
C Volumetric:	500 🔽 μL	Cancel
🔿 Total Count: 🔤	10000 particles	Save
C Modal Count:	1000 particles	
C Manual		
<u>N</u> umber of Runs: 3 ✓ Save Average of <i>i</i>	<u>W</u> ait Between Ru All <u>R</u> uns <b>I</b> ✓ Print /	ns: 60 seconds Average Report
Aperture: 100 μm	Kd: 111	
Sizing Threshold: 1.6 Noise Level: 1.6		<u>T</u> hreshold
Current: 1600 µA	Gain: 2	Current & <u>G</u> ain
ISOTO	IN II	<u>E</u> lectrolyte
After Each Run		
Flush Aperture Tube	e	
🔽 Sa <u>v</u> e File 🔽	Include Pulse Data	
<	G10>_ <s10>_<u2>.<x< td=""><td>File Name</td></x<></u2></s10>	File Name
Export Data	CV	S Export
I Print Report	Acrobat Distille	er Report
	(not specified	<u>D</u> irectory

#### 6.3 <u>Convert Pulses to Size</u> <u>Settings</u>

Although these settings can be changed during or after the analysis, it is convenient to select these settings before the analysis. In this way, they can be saved with the Standard Operating Method.

1. <sup></sup> **⊕**L→Settings, **⊕**L→Convert Pulses to Size Settings.

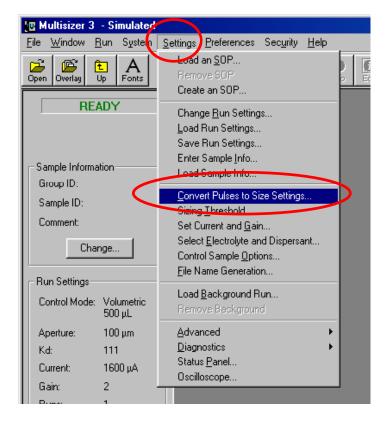

Convert Pulses to Size Settings	×
Sizing Threshold: 1.63 μm Aperture: 100 μm 1.6% of aperture diameter	OK
300 💌 Size Bins from 2 μm to 60 μm	Cancel
Bin Spacing C Linear Diameter	<u>2</u> % to 60%
<ul> <li>Log Diameter</li> <li>Linear Volume</li> </ul>	<u>D</u> efaults
C Log Volume	

2. Select the settings

• **Size Bins.** The number of size bins can be selected from 4 up to 300. Higher resolution is achieved when a larger number of channels is selected.

• **Bin Spacing.** Use Logarithmic when the X axis units are logarithmic, use Linear

when the X axis units are linear. Use Diameter to have the X-axis is in diameter units, use volume to have the X axis in volume units.

• From – To. Enter the desired range for the analysis. The range can be selected from the threshold size up to the size of the aperture.


The size distribution is guaranteed to be linear from 2 to 60% of the aperture diameter.

 Multisizer II Edit. This function is similar as in the Multisizer IIe. It is used to correct for non-axial pass of particles through the aperture. It can be used only with narrow size distributions.


#### 6.4 Saving Standard Operating Method

All the settings for the analysis can be saved for future use. The saved file containing the Standard Operating Method can also be used to create a **Standard Operation Procedure (SOP)** (see <u>Section 9</u>).

To save the Standard Operating Method  $\mathcal{D}$ -Settings on the Main Menu bar and then  $\mathcal{D}$ -Save SOM.

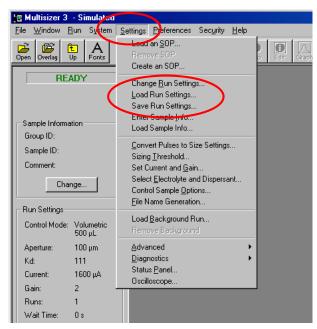

The **SOM** file can be saved in the **SOM** folder located in C:\MS\SOP. Any number of **SOM** files can be saved; they can be used at a later time for creating a **Standard Operating Procedure** (see <u>Section 9</u>).

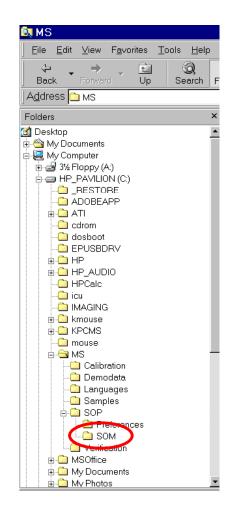

To use a Standard Operating Method previously saved,  $\mathcal{D}$ -Settings on the Main Menu bar, then  $\mathcal{D}$ -Load SOM. To change the Standard Operating Method,  $\mathcal{D}$ -Change SOM.

#### 6.5 Entering Sample Information

By entering the sample information, it is possible to save all the information regarding the sample in the same file for the analysis. When the determination of concentration is needed, (counts/ml or counts/g), it is necessary to enter additional information in order for the software to do the calculation.

1. ∽<sup>®</sup>→Change on the Sample Information sub-panel in the Control Panel,

or ↔ System on the Main Menu bar ↔ Enter Sample Info...

Multisizer 3	- Simulated		
<u>File Window F</u>	<u>R</u> un System	<u>S</u> ettings <mark>P</mark> eferences Sec <u>u</u> rity <u>H</u> elp	
	L A Jp Fonts	Lead an <u>S</u> OP Remove SOP Create an SOP	
Sample Informa Group ID:	ADY Ition	Change <u>B</u> un Settings Load Run Settings Save Run Settings Enter Sample Info Load Sample Info	
Sample ID: Comment Char Run Settings Control Mode:	Volumetric 500 μL	<u>Convert</u> Pulses to Size Settings Sizing <u>T</u> hreshold Set Current and <u>G</u> ain Select <u>E</u> lectrolyte and Dispersant Control Sample <u>D</u> ptions <u>F</u> ile Name Generation	
Aperture: Kd: Current: Gain:	100 μm 111 1600 μΑ 2	Load <u>B</u> ackground Run Remove Background	
Runs: Wait Time: Save: Print:	1 Os No No	Advanced ► Diagnostics ► Status Panel Oscilloscope	
Export: Chai	No nge		

2. In the dialog box, enter all the required information.

Enter Sample Info		X
Group ID:	Clear <u>A</u> ll	ОК
Sample ID:		Cancel
Operator:		
Bar Code:		
Comment:	Variable <u>2</u> :	0
Run number: 1	Shape <u>F</u> actor:	0
Sample	Electrolyte <u>V</u> olume:	0 mL
Vojume:	Analytic Volume:	0 μL
C Mass: 0 mL	Use Dilution Factor	0
De <u>n</u> sity: 0 g/mL	Control Sample Sample	Statistics
Template: <g10>_<s10>_<r2>.&lt;</r2></s10></g10>	* [	File Name
File Name:01.#M3		

Group ID, Sample ID, Operator, Bar Code and Comment are optional fields and their only purpose is to identify and classify the analyses. The run number will track the analyses sequence of each Group ID.

If it is necessary to determine the concentration of particles in the sample (counts), the followings fields must be filled in.

Sample volume or mass: This is the weight or volume of sample used for the analysis.

**Electrolyte volume:** This is the volume of electrolyte used for the analysis.

Analytical volume: This is the volume of sample analyzed.

Enter Sample Info	×
<u>G</u> roup ID:	Clear <u>A</u> ll OK
<u>S</u> ample ID:	Cancel
Operator:	
Bar Code:	Variable <u>1</u> : 0
	Variable <u>2</u> : 0
<u>C</u> omment:	
Run number: 1	Shape Eactor: 0
Sample	Electrolyte Volume: 0 mL
Volume:	Analytic Volume: Ο μ
C Mass: 0 mL	Use Dilution Factor,
De <u>n</u> sity: 0 g/mL	Control Sample Sample Statistics
Tamalaka (C10) (C10) (D2) (A)	File Name
Template: <g10>_<s10>_<r2>.<x> File Name: 01.#M3</x></r2></s10></g10>	

Other fields in the Enter Sample Info dialogue box.

Variable 1 and Variable 2: These fields won't affect the results and are just to enter any variable you would like to save in the sample information.

**Shape Factor:** If a constant correction is needed, a factor can be entered and it will affect the results. For example, if a fibrous material is analyzed and instead of the equivalent spherical diameter, the length is required, a conversion factor can be entered to compensate for the difference in shape.

**File Name:** This button has the same function as described in <u>Section 6.2.8</u> **Control Sample:** By checking this box, you indicate that the sample to be analyzed is a control sample.

**Density:** This information needs to be entered if the mass size distribution is required.

Enter Sample Info	×
<u>G</u> roup ID:	Clear <u>A</u> ll OK
<u>S</u> ample ID:	Cancel
Operator:	Variable <u>1</u> : 0
<u>B</u> ar Code:	
Comment:	Variable <u>2</u> : 0
Run number: 1	Shape <u>F</u> actor: 0
Sample	Electrolyte Volume: 0 mL
Vojume:	Analytic Volume: 0 µL
C Mass: 0 mL	Use Dilution Factor:
De <u>n</u> sity: 0 g/m	Control Sample Sample Statistics
Template: <g10>_<s10>_<r2>.&lt;</r2></s10></g10>	× File Name
File Name:01.#M3	

Enter Sample Info	× • • • • • • • • • • • • • • • • • • •
<u>G</u> roup ID:	Clear <u>A</u> ll OK
<u>S</u> ample ID:	Cancel
<u>O</u> perator:	
Bar Code:	Variable <u>1</u> : 0
	Variable <u>2</u> : 0
<u>C</u> omment:	Shape <u>F</u> actor: 0
Run number: 1	
Sample	Electrolyte Volume: 0 mL
• Volume:	Analytic Volume: 0 μL
C Mass: 0 mL	Use Dilution Factor:
De <u>n</u> sity: 0 g/mL	Control Sample Sample Statistics
Template: <g10>_<s10>_<r2>.<x> File Name: 01 #M3</x></r2></s10></g10>	File Name

**Use Dilution Factor:** By checking this box, it is possible to use a factor if dilution has been used when preparing the sample. The dilution factor will affect the results.

**Sample Statistics:** When a sample is selected as "Control Sample", its statistical parameters and acceptable limits can be entered here.

To use the Sample Statistics as a reference, save this control sample under an easily identifiable name.

Enter Sample Info		×
<u>G</u> roup ID:	Clear <u>A</u> ll	OK
<u>S</u> ample ID:		Cancel
Operator:		
	Variable <u>1</u> : 0	
<u>B</u> ar Code:	Variable <u>2</u> : 0	
<u>C</u> omment:		_
Run number: 1	Shape <u>F</u> actor: 0	
	Electrolyte ⊻olume: 0	mL
Sample	Analytic Volume:	μL
⊙ Vojume: O mL	· · · · · · · · · · · · · · · · · · ·	με
	Use Dilution Factor:	
De <u>n</u> sity: 0 g/mL	Control Sample Sample S	tatistics
Template: <g10>_<s10>_<r2>.&lt;&gt;</r2></s10></g10>	⇔Fil	e Name
File Name:01.#M3		

	Edit San	nple Statistic	50		
		Nominal Value	Lower Limit	Upper Limi	ОК
	Mean:	0	0	0	Carcel
1	Median:	0	0	0	
	Mode:	0	0	0	
	S.D.:	0	0	0	
	C.V.:	0	0	0 %	


This information can be use as reference or specification for other samples. When done,  $\textcircled{} \rightarrow OK$ .

## 6.6 <u>Running a Background.</u>

The background is an analysis of the electrolyte without sample material; it is also called a "Blank Analysis". It is advisable to run a background for every batch of analyses in a single day using the same electrolyte.

When analyzing low concentration samples, it is absolutely necessary to run a background.


- 1. Select the same settings and parameters that will be used for the sample analyses.
- 2. Place the beaker with clean electrolyte in the analyzer unit.
- 3. **A→Start** in the Status Panel.

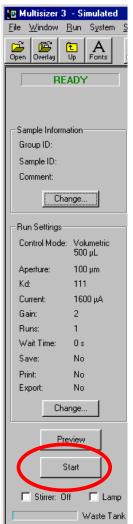

After the run is completed and saved, it can be automatically subtracted from each subsequent analysis.

## 6.7 Subtracting the Background


1. On the main menu, <sup>A</sup>→Settings, <sup>A</sup>→Load Background Run.

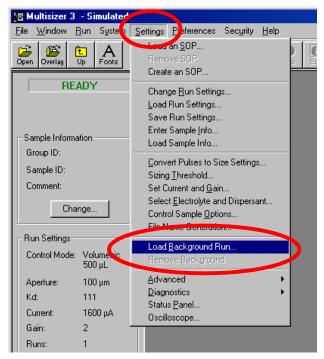

2. Select the background run.


Set Background			? ×
Look jn: 🔂 Msu		💽 🗈 🛃	<b></b>
BCKGRND_21.#M	3 🛋 BstlB.#M3	🛋 Msu.#14	🔊 MS
🛛 🛋 Bstl.#01 🛛 🛖	🛋 Msu.#07	폐 Msu.#15	🛋 MS
🛋 Bstl.#02	폐 Msu.#08	폐 Msu.#16	🔊 MS
📕 🖬 Bstl.#03 📃	폐 Msu.#10	폐 Msu.#17	🔊 MS
폐 Bstl.#04	폐 Msu.#11	폐 Msu.#18	🔊 MS
📓 BstlA.#M3	폐 Msu.#12	폐 MSU100_13.#M3	🔊 MS
			F
File <u>n</u> ame: BCKG	RND_21.#M3		<u>]</u> pen
Files of type: Multisi	zer (*.#*)		ancel
Recent <u>D</u> irectories: C:\MS\	Samples\MSU		•
		<u>G</u> et Info	

3. Double click on the background file or select the file and **<sup>A</sup>→Open**. From this point on, the background run will be automatically subtracted from each

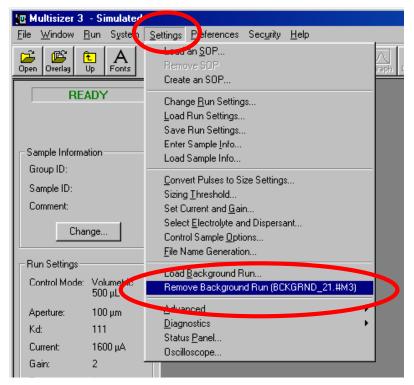

#### 6.8 <u>Running the Sample</u>

- Place the beaker with the sample in the analysis unit. <sup>⊕</sup>→Preview in the Status Panel, or <sup>⊕</sup>→Run on the Main Menu bar <sup>⊕</sup>→Preview.

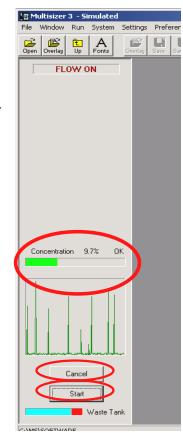

#### 6.9 Determination of Particle Concentration Using a 400µm Aperture or larger

When using a large aperture ( $400\mu$ m and larger) to determine particle concentration, it is not practical to use the volumetric control mode. If the volumetric control mode is used, the maximum volume allowed by the metering pump is 2 mL. Using a large aperture takes a very short time to pass 2 mL through the aperture. Since there is a limit for the appropriate concentration of sample, the numbers of particles counted are not enough to obtain good statistical results and therefore a poor reproducibility will achieved. The electrolyte used with large apertures should be thickened to avoid noise due to a high flow rate. It will also help in suspending heavy particles. As thickening agents, glycerol, sucrose and others can be used.

# Measure exactly the volume of electrolyte needed to fill the beaker to about 9/10 of its capacity.

- 1. Set the instrument in Time Mode to 60 seconds.
- 2. Run the electrolyte.
- 3. When the run is completed, measure accurately the volume of electrolyte left in the beaker.
- 4. Subtract the volume left from the original electrolyte volume.
- 5. Divide the resulting volume by 60. The result is the flow rate in ml/sec for this specific combination of aperture and electrolyte.
- 6. Knowing the flow rate, the Time Mode can be used to determine particle concentration. The volume of sample for the analysis is calculated using the flow rate. This volume is the "Analytical Volume".


#### 6.10 Cleaning the System

It is advisable to clean the system at the end of the day if the Multisizer 3 has been under any of the following working conditions.

- Analysis of biological samples
- Analysis of foods
- Use of corrosive fluids

#### 6.10.1 Cleaning Procedure

1. Disconnect the electrolyte reservoir from the instrument.

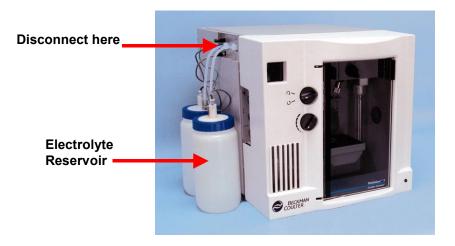

2. On the main menu, ↔→System, ↔→Drain System.

📴 Multisizer 3 - St	mulate
<u>File W</u> indow <u>R</u> un	System Settings Preferences Security Help
Open Overlay Up	Empty Waste Tank Unblock Aperture Elucit Aperture Tube
READ	<u>Fill System</u> Drain System
Group ID: Leti	<u>S</u> tirrer (on/off) Lamp (on/off)
Sample ID: Cefalo Comment:	fina Lot
Change	

- **3.** After the drain of the system is completed, in place of the electrolyte reservoir, connect a reservoir dedicated to holding the cleaning agent. The recommended cleaning agents for biological or food applications are Coulter Clenz or 2% bleach solution. For all other applications, DI water is suitable.
- **4.** On the main menu, *A*→**System**, *A*→**Fill System**.

📴 Multisizer 3 - Simulated						
<u>F</u> ile	<u>W</u> indow	<u>R</u> un	System	<u>S</u> ettings	<u>P</u> referen	nces
Dpen	Overlay	€ Up	<u>U</u> nbl	y <u>W</u> aste T ock Apertu	ire	2 e as
	R	EAD		n <u>A</u> perture ystem	Tubb	
			<u>D</u> rain	i System		
			<u>T</u> roul	bleshooting	g 🕨	
_ Sar	mple Infor	mation	<u>S</u> tirre	r (on/off)		

- 5. Repeat step 3.
- 6. Leave the system filled with the cleaning agent until the next working day.
- **7.** Before start using the instrument on the next working day, follow steps 1 trough 4, but in this case, switching back to the electrolyte.

#### Back to TOC


# SECTION 7 REVIEWING AND SAVING THE RESULTS

## 7.1 Introduction

While an analysis file is open it can be viewed, changed and saved in many different formats by using the commands and options described in this section.

## 7.2 Opening a File

To open a previously saved file, on the main menu,  $\textcircled{} \rightarrow File$ ,  $\textcircled{} \rightarrow Open$ .

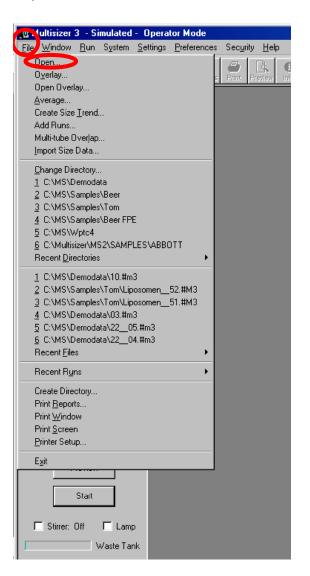

Or on the tool bar  $\mathcal{D}$  the **Open** icon.

😈 Multisizer 3 - Simulated								
<u>File W</u> indow <u>R</u> un	System	<u>S</u> ettings	<u>P</u> refer	rences	Sec <u>u</u> ri	ity <u>H</u> el	P	
Open Verlay Up	A	Cverlag	Save	<b>R</b> Save as	E Print	Preview	<b>()</b> Info	

Open			?×
Look in: 🔂	Ms	💌 🗈 💆	🖻 🔳
Demodata	_		
File <u>n</u> ame:			<u>O</u> pen
Files of type:	Multisizer (*.#*)	<b>T</b>	Cancel
	C Open as read-only		
Recent	C:\MS\MS		<b></b>
<u>D</u> irectories:			
		<u>Gi</u> et Info	0 <u>v</u> erlay
		🔽 Quick View	Size 💌

The Open command automatically displays the Open dialog box. To open the **Demodata** file,  $\textcircled{}{} \rightarrow$  the file and,  $\textcircled{}{} \rightarrow$  **Open.** 

Open			? ×
Look jn: 🔂	Demodata	🗖 🖻 🗾	<u>*</u>
🖻 01.#m3	🇃 07.#m3	폐 13.#m3	🔊 19.#mi
02.#m3 ┥	📕 📄 08.#m3	폐 14.#m3	📄 20.#m:
🛃 03.#m3	📄 09.#m3	🗾 15.#m3	🗾 21.#m:
🇃 04.#m3	🔊 10.#m3	🛋 16.#m3	🗃 22.#m:
🔄 05.#m3	🗾 11.#m3	🌌 17.#m3	🗃 23.#m:
🄊 06.#m3	🔊 12.#m3	🗃 18.#m3	🗃 24.#m:
			F
File <u>n</u> ame:	02.#m3		<u>O</u> pen
Files of <u>type</u> :	Multisizer (*.#*)	•	Cancel
- I	Open as <u>r</u> ead-only		
Recent <u>D</u> irectories:	C:\MS\Demodata		•
www.www.www.	M	<u>G</u> et Info I <u>Q</u> uick View	0 <u>v</u> erlay Size ▼

				al 📻 💶 I	
		Demodata	- <u>-</u> • <u>-</u>		
	■ 01.#m3	≥ 07.#m3 ≤ 308.#m3	i 13.#m3	)∎ 19.#mi ■ 20.#mi	
	<mark>)                                    </mark>	Image: 100.4 m3 Image: 100.4 m3 Image: 100.4 m3 Image: 100.4 m3	⊠a14.#m3 ⊛a15.#m3	i≇i 20.#m. i≇i 21.#m:	
	04.#m3	10.#m3	16.#m3	22.#m:	
	🍺 05.#m3	🍺 11.#m3	🍓 17.#m3	🗾 23.#m:	
	폐 06.#m3	🗃 12.#m3	🛋 18.#m3	🛋 24.#m:	
	•				
	File <u>n</u> ame:	02.#m3	(	<u>O</u> pen	
	Files of <u>t</u> ype:	Multisizer (*.#*)		Cancel	
	Thes of type.	· ·			
	Description	Open as read-only			
	Recent <u>D</u> irectories:	C:\MS\Demodata		•	
		MA	<u>G</u> et Info	O <u>v</u> erlay	
		y <sup>n</sup> W			
		/ \	🔽 Quick View	Size 💌	
	Whenhow when we will	W. March			
			Differential Number		
	2-		Differential Number		— 08.#m3
			Differential Number		— 08.#m3
	1.8 -		Différential Number		— 08.#m3
			Differential Number		— 08.#m3
	1.8 -		Differential Number		— 08.#m3
	1.8 - 1.6 - 1.4 -		Differential Number		— 08.#m3
[ (w)	1.8 - 1.6 - 1.4 - 1.2 -		Differential Number		— 08.#m3
	1.8 - 1.6 - 1.4 - 1.2 - 1 -		Differential Number		— 08.#m3
	1.8 - 1.6 - 1.4 - 1.2 -		Differential Number		— 08.#m3
	1.8 - 1.6 - 1.4 - 1.2 - 1 -		Differential Number		— 08.#m3
	1.8 - 1.6 - 1.4 - 1.2 - 1 - 0.8 -		Differential Number		08.#m3
	1.8 - 1.6 - 1.4 - 1.2 - 1 - 0.8 - 0.6 - 0.4 -		Differential Number		— 08.#m3
	1.8 - 1.6 - 1.4 - 1.2 - 1 - 0.8 - 0.6 -		Differential Number		08.#m3

To open any of the files,  $\Im \rightarrow$  the file name and  $\Im \rightarrow Open$ .

## 7.3 Overlaying Files

The Multisizer 3 software allows you to view more than one data file on the same graph. This is useful when you want to view point-by-point data comparisons using the same graph.

To open the data files for overlay,  $\mathcal{T} \rightarrow \mathbf{File}$  then  $\mathcal{T} \rightarrow \mathbf{Overlay}$  on the main menu, or on the tool bar,  $\mathcal{T} \rightarrow \mathbf{the Overlay}$  icon.


	🚾 Multisizer 3 - Simulated - Operator Mod	e	
	File Window Run System Settings Prefere	nces	Sec
I	Open		E,
	0 <u>v</u> erlay	s	Print
	Open a verlay	F	
	Average	- 1	
	Create Size <u>T</u> rend	- 1	
	Add Runs	- 1	
	Multi-tube Overlap Import Size Data	- 1	
	Import Size Data	- 1	
	Change Directory	- 1	
	1 C:\MS\Demodata	- 1	
	2 C:\MS\Samples\Beer	- 1	
	3 C:\MS\Samples\Tom	- 1	
	4 C:\MS\Samples\Beer FPE	- 1	
	5 C:\MS\Wptc4	- 1	
	6 C:\Multisizer\MS2\SAMPLES\ABBOTT		
	Recent <u>D</u> irectories	-	
	1 C:\MS\Demodata\10.#m3	- 1	
	<u>2</u> C:\MS\Samples\Tom\Liposomen52.#M3	- 1	
	<u>3</u> C:\MS\Samples\Tom\Liposomen51.#M3	- 1	
	4 C:\MS\Demodata\03.#m3	- 1	
	5 C:\MS\Demodata\2205.#m3	- 1	
	6 C:\MS\Demodata\22_04.#m3		
	Recent <u>F</u> iles		
	Recent R <u>u</u> ns		
	Create Directory		
	Print <u>B</u> eports		
	Print <u>W</u> indow		
	Print <u>S</u> creen		
	Printer Setup		
	E <u>x</u> it		

📴 Multisizer 3 - Simulated										
<u>F</u> ile <u>W</u>	indow	<u>B</u> un	System	<u>S</u> ettings	<u>P</u> refei	rences	Sec <u>u</u> ri	ty <u>H</u> el	P	
Open O	)verlay	Ê Up	A	Cverlag	Save	N. Save as	E Print	Preview	<b>f</b> Info	


The **Overlay** dialog box will be displayed.

Overlay			×
<u>R</u> ecent Directories:	C:\MS\Demodata		•
File <u>N</u> ame: <b>****</b> Directories:	Multisizer (*.#*) 28 Files:	Open as Read- <u>o</u> nly     O Selected Files:	
[.] [-a-] [-d-]	01. #m3 02. #m3 03. #m3 04. #m3 05. #m3 06. #m3 07. #m3 08. #m3 09. #m3 10. #m3 11. #m3 11. #m3 13. #m3		
<u>U</u> pdate	Giet [nfo >>> ≜dd >	> << Remo <u>v</u> e << Car	ncel OK


To **Overlay** files,  $\neg \ominus \rightarrow$  to highlight the file name.  $\neg \ominus \rightarrow Add$  in the bottom of the dialog box. Repeat this step to select all the files to overlay. When you have finished choosing the files,  $\neg \ominus \rightarrow OK$ . To overlay consecutive files,  $\neg \ominus \rightarrow to$  highlight the first file name, hold down the left mouse button and drag to highlight all the required files then  $\neg \ominus \rightarrow Add$ , and  $\neg \ominus \rightarrow OK$ .

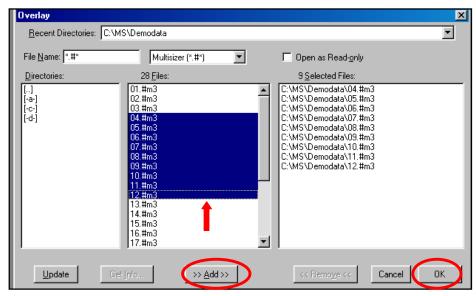

A graph containing data from all selected files is displayed. The legend that appears in the upper right-hand corner of the graph displays the names of the data files used in the overlay.

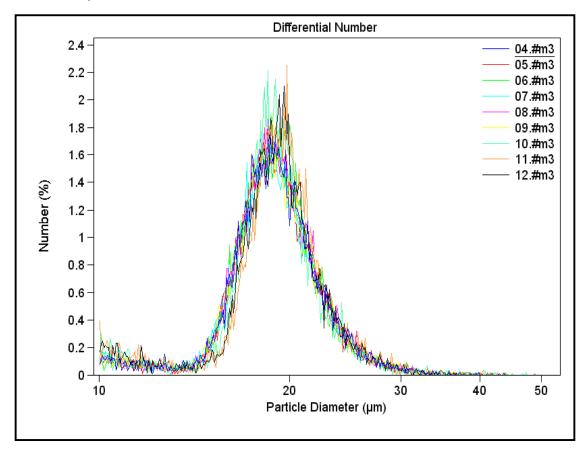

To add an additional file to an already existing overlay,  $\mathcal{D} \rightarrow \mathbf{Run}$  File, then  $\mathcal{D} \rightarrow \mathbf{Open}$  for **Overlay** on the second level menu, or  $\mathcal{D} \rightarrow \mathbf{Overlay}$  (second icon) on the tool bar.

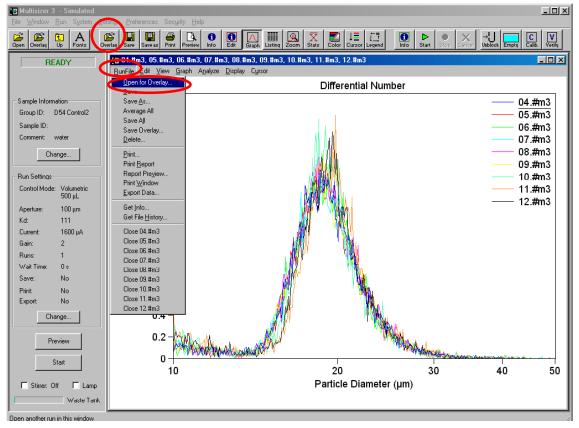


The **Open for Overlay** dialog box will appear automatically. Select the file to be added to the existing overlay and  $\Im \rightarrow Open$ .

Open for Overla	y.		? ×
Look jn: 🔂 D	emodata	- 🗈 💋	* 📰 🖿
🔊 01.#m3	🛋 07.#m3	폐 13.#m3	🔊 19.#m:
🗃 02.#m3	🍺 08.#m3	🍺 14.#m3	🙍 20.#m:
03.#m3 🔶	💻 🛋 09.#m3	連 15.#m3	🔊 21.#m
04.#m3	🎑 10.#m3	🎑 16.#m3	🧃 22.#m:
05.#m3	連 11.#m3	🎑 17.#m3	🔊 23.#mi
i 06.#m3	🔊 12.#m3	🄊 18.#m3	🔊 24.#mi
•			Þ
File <u>n</u> ame:	3.#m3	(	<u>O</u> pen
Files of type:	/ultisizer (*.#*)		Lancel
Γ	Open as <u>r</u> ead-only		
Recent C: Directories:	\MS\Demodata		•
hd.	Mu	<u>G</u> et Info	Add to <u>W</u> indow
Ĵ	<b>N</b>	🔽 Quick View	Size 💌
Merson and a stand	My and a hard a second		

The selected file will be added to the existing overlay.

The **Overlay** can be saved as an **Overlay File**. To save an **Overlay**, on the second level menu,  $\mathcal{P}$ -**Run File**,  $\mathcal{P}$ -**Save Overlay**.

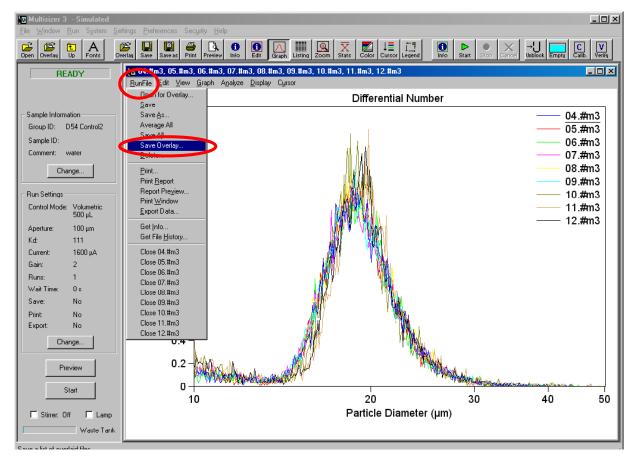

The Save Overlay dialog box will appear automatically. Assign a name for the Overlay File and  ${}^{\circ}$   $\rightarrow$  Save.

Save a list of o	verlaid files			? ×
Save in: 🔂 D	emodata	•		* 🔳
🔊 Overlay4-12.0	IVO			
File <u>n</u> ame:	)verlay4-12.ovr		_(	<u>S</u> ave
Save as <u>t</u> ype:	)verlay File (*.ovr)			Cancel
Recent Directories:	\MS\Demodata			•
Directones: )				

🚾 Multisizer	3 - Si	mulated	- Opera	tor Mode	
<u>File</u> indow	<u>R</u> un	System	<u>S</u> ettings	<u>P</u> reference	s Sec <u>i</u>
<u>0</u> pen					
Overlau					s Print
Open Over	lay				
<u>Average</u>					
Create Size	_				
Add Runs.					
Multi-tube 0					
Import Size	Data				
<u>C</u> hange Dir	ectory				
<u>1</u> C:\MS\D	emoda	ita			
<u>2</u> C:\MS\S					
<u>3</u> C:\MS\S					
<u>4</u> C:\MS\S		NBeer FP	E		
<u>5</u> C:\MS\V					
<u>6</u> C:\Multis			LESVABBO	TTC	
Recent <u>D</u> ire	ectories	;		<u> </u>	
<u>1</u> C:\MS\D	emoda)	ita\10.#m	3		
2 C:\MS\S		•			
<u>3</u> C:\MS\S		•		51.#M3	
<u>4</u> C:\MS\D					
5 C:\MS\D					
6 C:\MS\D		ita\220	4. <b>⊞</b> m3		
Recent <u>F</u> ile	s			•	
Recent R <u>u</u>	ns			•	
Create Dire	ctory				
Print <u>R</u> epor	ts				
Print <u>W</u> indo	W				
Print <u>S</u> cree	n				
<u>P</u> rinter Setu	ю				
E <u>x</u> it					

To open an **Overlay File**, on the main menu, P **File**, P **Open Overlay**.

The **Open Overlay** dialog box will appear automatically. Select the file and  $\textcircled{} \rightarrow \textbf{Open}$ . The file will be added to the **Overlay**.

Open Overla	у				? ×
Look jn: 🧲	Demodata		-	<u></u>	<u>r i i i i i i i i i i i i i i i i i i i</u>
Overlay4-	12.ovr				
File <u>n</u> ame:	Overlay4-12.ovr				<u>O</u> pen
Files of type:	Overlay File (*.ov	/r)		⊡	Lancel
Recent Directories:	C:\MS\Demodata	1			•
Directones.					

This file is linked to the source files. It means that the file may only be opened when the original files are in the location from where the overlay file was saved.

## 7.4 Averaging Files

The averaging function allows data from multiple data files to be averaged together. Data from individual files are combined, channel contents are averaged, sample statistics for each channel are calculated and the distribution is plotted on a single graph. The resultant file provides a graphical representation of the averaged contents of each channel.


i a							_
	u Multisizer	3 - Si	imulated	- Opera	tor Mode		
	<u>F</u> ile <u>W</u> indow	<u>R</u> un	System	<u>S</u> ettings	Preference	es S	ec
Ì	<u>0</u> pen						e.
l	O <u>v</u> erlay					s P	
I	Opt., Over	ay				H	_
	Average						
	Create Size	<u>T</u> rend					
	Add Runs						
	Multi-tube 0	)verlap					
	Import Size	Diata					
	<u>C</u> hange Dir	ectory.					
	1 C:\MS\D						
	2 C:\MS\S	amples	s∖Beer				
	3 C:\MS\S	amples	NT om				
	4 C:\MS\S	amples	NBeer FP	E			
	5 C:\MS\V	Vptc4					
	<u>6</u> C:\Multis	izer\M	52\SAMP	LESVABBO	TT		
	Recent <u>D</u> ire	ectories	;		•		
	1 C:\MS\D	)emoda	ata\10.#m	3			
	2 C:\MS\S	amples	NTom\Lip	osomen	52.#M3		
	3 C:\MS\S	amples	NTom\Lip	osomen	51.#M3		
	4 C:\MS\D	emoda	ata\03.#m	3			
	5 C:\MS\D	)emoda	ata\220	5. <b>#</b> m3			
	<u>6</u> C:\MS\D	emoda	ata\220	I4.#m3			
	Recent <u>F</u> ile	s			+		
	Recent R <u>u</u>	ns			•		
	Create Dire	ctory					
	Print <u>R</u> epor	ts					
	Print <u>W</u> indo	W					
	Print <u>S</u> cree	n					
	<u>P</u> rinter Setu	ю					
	E <u>x</u> it						


To average files, on the main menu,  $\textcircled{} \rightarrow File$ ,  $\textcircled{} \rightarrow Average$ .

The Average dialog box will appear automatically.


Average			×
<u>R</u> ecent Directories:	C:\MS\Demodata		•
File <u>N</u> ame: .#*	Multisizer (*.#*)	Dpen as Read-only	
Directories:	28 <u>F</u> iles:	0 <u>S</u> elected Files:	
[] [-a-] [-d-] [-d-] [-h-] [-h-] [-h-] [-l-] [-y-]	<ul> <li>■ 01.#m3 02.#m3 03.#m3 04.#m3 05.#m3 06.#m3 06.#m3 07.#m3 08.#m3 09.#m3</li> </ul>		
<u>U</u> pdate	Gret Info >> Add >>	<< Remo <u>v</u> e << 1	Cancel OK

For averaging files,  $\textcircled{} \rightarrow$  file name.  $\textcircled{} \rightarrow$  Add in the bottom of the dialog box. Repeat this step to select all the files to overlay. When you have finished choosing the files,  $\textcircled{} \rightarrow$ OK. To **Average** consecutive files,  $\textcircled{} \rightarrow$ OK highlight the first file name, hold down the mouse left mouse button and drag to highlight all the required files,  $\textcircled{} \rightarrow$ Add and then  $\textcircled{} \rightarrow$ OK.

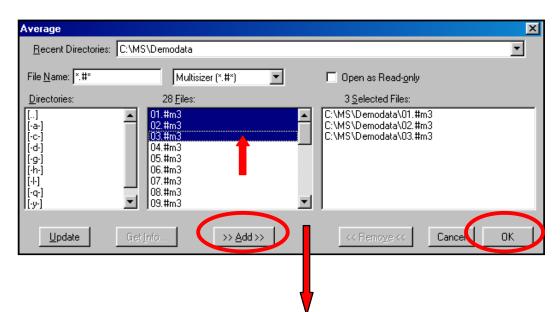

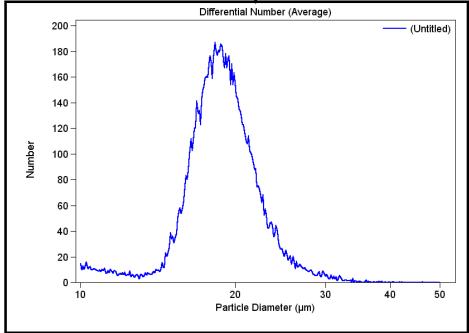


To add an additional file to an already existing **Average**, on the second level menu,  $\mathcal{D} \rightarrow \mathbf{Run \ File}$ ,  $\mathcal{D} \rightarrow \mathbf{Open \ and \ Add \ to \ Average}$ .

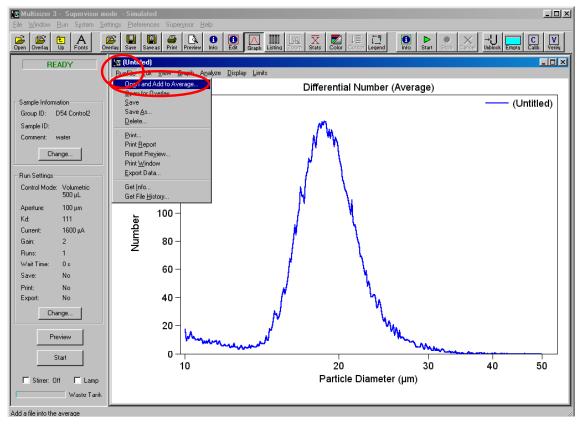


The **Open** dialog box will appear automatically. Select the file and  ${}^{\circ} \rightarrow$ **Open**. The file will be added to the **Average**.

Open			? ×
Look jn: 🔂 Dem	odata	- 🗈 💋	* 📰 📰
🛋 01.#m3	🛋 07.#m3	🛋 13.#m3	🛋 19.#m:
🗃 02.#m3	폐 08.#m3	🍙 14.#m3	🗃 20.#m:
🔊 03.#m3	폐 09.#m3	폐 15.#m3	폐 21.#m:
04.#m3	폐 10.#m3	폐 16.#m3	🛋 22.#m:
🗃 05.#m3	폐 11.#m3	폐 17.#m3	폐 23.#m:
폐 06.#m3	폐 12.#m3	폐 18.#m3	폐 24.#m:
Files of type: Mul	#m3 ltisizer (*.#*)		Dpen Cancel
Directories: C:\M	S\Demodata		<u> </u>
and the second of the second o	My how we want the second seco	<u>G</u> et Info I Quick View	Size 💌

The Average file can be saved. On the second level menu, A→Run File, A→Save.

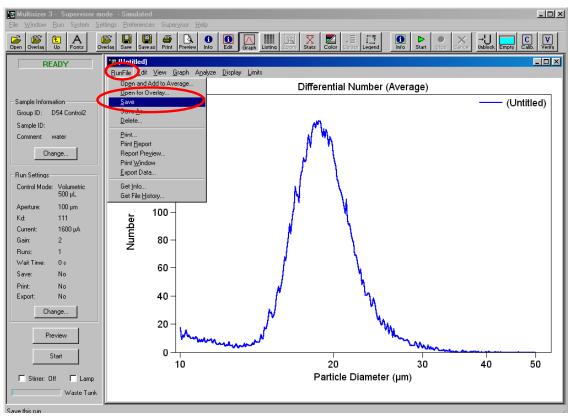

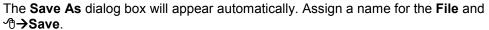

Save As			? ×
Savejn: 🔂 Der	nodata	💿 🖻 💆 🖸	* 🔳
🔊 01.#m3	🔊 07.#m3	폐 13.#m3	🔊 19.#mi
🙍 02.#m3	폐 08.#m3	🐋 14.#m3	🙍 20.#m:
🛋 03.#m3	폐 09.#m3	🐋 15.#m3	폐 21.#mi
🔊 04.#m3	🎽 10.#m3	連 16.#m3	🗾 22.#m:
🔊 05.#m3	🎽 11.#m3	🌌 17.#m3	🔊 23.#mi
🔊 06.#m3	🄊 12.#m3	폐 18.#m3	🄊 24.#m
•			Þ
File <u>n</u> ame:		>	<u>S</u> ave
Save as type: Mu	ultisizer (*.#*)	•	Cancel
Γ	Save as <u>r</u> ead-only		
Recent <u>D</u> irectories:	/IS\Demodata		•
	No.	<u>G</u> et Info I Quick View	Size 💌
Mary and and a second s	· · · · · · · · · · · · · · · · · · ·		

## 7.5 Creating a Size Trend

**Size Trend** is most often used to show statistics from a number of sample runs on one graph or report. This would be useful, for example, in a manufacturing process where a sample is measured with the Multisizer at specific time intervals, (minutes, hours, days etc.). These data files can be used to create a trend graph or listing at the end of a chosen period of time to track changes of selected statistical values. User-selectable pass/fail limits can also be shown on trend graphs and listings. The Size Trend can be updated at any time by adding new points.


To create a Size Trend, on the main menu, <sup>⊕</sup>→File, <sup>⊕</sup>→Create Size Trend

🚾 Multisizer 3 - Simulated - Operator Mode		
<u>File W</u> indow <u>R</u> un System <u>S</u> ettings <u>P</u> referen	ces	Sec <u>u</u>
Dpen O⊻erlay Open Overlay Average	s	Print
Create Size <u>T</u> rend		
Add Bure Multi-tube Over]ap Import Size Data		
Change Directory 1 C:\MS\Demodata 2 C:\MS\Samples\Beer 3 C:\MS\Samples\Tom 4 C:\MS\Samples\Beer FPE 5 C:\MS\Wptc4		
<u>6</u> C:\Multisizer\MS2\SAMPLES\ABBOTT Recent <u>D</u> irectories		
C:\MS\Demodata\10.#m3     C:\MS\Samples\Tom\Liposomen_52.#M3     C:\MS\Samples\Tom\Liposomen_51.#M3     C:\MS\Demodata\03.#m3     C:\MS\Demodata\22_05.#m3     C:\MS\Demodata\22_04.#m3		
Recent <u>F</u> iles		
Recent R <u>u</u> ns		
Create Directory Print <u>B</u> eports Print <u>W</u> indow Print <u>S</u> creen <u>P</u> rinter Setup		
Exit		


The Create Size Trend dialog box will appear automatically.

Create Size Trend			X
Recent Directories:	C:\MS\Demodata		•
File <u>N</u> ame: <b>*:#*</b>	Multisizer (*.#*)	Open as Read-only	
<u>D</u> irectories:	28 <u>F</u> iles:	0 <u>S</u> elected Files:	
[] [-a-] [-c-] [-d-] [-g-] [-h-] [-h-] [-h-] [-y-]	<ul> <li>■ 01. #m3 02. #m3 03. #m3 04. #m3 05. #m3 06. #m3 06. #m3 07. #m3 08. #m3</li> <li>■ 09. #m3</li> </ul>		
<u>U</u> pdate	Get Info >> Add >>	<< Remove <<	Cancel OK.


To select the files to be included in the **Size Trend**,  $\neg \ominus \rightarrow$  the file name. Once you have highlighted the file name,  $\neg \ominus \rightarrow Add$  in the bottom of the dialog box. Repeat this step to select all the files for the **Size Trend**. When you have finished selecting the files,  $\neg \ominus \rightarrow OK$ . To select consecutive files,  $\neg \ominus \rightarrow$  to highlight the first file name, hold down the left mouse button and drag to highlight all the required files.  $\neg \ominus \rightarrow Add$ , and  $\neg \ominus \rightarrow OK$ .

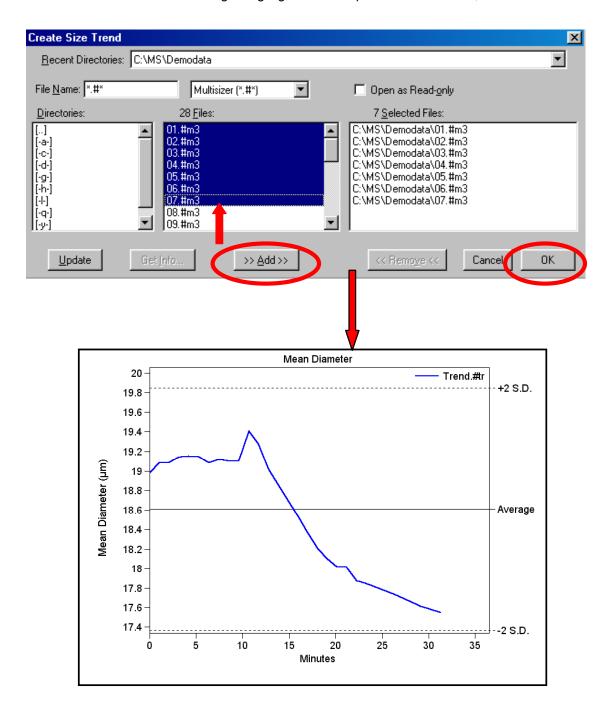


Additional files may be added to an already existing **Size Trend**. On the second level menu,  $\mathcal{T} \rightarrow \mathbf{Run}$  **File**,  $\mathcal{T} \rightarrow \mathbf{Open}$  and **Add to Trend**.

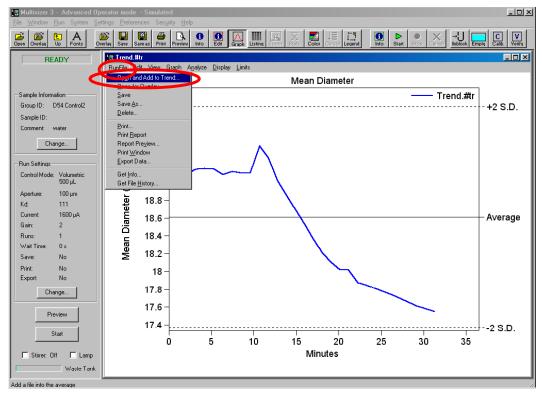

The **Open** dialog box will appear automatically. Select the file and  $\textcircled{O} \rightarrow Open$ . The file will be added to the **Size Trend**.

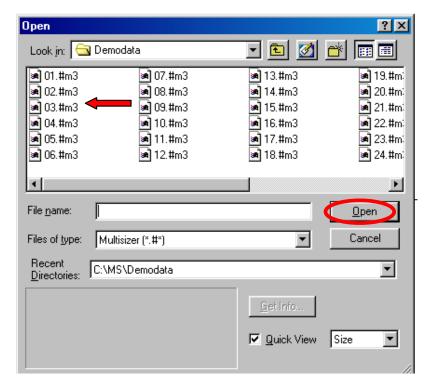

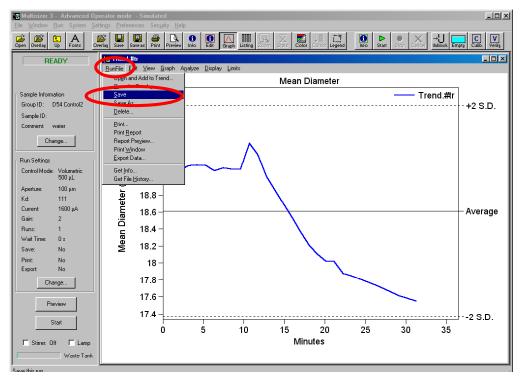


The **Trend** file can be saved. On the second level menu,  $\mathcal{D} \rightarrow \mathbf{Run}$  File,  $\mathcal{D} \rightarrow \mathbf{Save}$ .

Save As			? ×
Savejn: 🔂 De	modata	- 🗈 🗹 c	* 🔳
🛋 01.#m3	🛋 07.#m3	🛋 13.#m3	🛋 19.#mi
🛃 02.#m3	폐 08.#m3	🍺 14.#m3	폐 20.#m:
🛃 03.#m3	폐 09.#m3	폐 15.#m3	폐 21.#m:
🛋 04.#m3	폐 10.#m3	🛋 16.#m3	🛋 22.#m:
🛋 05.#m3	🛋 11.#m3	🛋 17.#m3	🛋 23.#mi
🛋 06.#m3	폐 12.#m3	🛋 18.#m3	🛋 24.#m:
•			
File <u>n</u> ame:	end.#tr		<u>S</u> ave
Save as type: M	ultisizer (*.#*) Save as <u>r</u> ead-only		Lancel
Recent Directories:	MS\Demodata		•
		<u>G</u> et Info I▼ <u>Q</u> uick View	Size 💌

## 7.6 Overlapping Files from Different Aperture Sizes.

The **Multi-Tube Overlap** procedure is used to combine up to 5 distributions into a single distribution with a new file name. This merging routine is able to create particle size distributions outside the range of any single aperture by combining 2 to 5 aperture size analyses from the same sample.

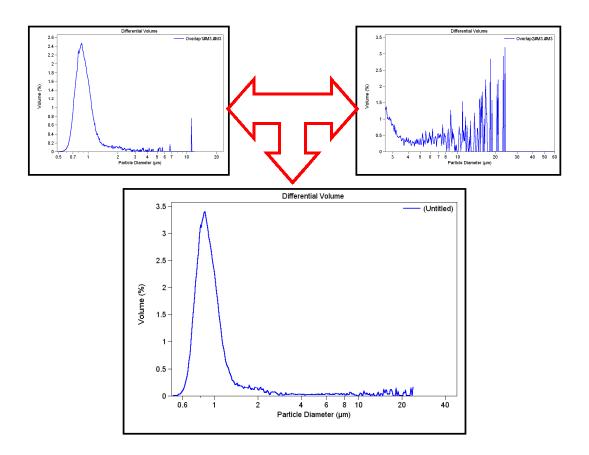

Obtain a size distribution on the Multisizer analyzer with an aperture tube suitable for the coarsest material in the sample. After analysis, sieve, filter or use any other suitable method to remove particles from the remaining suspension which are likely to block the smaller aperture. Continue the analysis with the next smaller aperture.


To find the combined distribution, find the point or range where the plots are most "parallel". This point, or range, is called the "offset point". At this point the vertical displacement (or offset), between the two graphs is measured. For example, if at the chosen point the cumulative number oversize is 1,000 from the small aperture and 10,000 from the large aperture, then the dilution is exactly 10:1.To combine the two plots the Multisizer 3 program multiplies the sample data from the small aperture by 10 and overplots it on the large aperture plot. The program will then exclude the small aperture data with a size larger than the offset point. The result is the combined distribution.


To create a Multi-Tube Overlap file, on the main menu,  $\mathcal{D} \rightarrow File$ ,  $\mathcal{D} \rightarrow Multi-Tube Overlap$ .

The **Multi-Tube Overlap** dialog box will appear automatically. Up to five files may be selected to be included in the **Multi-Tube Overlap**. To select files  $\textcircled{O} \rightarrow$  the file name. Once you have highlighted the file name,  $\textcircled{O} \rightarrow Add$  in the bottom of the dialog box. Repeat this step to select all the files for the **Multi-Tube Overlap**. When you have finished selecting the files,  $\textcircled{O} \rightarrow OK$ . To select consecutive files,  $\textcircled{O} \rightarrow to$  highlight the first file name, hold down the left mouse button and drag to highlight all the required files,  $\textcircled{O} \rightarrow Add$ , and  $\textcircled{O} \rightarrow OK$ .

Multi-tube Overlap		×
<u>R</u> ecent Directories:	C:\MS\Demodata	<b>•</b>
File <u>N</u> ame: <b>*.</b> #*	Multisizer (*.#*)	Open as Read- <u>o</u> nly
<u>D</u> irectories:	28 <u>F</u> iles:	2 <u>S</u> elected Files:
[] [-a-] [-c-] [-d-]	20.#m3 21.#m3 22.#m3 23.#m3 24.#m3 Overlap1#M3.#M3 Overlap2#M3.#M3 Ovp#m3.#m3 Trend.#tr	C:\MS\Demodata\0verlap1#M3.#M3 C:\MS\Demodata\0verlap2#M3.#M3
<u>U</u> pdate	Giet Info	K Remove K Cancel OK




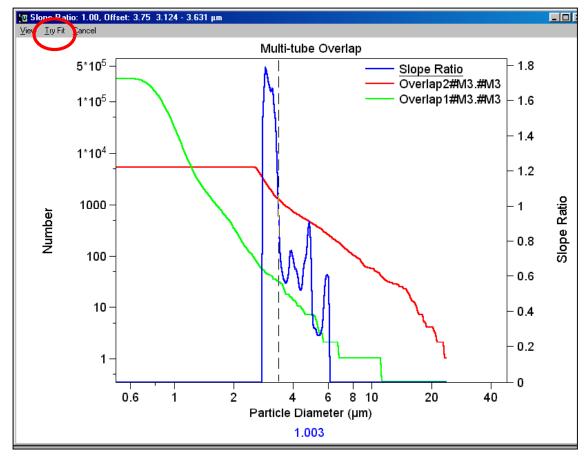


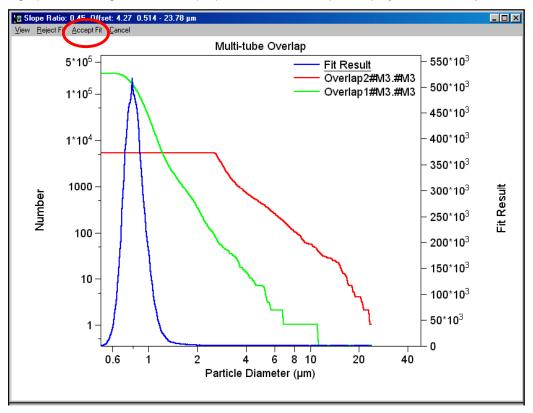

A graph containing the software proposal for the overlap is displayed. ↔ Accept Fit.

The size distribution graph for the resultant file will appear.

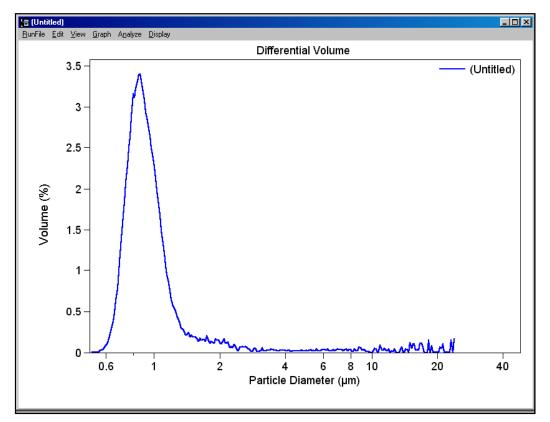

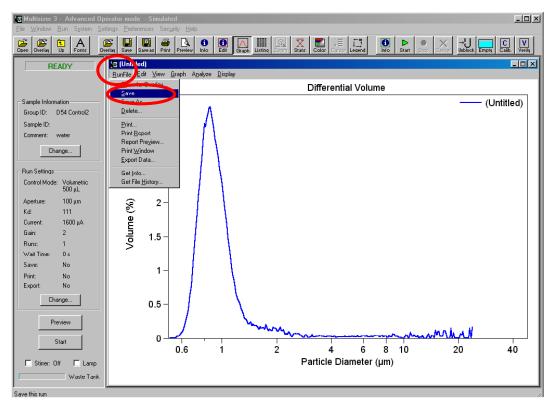

To save the file, on the second level menu,  $\mathcal{D} \rightarrow \mathbf{Run \ File}$ ,  $\mathcal{D} \rightarrow \mathbf{Save}$ .

Save As			? ×
Save in: 🔁	Demodata	- 🗈 💆 🕻	
🍺 01.#m3	폐 07.#m3	🍺 13.#m3	🍺 19.#m:
🔊 02.#m3	🍺 08.#m3	🍺 14.#m3	🍺 20.#m:
🍺 03.#m3	폐 09.#m3	폐 15.#m3	폐 21.#m:
🍺 04.#m3	폐 10.#m3	폐 16.#m3	폐 22.#m:
🍺 05.#m3	🍺 11.#m3	폐 17.#m3	폐 23.#m:
🍺 06.#m3	폐 12.#m3	폐 18.#m3	폐 24.#m:
			ľ
File <u>n</u> ame	Ovp#m3.#m3		<u>S</u> ave
Save as <u>t</u> ype:	Multisizer (*.#*)	•	Cancel
	Save as <u>r</u> ead-only		
Recent <u>D</u> irectories:	C:\MS\Demodata		•
$\square$		<u>G</u> et Info	
		☑ Quick View	Size 💌

**Back to TOC** 

## **SECTION 8 PRINTING A REPORT**

#### 8.1 Introduction

This section describes how to set up the **Preferences Menu** in order to customize the printed report from the analysis. **The Preferences Menu** is found on the main menu, and after selecting the desired options, they can be saved as a preference file or as the **Default Preferences** file. Saving these files allows the user to recall them in the future and also to be used in the creation of a **Standard Operation Procedure (SOP)** (see <u>Section 9</u>).

The **Default Preferences** file will determine how the format of the on-screen graph and data is displayed when the program is started at the beginning of a working session.


The **Preferences File** in use will determine the format of the on-screen graph and data displayed in the following situations:

- During the length of the analysis.
- After the analysis is completed.
- When a file is opened

Samples of data files, chosen to illustrate how to work with them, are included in the software's **Demodata** directory and are installed in **C:/MS/Demodata**.

The way the on-screen results are displayed can be changed at any time while the file for the run is open, but the printed report will be changed only if the Preferences File is altered.

To access the **Preferences Menu**, go to the main menu and *<sup>o</sup>*→**Preferences**.

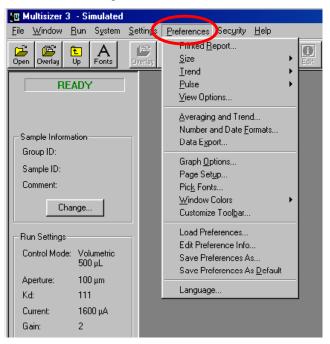

## 8.2 Selecting the Content of the Printed Report

The **Printed Report** submenu is where the choices for the content of the report are found. Once the selection of the content is done, they can be customized on the subsequent submenus.

In the **Preferences Menu**, *A*→**Printed Report.** 

## 8.2.1 Sample Info.

Checking the **Sample Info** box allows you to place basic information about the analysis on the report. The sample information enterd by the operator in <u>Section 6.5</u> can be reported in two formats, standard and short.

	Printed Report	×
Printed Report         Size         Irend         Pulse         View Options         Averaging and Trend         Number and Date Formats         Data Export	Sample Into Short Size Graphs Select Statistics Averaged Statistics Compare Statistics	Clear All       OK         Cancel         Size Statistics         Mass       Select         Number       ℤ % □ /mL
Graph Options Page Setup Pick Fonts Window Colors Customize Toolbar Load Preferences Edit Preferences As Save Preferences As Save Preferences As	Listing Select Listing Select Pulses Graph Select Listing Select Printer	Volume     %     /mL       Surface Area     %     /mL       Irend     %     /mL       Graph     Select       Listing     Select
Language	<ul> <li>Use Default Printer</li> <li>Use Specified Printer:</li> </ul>	Select

Example for standard Sample Info:

Multisizer 33.5120 May 2001File name: 14931s1.#01Group ID: 14931s1Sample ID: A11Operator: ALRun number: 1
File name: 14931s1.#01 Group ID: 14931s1 Sample ID: A11 Operator: AL
Group ID: 14931s1 Sample ID: A11 Operator: AL
Sample ID: A11 Operator: AL
Operator: AL
•
Dup pumber: 1
Electrolyte: 4% NaCl IN H2O
Dispersant: None
Aperture diameter: 30 µm Kd: 42.393
Aperture current: -800 µA Gain: 4
Size bins: 256 from 0.6 µm to 18 µm
Sigma: 7087 (Coincidence corrected)
Count > 0 $\mu$ m: 0 Coincidence corrected: 0
Coincidence correction: 0.0%
Control method: Volumetric, 500 µL
Acquired: 9:19 30 Mar 2000
Electrolyte volume: 20 mL
Analytic volume: 500 µL
Sample: 2.5 mL
Background run: 14931.#01

Example for short **Sample Info**:

 Multisizer 3
 3.51

 20 May 2001
 File name: 14931s1.#01

 Group ID: 14931s1
 Sample ID: A11

 Size bins: 256 from 0.6 μm to 18 μm
 Sigma: 7087 (Coincidence corrected)

 Acquired: 9:19 30 Mar 2000
 Background run: 14931.#01

## 8.2.2 Size Section of the Printed Report

#### 1. Size Graph

By checking the Graph box, the size distribution histogram for the samples will appear on the Printed Report. Different options are available for the size distribution **Graph**. To select these options,  $\neg \ominus \rightarrow$  **Select** and a dialog box containing all the options for the Graph will pop up. After selecting the format of the **Graph**,  $\neg \ominus \rightarrow OK$ .

Printed Report	×					
Sample In <u>f</u> o	Clear <u>A</u> ll OK					
Select	Cancel					
	Mass Select					
	Number 🔽 % 🗖 /mL					
	/olume □% □ /mL					
Interpolation Select	Surface Area 🗖 % 🗖 /mL					
Pulses	d					
Graph Select	Graph Select					
Listing Select	_istingSelect					
Printer						
<ul> <li>Use Default Printer</li> </ul>	Select					
C Use Specified Printer:	Size Graphs					×
	Distribution Types	(Select up t	o 12)———			
		Mass N	umber	Volume	Surface Area	OK
	Differential:		% /mL ☑ □			Cancel
	_					
	<u>C</u> umulative <:					
	C <u>u</u> mulative >:					Clear <u>A</u> ll
	Djff. + Cum. <:					
	Diff. + Cum. >:					
	Graph Printing Or	der				
	Diff. Number %		- <sub>E</sub> XA	xis	🔲 Log Y Axis	
	Diff. Volume 炎			Linear	<u>Multisizer Curs</u>	ore
				Log		
				As Accumulated	Y/dX Diff. Grap	
				As Accumulated	🔲 Use Sample ID	) as Title

#### 2. Size Statistics

The **Statistics** for the samples will be contained in the Printed Report by checking the corresponding box. In Section 8.3 is described how to select the statistics parameters to be shown on the report.

Printed Report	×
I Sample In <u>f</u> o I Short	Clear <u>A</u> ll
Size	Cancel
Statistics	Size Statistics
Averaged Statistics	Mass Select
Compare Statistics	🗖 Number 🛛 🖉 % 🗖 /mL
Listing Select	□ Volume □ % □ /mL
Interpolation Select	🗖 Surface Area 🗖 % 🗖 /mL
P <u>u</u> lses	_ <u>I</u> rend
Graph Select	Graph Select
Listing Select	Listing Select
<u>P</u> rinter	
Use Default Printer	Select
C Use Specified Printer:	

#### 3. Averaged Statistics

When several runs are averaged, the statistics of the resultant run can be obtained on the printed report by selecting the **Averaged Statistics** option.


Printed Report	<u>×</u>
Sample In <u>f</u> o	Clear <u>A</u> ll
Size Graphs Select	Cancel
Averaged Statistics	Mass Select
Compare Statistics	🗖 Number 🛛 🖉 % 🗖 /mL
Listing Select	□ Volume □ % □ /mL
☐ InterpolationSelect	🗂 Surface Area 🗖 🖏 🗂 /mL
P <u>u</u> lses	_ <u>I</u> rend
Graph Select	Graph Select
Listing Select	Listing Select
_ <u>P</u> rinter	
O Use Default Printer	Select
C Use Specified Printer:	

## 4. Compare Statistics

The **Multisizer 3** software allows one data file to be used as a statistical standard to which other data files may be compared. Limits from this standard file can also be applied to other data files.

To use the **Compare Statistics** function set up, as described below, a data file to serve as a standard file against which other files will be compared. When a sample is selected as control and its statistics are entered in the **Sample Info** box, this run can be used as a reference to compare to other runs (see <u>Section 6.5</u>). If a file from a sample was not selected as **Control Sample** at the time of the analysis, and it is required to **Compare Statistics**, the following procedure can be used.

- 1. Use the **Open** command as described in <u>Section 7.2</u> to open the required file.
- 2. On the second level menu,  $\mathcal{T} \rightarrow \mathsf{Edit}$ ,  $\mathcal{T} \rightarrow \mathsf{Edit}$  Sample Statistics.

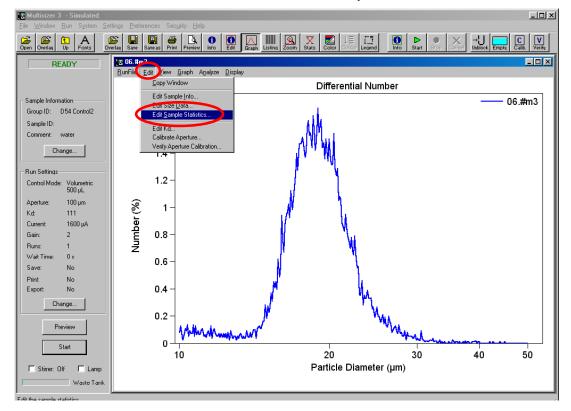

3. In the **Edit Sample Statistics** dialog box, enter the statistics information: *√*θ→**OK**. Save the file.

Edit San	ple Statistic	:\$		×
	Nominal Value	Lower Limit	Upper Limit	ОК
Mean:	Ē.	0	0	Cancel
Median:	0	0	0	
Mode:	0	0	0	
S.D.:	0	0	0	
C.V.:	0	0	0	%

# To use the Sample Statistics as a reference, save this file under an easily identifiable name.

To display on the **Printed Report** the **Compare Statistics**, check the appropriate box in the **Printed Report** dialog box.

Printed Report	<u>&gt;</u>
✓ Sample In <u>f</u> o ✓ Short	Clear <u>A</u> ll
Size	Cancel
🔽 Statistics	Size Statistics
Averaged Statistics	Mass Select
Compare Statistics	🗖 Number 🔽 🛠 🗖 /mL
Listing Select	□ Volume □ % □ /mL
Interpolation Select	🗖 Surface Area 🗖 % 🗖 /mL
P <u>u</u> lses	_ <u>I</u> rend
Graph Select	Graph Select
Listing Select	Listing Select
Printer	
Use Default Printer	Select
C Use Specified Printer:	

#### 5. Size Listing

By selecting the **Listing** option, a section containing a list of the data for each channel of the size distribution will be placed on the **Printed Report**.

Printed Report	×
Sample In <u>f</u> o	Clear <u>A</u> ll
<u>S</u> ize ✓ GraphsSelect	Cancel
✓ Statistics	Size Statistics
Averaged Statistics	Mass Select
Compare Statistics	🗖 Number 🛛 🖉 % 🗖 /mL
Listing Select	□ Volume □ % □ /mL
Interpolation Select	🗖 Surface Area 🗖 % 🗖 /mL
P <u>u</u> lses	_ <u>I</u> rend
Graph Select	Graph Select
Listing Select	Listing Select
Printer	
Use Default Printer	Select
C Use Specified Printer:	

The format and content of the Listing can be selected by →Select next to the Listing box in the **Printed Report** dialog box.

The Size Listing dialog box appears automatically. Select the desired options, then  $\textcircled{} \rightarrow OK$ .

Size Listing					X
- Columns (Select up to 7)					
☑ Bin <u>N</u> umber	Mass	Number %/mL	Volume %/mL	Surface Area % /mL	
🔽 Lower Edge 🛛 💿 Diamete	er <u>D</u> iff.: 🗖				Cancel
🗌 🗖 Center 🛛 🔿 Volume	<u>C</u> um. <: 厂				
🔲 Upper Edge 🛛 Surface	Area C <u>u</u> m.≻: ⊫			ГГГ	Clear <u>A</u> ll
Bin Number Bin Diameter (Lower) µm	Diff. Diff. Number Volum え え	e			
Bin Grouping	<b>.</b>				
<u>G</u> roup Size: 1	C All Bins				
O First Bin in Group	Between Cursors     Between Cursors     Addition     Section     Secti				
Sum Bins in Group	● From 1 to	1			

#### 6. Size Interpolation

The **Multisizer 3** software can be used to create sample data reports that will arrange data to interpolate data to conform to a set of user defined boundaries, also called "channels", "bins", or "size classes". (This is similar to how particle data points are reported when different size sieves are used.) The Multisizer 3 software refers to this process as **Size Interpolation**.

To include on the **Printed Report** the **Interpolation Points**, check the appropriate box in the **Printed Report** dialog box.

Printed Report	×
I Sample In <u>f</u> o I Short	Clear <u>A</u> ll
Size	Cancel
✓ Statistics	Size Statistics
Averaged Statistics	Mass Select
Compare Statistics	🗖 Number 🔽 🎖 🗖 /mL
Listing Select	□ Volume □ % □ /mL
Interpolation Select	🗖 Surface Area 🗖 % 🗖 /mL
P <u>u</u> lses	Irend
Graph Select	Graph Select
Listing Select	Listing Select
_ <u>P</u> rinter	
Use Default Printer	Select
O Use Specified Printer:	

The Size Interpolation dialog box appears automatically. When finished selecting the options,  $\neg \oplus \rightarrow OK$ .

Size Interpolation	×
Interpolation <u>Type</u> ⓒ <u>Size to</u> ] ☐ <u>N</u> umber ☐ % ☑ /mL ☐ <u>V</u> olume ☐ % ☐ /mL ☐ <u>S</u> urface Area ☐ % ☐ /mL ⓒ to Size	Interpolation Points ↓ Add > 10 25 50 75 90 Delete
<ul> <li>Under (&lt;)</li> <li>Over (&gt;)</li> <li>Differential</li> </ul>	Bange         C All         Setween Cursors         C From       0         μm       to

## 8.2.3 Size Statistic Section of the Printed Report

This section of the Printed Report dialog box allows a choice of the statistic parameters to be printed on the report. Here the formats of the size distributions from which the statistic is calculated can be selected as well.

Printed Report	X
Sample In <u>f</u> o	Clear <u>A</u> ll
Size Graphs Select	Cancel
🔽 Statistics	Size Statistics
C Averaged Statistics	Mass Select
Compare Statistics	□ Number 🔽 🕷 🗆 /mL
🗖 Listing 🛛 Select	□ Volume □ % □ /mL
Interpolation Select	🗖 Surface Area 🗖 🖏 🗖 /ml
P <u>u</u> lses	<u>I</u> rend
Graph Select	Graph Select
Listing Select	Listing Select
Use Default Printer	Select
C Use Specified Printer:	

The kind of size distribution can be selected by checking the appropriate box.

Volume % Mass Number Volume/mL Number % Surface Area Number/mL Surface Area % Volume Surface Area/mL 

The Size Statistics dialog box appears when  $\mathcal{D}$ -Select inside the Size Section. Choose the parameters needed on the Printed Report and  $\mathcal{D}$ -OK.

Size Statistics				×
	> 🔽 <	× کا	Results to Print <u>M</u> ean Median	Set All OK Clear <u>A</u> ll Cancel
10       %       10         25       %       25         50       %       50         75       %       75         90       %       90         %           Label Percent	%         2           %         10           %         20           %         40           %         50           %         -           %         -           %         -           %         -           %         -           %         -           %         -           %         -           %         -           %         -           %         -           %         -	μm 2 μm μm 10 μm μm 20 μm μm 40 μm μm 50 μm μm	<ul> <li>□ D(p,q) p 0 q 0</li> <li>□ Mean/Median ratio</li> <li>□ Mode</li> <li>□ S.D.</li> <li>□ Variance</li> <li>□ Q.V.</li> <li>□ Skewness</li> <li>□ Kurtosis</li> <li>□ 95% Conf. Limits</li> <li>□ Specific Surface Area</li> </ul>	Clear <u>A</u> ll Cancel Bange ○ All ○ Between Cursors ○ From 0 µm to 0 µm Statistic <u>Lype</u> ○ Arithmetic ○ Geometric
			J □ d <u>1</u> 0 □ d <u>5</u> 0 □ d <u>9</u> 0	C As Accumulated


## 8.2.4 Pulses Section of the Printed Report

#### 1. Pulse Graph

If Pulse Data is saved as described in **Section 6.2.7**, the options to display and print this information can be selected now. Check the appropriate box for including the **Pulse Graph** in the **Printed Report** and select options by **O**→**Select**.

Printed Report	x
I Sample In <u>f</u> o I Short	Clear <u>A</u> II
Size	Cancel
✓ Statistics	Size Statistics
Averaged Statistics	Mass Select
Compare Statistics	🗖 Number 🔽 🕷 🗖 /mL
Listing Select	□ Volume □ % □ /mL
Interpolation Select	🗖 Surface Area 🗖 % 🗖 /mL
Pulses	<u>I</u> rend
Graph Select	🗖 Graph 🛛 Select
Listing Select	Listing Select
Printer	
Use Default Printer	Select
C Use Specified Printer:	

The **Pulse Graph** dialog box will appear. After selecting the options, *A*→**OK**.

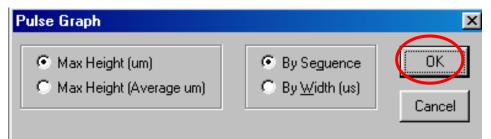

## 2. Pulse Listing

Check the appropriate box for including the **Pulse Listing** in the **Printed Report** and select options by  $\mathcal{D}$ -**Select**.

Printed Report	×
Sample In <u>f</u> o	Clear <u>A</u> ll
<u>S</u> ize	Cancel
Graphs Select	
V Statistics	Size Statistics
Averaged Statistics	Mass Select
Compare Statistics	🗖 Number 🔽 🎖 🗖 /mL
Listing Select	🗖 Volume 🗖 % 🗖 /mL
☐ InterpolationSelect	🗖 Surface Area 🗖 % 🗖 /mL
P <u>u</u> lses	_ <u>I</u> rend
Graph Select	🗖 Graph 🥂 Select
Listing Select	Listing Select
_ <u>P</u> rinter	
Use Default Printer	Select
C Use Specified Printer:	

The **Pulse Listing** dialog box will appear. After selecting the options,  $\mathcal{D} \rightarrow OK$ .

Pulse Listing	×
Columns Seguence	
Max Height (um)	Max Height (average)(um)
Max Height (V)	
□ <u>W</u> idth (us)	
Sequence Max Pulse Height	Max Pulse Height μm
Pulse Grouping	
<u>G</u> roup Size: 10	C All Pulses
C First Pulse in Group	Between Cursors
Sum Pulses in Grou	p C From 1 to 1


## 8.2.5 Trend Section of the Printed Report

#### 1. Trend Graph

Check the appropriate box for including the **Trend Graph** in the **Printed Report** and select options by  $\mathcal{A}$ ->Select.

Printed Report	x
Sample In <u>f</u> o	Clear <u>A</u> ll OK
Size	Cancel
✓ Statistics	Size Statistics
Averaged Statistics	Mass Select
Compare Statistics	T Number 🔽 % T /mL
Listing Select	□ Volume □ % □ /mL
☐ InterpolationSelect	☐ Surface Area ☐ % ☐ /mL
P <u>u</u> lses	Trend
Graph Select	Graph Select
Listing Select	Listing Select
Printer-	
O Use Default Printer	Select
C Use Specified Printer:	

The **Trend Graph** dialog box will appear. After selecting the options, *A*→**OK**.

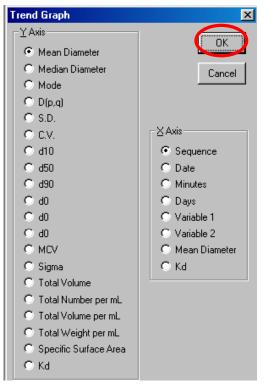

## 2. Trend Listing

Check the appropriate box for including the **Trend Listing** in the **Printed Report** and select options by <sup>-</sup>∂**-Select**.

Printed Report	X
Sample In <u>f</u> o	Clear <u>A</u> ll
Size Graphs Select	Cancel
Statistics	Size Statistics
Averaged Statistics	Mass Select
Compare Statistics	🗖 Number 🛛 🖉 🎖 🗖 /mL
Listing Select	□ Volume □ % □ /mL
☐ Interpolation Select	🗖 Surface Area 🗖 % 🗖 /mL
P <u>u</u> lses	Irend
Graph Select	Graph Select
Listing Select	Listing Select
_ <u>P</u> rinter	
Ose Default Printer	Select
C Use Specified Printer:	

The **Trend Listing** dialog box will appear. After selecting the options, *A*→**OK**.

Trend Listing		×
Columns (Select up to ☐ Seguence ☐ Eile Name ☑ Mean ☐ Median ☐ Mode ☐ D(p,q) ☐ S.D.	☐ Date ☐ Time ☐ Days ☑ Minutes ☐ MC⊻ ☐ Sigma ☐ Total Volume	Cancel
「 <u>C</u> .V. 「d <u>1</u> 0 「d <u>5</u> 0 「 <u>d<u>9</u>0 「<u>K</u>d</u>	<ul> <li>Total Number per mL</li> <li>Total Volume per mL</li> <li>Total Weight per mL</li> <li>Variable 1</li> <li>Variable 2</li> <li>Specific Surface Area</li> </ul>	☐ <u>U</u> se Trend Limits
Mean Minutes Diameter μm		

#### 8.2.6 Printer Section

The default printer can be used to print the report or another printer attached to the computer can be selected.

Printed Report	
I Sample In <u>f</u> o I Short	Clear <u>A</u> ll OK
_ <u>S</u> ize ✓ GraphsSelect	Cancel
✓ Statistics	Size Statistics
Averaged Statistics	Mass Select
Compare Statistics	🗖 Number 🔽 🕷 🗖 /mL
Listing Select	□ Volume □ % □ /mL
Interpolation Select	🗖 Surface Area 🗖 % 🗖 /mL
P <u>u</u> lses	_ <u>T</u> rend
Graph Select	🗖 Graph 🛛 Select
Listing Select	Listing Select
Dista	
Printer C Use Default Printer	Select
Use Specified Printer:	

## 8.2.7 Closing the Printing Report Dialog Box

At this moment all the selections for the Printed Report have been made.  $\textcircled{} \rightarrow OK$ .

Printed Report	<u>×</u>
I Sample In <u>f</u> o I Short	Clear <u>A</u> ll OK
Size	Cancel
✓ Statistics	Size Statistics
Averaged Statistics	Mass Select
Compare Statistics	🗖 Number 🔽 🕱 🗖 /mL
Listing Select	□ Volume □ % □ /mL
Interpolation Select	🗖 Surface Area 🗖 % 🗖 /mL
P <u>u</u> lses	<u>I</u> rend
Graph Select	Graph Select
Listing Select	Listing Select
<u>Printer</u>	
O Use Default Printer	Select
Use Specified Printer:	

## 8.3 <u>Customizing the Preferences</u>

This section describes how to set up the preference file. **Section 8.2** described how to select the options for some of the preferences, in this section they are explained in more detail.

Size Graph	Trend Graph
Size Listing	Trend Listing
Size Statistics	Pulse Graph
Size Interpolation	Pulse Listing

If you already made these selections, it is not necessary to modify them again.

#### 8.3.1 Size

To set up the Size Preferences, on the main menu  $\mathcal{D} \rightarrow Preferences$  and  $\mathcal{D} \rightarrow Size$ . A submenu showing the options for Size will appear.

📴 Multisizer 3 - Simulated		
<u>File Window R</u> un System <u>S</u> ettings	<u>P</u> references Sec <u>u</u> rity <u>H</u> elp	
	Frinted <u>R</u> eport	
Open Overlay Up Fonts Overlay	( <u>S</u> ize)	Size <u>G</u> raphs
	<u>T</u> rend	Size Listing
READY	<u>P</u> ulse I	Size <u>S</u> tatistics
	⊻iew Options	Size Interpolation
	Averaging and Trend	
	Number and Date Formats	
Sample Information	Data E <u>x</u> port	
Group ID: D54 Control2	Creek Oeviews	-
Sample ID:	Graph <u>O</u> ptions Datas Catura	
Comment: water	Page Set <u>u</u> p Diele Ferrie	
	Pic <u>k</u> Fonts	
Change	<u>W</u> indow Colors	
	Customize Tool <u>b</u> ar	_
Run Settings	Load Preferences	
	Edit Preference Info	
Control Mode: Volumetric 500 μL	Save Preferences As	
	Save Preferences As <u>D</u> efault	
Aperture: 100 μm	1	
Kd: 111	Language	
Current: 1600 µA		
Gain: 2		


#### 1. Size Graph


In the **Size** submenu  $\mathcal{O}$  **Size Graph**. The dialog box containing the options to display and print the **Size Graph** will appear. A separate graph will be printed for every box checked in the **Distribution Types** section. These graphs will be printed in the order they are selected, and the printing order will be displayed in the lower left-hand corner of the dialog box.

Size Graphs										×
Distribution Type: <u>D</u> ifferential: <u>C</u> umulative <: C <u>u</u> mulative >: Djff. + Cum. <: Diff. + Cum. >:	Nu	mber %/m I▼ □ □□□□	     [] 				Surl	íace A % Г Г	rea /mL C C	OK Cancel Clear <u>A</u> ll
Graph Printing O			-⊻Axis ⊂Lir ∙Lo ⊂As	s near	umula	ated	⊡ Y	- 7/d× D	er Curs liff. Graj	


#### **Types of Distribution**


For each kind of **Size Distribution**; (mass, number, number %, number/mL, volume, volume %, volume/mL, surface area, surface are % and surface area/mL), there are five different types of graph to choose from.

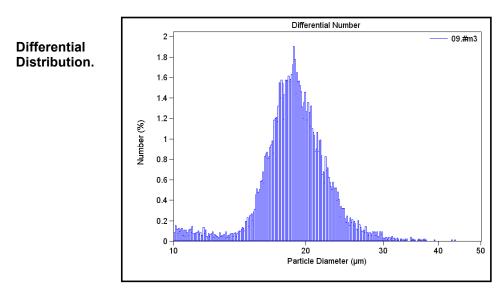

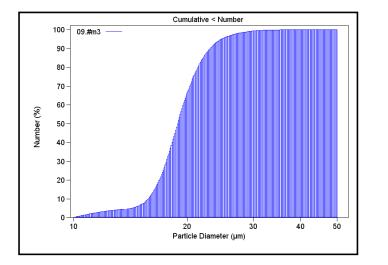


Cumulative <

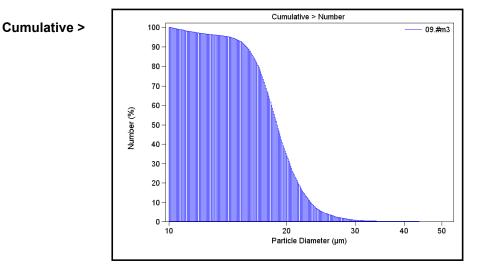

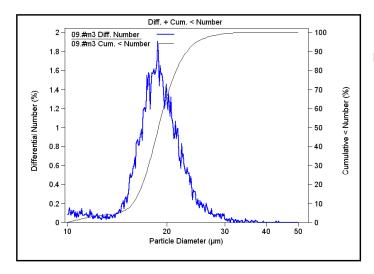


Differential + Cumulative <

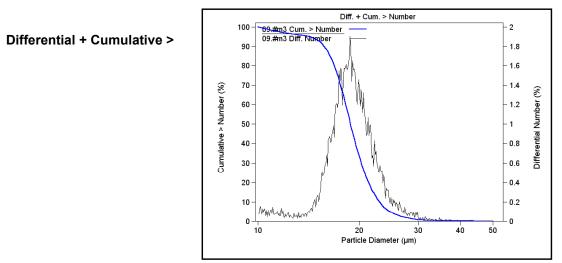


#### X Axis Scale

The X axis scale may be displayed in either linear or logarithmic. The logarithmic scale offers more detail for the lower side of the distribution. When a file acquired on a Multisizer IIe is open using the Multisizer 3 software, it is advisable to choose the option "as accumulated".

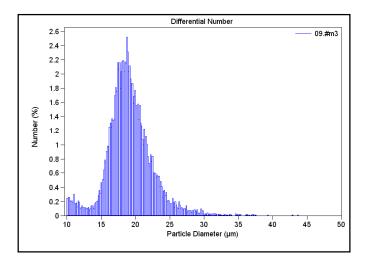

Linear X axis



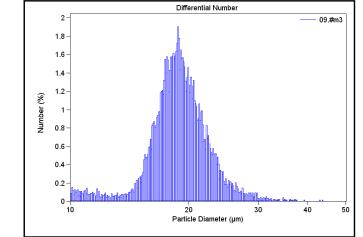

#### Other Options for the Size Graph

- □ Logarithmic Y axis
- □ **Multisizer Cursors.** This option shows where the cursors are placed when a file acquired on the Multisizer IIe is open using the Multisizer 3 Software.
- □ Y/dX Graph.
- □ Use Sample ID as Title. It places the sample ID entered in Step 6.5 as the graph title.

When done selecting the **Size Graph** options,  $\neg \ominus \rightarrow OK$  in the **Size Graph** dialog box.

#### 2. Size Listing

In the Size submenu, A→Size Listing, the dialog box containing the options to display and print the Size Listing will appear.

Size Listing							X
Columns (Select	: up to 7)						OK
Bin Number			Mass N	umber %/mL	Volume %/mL	Surface Area % /mL	
🔽 Lower Edge	💿 Diamete	r <u>D</u> iff.:					Cancel
Center	C Volume	<u>C</u> um. k:		ГГ	ГГГ		
🔲 Upper Edge	O Surface	Area ⊂C <u>u</u> m. >:					Clear <u>A</u> ll
	Bin Diameter (Lower) μm	Diff. Number %	Diff. Volume %				
Bin Grouping							
<u>G</u> roup Size:	1	🔿 All Bins					
C First Bin in (	àroup	🖲 <u>B</u> etween 0	Cursors				
Sum Bins in	Group	C From 1	to 1	]			

Selecting "Between Cursors" will display the Listing for only the channels within the range being used. When "All Bins" is selected, all the channels for the entire range of the aperture (2% to 60%), will be listed. "From" allows selection of the channel range for the Listing.

When done selecting the **Size Listing** options,  $\neg \ominus \rightarrow OK$  in the **Size Listing** dialog box.

#### 3. Size Statistics

In the Size submenu,  $\mathcal{D}$   $\rightarrow$  Size Statistics. The dialog box containing the options to display and print the Size Statistics will appear.

Selecting "Between Cursors" will calculate the Statistics for only the size data within the range selected with the cursors on the graph. When "All Bins" is selected, all the channels for the entire range of the aperture (2% to 60%), will be used in the calculations. "From" allows selection of the size range for the calculations.

When done selecting the Size Statistics options,  $\mathcal{A} \rightarrow OK$  in the Size Statistics dialog box.

#### 4. Size Interpolation

The Multisizer 3 software may be used to create sample data reports that will arrange data to interpolate in a set of user defined boundaries. This is similar to how particle data points are reported when different size sieves are used.

Two types of interpolation format may be used. In the first format, the boundaries represent size classes. In the second format, the particle data are contained in percentage classes. There are 3 choices: "Over", "Under", and "Differential". You can redefine both the number of boundaries and the value assigned to each boundary.

In the Size submenu,  $\mathcal{D}$  -> Size Interpolation. The dialog box containing the options to display and print the Size Interpolation will appear.

Size Interpolation		×
Interpolation <u>Iype</u> ○ Size to:         □ <u>N</u> umber       2 □ /mL         □ <u>V</u> olume       2 □ /mL         □ <u>S</u> urface Area       2 □ /mL         ⓒ to Size       - /mL         ○ <u>U</u> nder (<)	Interpolation Points         /mL         10.5         15.3         20.2         Delete         Bange         C         All         Image         C         From         Image         Image	OK Cancel

Selecting "Between Cursors" will calculate the Interpolation points for only the size data within the range selected with the cursors on the graph. When "All Bins" is selected, all the data for the entire range of the aperture (2% to 60%) will be used in the calculations. "From" allows selection of the size data range for the calculation of the Interpolation points.

When done selecting the Size Interpolation options,  $\mathcal{D} \rightarrow OK$  in the Size Interpolation dialog box.

#### 8.3.2 Trend

🔟 Multisizer 3 - Simulated - Operator mode				
<u>File W</u> indow <u>R</u> un System <u>S</u> etting	Preferences ec <u>u</u> rity <u>H</u> elp			
Open Overlay Up Fonts	Printed <u>S</u> eport			
READY	Irend Pulse	Trend <u>G</u> raph Trend Listing		
READT	View Options	<u>Custom Trend Points</u>		
Sample Information	<u>A</u> veraging and Trend Number and Date <u>F</u> ormats Data E <u>x</u> port			
Group ID: 22 Sample ID: Comment:	Graph <u>O</u> ptions Page Set <u>u</u> p Pick Fonts			
Change	Window Colors ► Customize Tool <u>b</u> ar			
SOM	Load Preferences Preference Info			
Control Mode: Volumetric 500 µL	Save Preferences As			
Aperture: 100 μm	Save Preferences As <u>D</u> efault			
Kd: 111	Language			
Current: 1600 µA				


#### 1. Trend Graph

## In the **Trend** submenu, <sup>A</sup>→**Trend Graph.**

The parameters in the larger section at the left side of the dialog box can be selected for the Y-axis. The parameters in the smaller section can be used on the Xaxis. When done selecting the **Trend Graph** options, →**OK** in the **Trend Graph** dialog box.

#### 2. Custom Points

Up to three points can be customized.

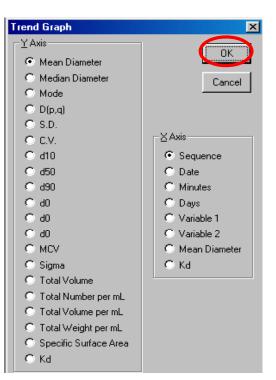

Mean           Mean           D(p,q)           D(p,q)	Date     Clear All       Time     Days       Days     Minutes       MC⊻     Sigma       Total Volume     Total Volume       Total Volume per mL     Total Volume per mL       Total Volume per mL     Variable 1       Variable 1     Variable 2       Specific Surface Area	Cancel	Trend Listing In the Trend submenu, ♪⊕→Trend Listing. When finished selecting the Trend Listing options, ♪⊕→OK in the Trend Listing dialog box.
Mean Minutes Diameter μm			

#### 8.3.3 Pulses

If Pulse Data is saved as described in <u>Section 6.2.7</u>, the options to display and print this information may be selected at this time. On the main menu  $\textcircled{} \rightarrow \mathbf{Preferences}$  and  $\textcircled{} \rightarrow \mathbf{Pulse}$ . A submenu showing the options for **Pulse** will appear.

🖪 Multisizer 3 - Simulated		
<u>File W</u> indow <u>R</u> un System <u>S</u> ettings	Preferences Security Help	
Open Overlay Up Fonts Overlay	Printed <u>R</u> eport Size ▶	Edit Graph Listing Zoom
READY	Pulse	Pulse <u>G</u> raph Pulse <u>L</u> isting
Sample Information Group ID: D54 Control2 Sample ID: Comment: water Change	Averaging and Trend Number and Date <u>F</u> ormats Data E <u>x</u> port Graph <u>O</u> ptions Page Set <u>up</u> Pic <u>k</u> Fonts <u>W</u> indow Colors Customize Tool <u>b</u> ar	
Run Settings Control Mode: Volumetric 500 μL Aperture: 100 μm Kd: 111 Current: 1600 μA	Load Preferences Edit Preference Info Save Preferences As Save Preferences As <u>D</u> efault Language	

1. Pulse Graph

In the **Pulse** submenu, **A**→**Pulse Graph**.

Pulse Graph		×
<ul> <li>Max Height (um)</li> <li>Max Height (Average um)</li> </ul>	<ul> <li>G By Seguence</li> <li>C By <u>W</u>idth (us)</li> </ul>	OK Cancel

Pulse Maximum High can be used for the Y-axis and Pulse Sequence or Pulse Width on the X-axis. When finished selecting the **Pulse Graph** options,  $\mathcal{O} \rightarrow OK$  in the **Pulse Graph** dialog box.

#### 2. Pulse Listing


In the **Pulse** submenu, **A**→**Pulse** Listing.

Pulse Listing	×
Columns Seguence Max Height (um) Max Height (V) Width (us)	Clear <u>A</u> ll OK Max Height (average)(um) Cancel
Sequence Max Pulse Height	Max Pulse Height μm
Pulse Grouping	
<u>G</u> roup Size: 10	C All Pulses
C First Pulse in Group	Between Cursors
Sum Pulses in Group	D C From 1 to 1

When finished selecting the **Pulse Listing** options,  $\mathcal{D} \rightarrow OK$  in the **Pulse Listing** dialog box.

#### 8.3.4 Other Options for the Preference File

On the main menu, → Preferences and → View Options. A dialog box will appear.

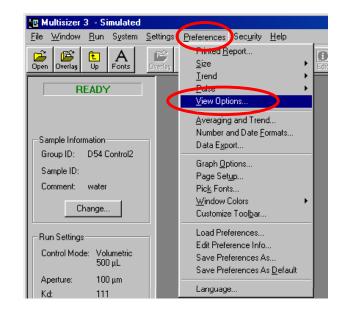

View Options	×
– Size ⊻ Axis	ОК
Diameter	
C Volume	Cancel
🔿 Surface Area	
C As Accumulated	
_ <u>M</u> easuring	
Particles	• μm <sup>3</sup>
🔿 Cells	C fL
Specific Surface <u>A</u> rea © cm²/mL © m²/mL	Units
 Size Results Represer	ntation
🔲 🗍 ISO 9276-1 (prelim	ninary)
Background Subtract	
Allow <u>n</u> egative res	ults
Multisizer 3 Pulse Data	
Show at most 52500	0 Pulses
Z2 Data	
Suppress last char	nnel

#### 1. X-Axis Section

The units for the X-axis can be expressed as diameter, volume or surface area.

**Diameter:** this is the most common use for the X-axis for industrial applications.

**Volume:** this is widely used in biological applications **Area:** this is used for specialized studies.

**As Accumulated:** this option is used when a file acquired on the Multisizer IIe is opened using the Multisizer 3 software.

#### 2. Measuring Section

Particles or Cells can be selected. This selection is used only for changing the X-axis units for displaying the data and /or printing the report. Where it is critical to select Cells is in <u>Step 6.2.4</u>.

#### 3. Volume Units Section

When Diameter is used for the X-axis, units can be  $\mu^3$  or fL. In cell biology, fL is more commonly used

#### 4. Specific Surface Area Units

When Surface Area is used for the X-axis, units can be selected as  $\mbox{cm}^2/\mbox{mL}$  or  $\mbox{m}^2/\mbox{mL}.$ 

#### 5. Multisizer 3 Pulse Data

In this box the number of pulses to be saved can be selected. Up to 525,000 pulses may be saved.

#### 6. Z2 Data


If a file acquired on the Z2 is opened using the Multisizer 3 software, the last channel can be suppressed by checking this box.


#### 7. ISO 9276 (Preliminary)

The choice of having all units represented in accordance with ISO 9276-1 (preliminary) is also included as an option for both graphs and listings.

#### 8.3.5 Averaging and Trending

The options for the weighting and type of distribution for the average and also the setting for the limits averages and trends are located under **Averaging and Trend**.

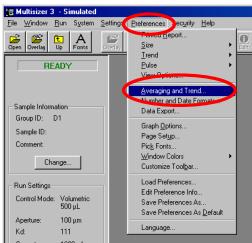

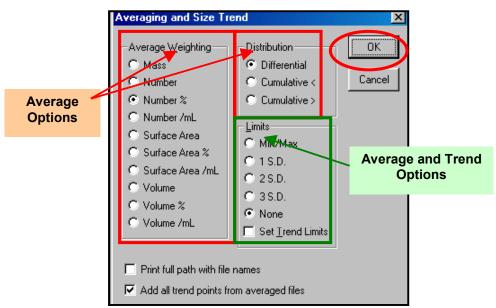

On the main menu  $\mathcal{T} \rightarrow \mathsf{Preferences}$  and  $\mathcal{T} \rightarrow \mathsf{Averaging}$  and Size Trend. The following dialog box will appear.

The two boxes at the bottom of the dialog allow file names to be printed with their full path, and another option to add to the report all data points from the averaged or trend files.

Once the options have been selected,  $\mathcal{P} \rightarrow OK$ .

#### 8.3.6 Number and Date Formats

The **Number and Date Format** dialog specifies what number format will be used for data files that are being exported to other applications. This function also determines how numbers are displayed in tabular data listings, and determines whether the Multisizer software represents the decimal point as a space, a comma or a period. The date can also be formatted under this **Preferenc**e option.


On the main menu,  $\mathcal{P} \rightarrow \mathsf{Preferences}$  and  $\mathcal{P} \rightarrow \mathsf{Number}$  and Date Format

🔟 Multisizer 3 - Simulated	
<u>File Window Run System Settings</u>	Preferences Security Help
Copen Overlag Up Fonts Corrier	Printed Beport       Size         Size       Image: Size size         Irend       Image: Size size size size size size size size s
Sample Information Group ID: D1 Sample ID: Comment: Change Run Settings Control Mode: Volumetric 500 µL Aperture: 100 µm Kd: 111 Current: 1600 µA Gain: 2 Runs: 1	Number and Date Export       Data Export         Graph Options       Page Setup         Pick Fonts       Window Colors         Window Colors       Cancel         Customize Toolbar       Cancel         Load Preferences       Call 23456.78         Edit Preferences       Call 23456.78         Save Preferences As       Cancel         Save Preferences As       Call 23456.78         Language       No Rounding

Once the options have been selected,  $\neg \ominus \rightarrow OK$ .

#### 8.3.7 Data Export


Options in the **Data Export** dialog allow the user to specify which subsections of the runfile (e.g., the file header, size data, pulse data, etc.) will be exported. An extension, such as .xls, may be entered so that another application will recognize the exported data. Formatting options are also included in the **Data Export** dialog. On the main menu,  $\textcircled{O} \rightarrow Preferences$  and  $\textcircled{O} \rightarrow Data Export$ .

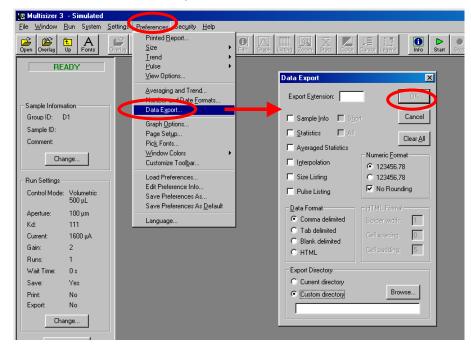

Once the options have been selected,  $\mathcal{D} \rightarrow OK$ .

#### 8.3.8 Graph Options

The **Graph Options** dialog box allows customizing the appearance of the graphs. In this dialog, separate color, line or dot styles may be specified for each data file, and the dialog also includes options for grid lines and legend position. The colors and styles for each file may be defined differently depending on whether the choices refer to a printed document or an on-screen display. On the main menu,  $- P \rightarrow Preferences$  and  $- P \rightarrow P P$ 

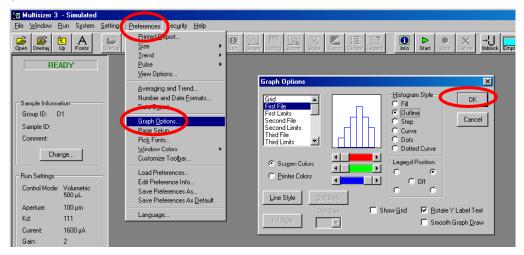

Once the options have been selected,  $\mathcal{D} \rightarrow \mathbf{OK}$ .

#### 8.3.9 Page Setup

Select this menu item to control printer functions, including printer set-up, margins, titles and page orientation of printed output. **Page Setup** also allows inserting a company logo as a page header by entering the file name of a **custom logo** in bitmap format using the **Custom Bitmap** option. This can be created using Microsoft Paintbrush to make a black and white or 16-color bitmap up to 600 pixels wide by 200 pixels high.

The full path to the bitmap must be entered, e.g. ("c:\windows\ls\mylogo.bmp",). This bitmap will show up at the top left of every page printed, replacing the Beckman Coulter logo. On the main menu,  $\textcircled{} \rightarrow \mathbf{Preferences}$  and  $\textcircled{} \rightarrow \mathbf{Page Setup.}$ 

📴 Multisizer 3 - Simulated		
<u>File W</u> indow <u>R</u> un System <u>S</u> etting:	Preferences ec <u>u</u> rity <u>H</u> elp	
Open Overlay Up Fonts Overlay	Printed Deport Size Trend	Eller Graph Listing Zoom Stats Color Largerd Legend
READY	<u>P</u> ulse <u>V</u> iew Options	Page Setup
Sample Information Group ID: D1 Sample ID: Comment: Change Run Settings Control Mode: Volumetric 500 µL Aperture: 100 µm Kd: 111 Current: 1600 µA Gain: 2 Runs: 1	Averaging and Trend Number and Date Eormats Data Egoort Page Setup Pick Fonts Window Colors Customize Toolbar Load Preferences Edit Preferences As Save Preferences As Save Preferences As Save Preferences As	Page Setup         Include Qustom Title         Print Graphs using Screen Colors         Landscape Qrientation         Include Page Number         Include Page Number         Include Date         Margin Settings (inches)         Include Date         Top:         Left:         O         Right:         OK         Printer Setup         Cancel         OK         Fill Page         Sixth Page         Fill Page

When the page setup has been selected,  $\mathcal{D} \rightarrow \mathbf{OK}$ .

#### 8.3.10 Pick Fonts

Use **Pick Fonts** to make changes in font type and font size. Separate settings are available for graphs, listings, statistics, and headers. The program allows previewing font size and font samples before making the choices. On the main menu  $\bigcirc \rightarrow$  **Preferences**, then  $\bigcirc \rightarrow$  **Pick Fonts**.

📴 Multisizer 3 - Simulated	
<u>File W</u> indow <u>R</u> un System <u>Settings</u> <u>P</u> referen	nces ecurity Help
Open Overlay Up Fonts Overlay Size	Lan arean ability additi orace additi
READY Pulse	e → v Dpions
Avera	raging and Trend Pick Fonts
	nber and Date Eormats a Export <u>G</u> raphs Arial, 12 OK
Sample ID: Grap	e Serban E Serban Listings Arial, 12
Change	dow P ofs
Lusto	tomize Toolbar Arial, 12
Run Settings Edit F Control Mode: Volumetric Save	d Preferences Preferences Info e Preferences As e Preferences As Default
Aperture: 100 μm	
Kd: 111 Lang	guage
Current: 1600 µA	

After making changes, ↔ → Apply to implement changes, then ↔ → OK.


#### 8.3.11 Windows Colors

Choices here will control the default colors for the text, warning text and the window background.  $\textcircled{} \rightarrow Preferences$ , and  $\textcircled{} \rightarrow Windows$  Colors. After selecting colors for each option,  $\textcircled{} \rightarrow OK$ .

✓ 🙂 These cho	pices will not affect any other Windows applications.
T Multisizer 3 - Simulated	
File     Window     Bun     System     Settin       Image: Comparison of the processing of the proc	Bised Beport     Size       Size     Fair       Irend       Pulse       Vew Options
Sample Information Group ID: D1 Sample ID: Comment: Change SOM Control Mode: Volumetric 500 µL Aperture: 100 µm	Averaging and Trend         Number and Date Eornats         Data Egont         Graph Options         Page Setup         Window Colors         Window Colors         Usad Preferences         Edit Preferences As         Save Preferences As         Save Preferences As         Save Preferences As         Save Preferences As
Kd: 111 Current: 1600 μA Gain: 2 Runs: 3 Wat Time: 60 s Save: Yes Print: No Export: No Chonge	Custom colors:

#### 8.3.12 Customize Toolbar


This option allows you to select wether to show or to hide the toolbar and the way the toolbar buttons will be displayed. On the main menu  $\neg \ominus \rightarrow Preferences$ , and,  $\neg \ominus \rightarrow Customize Toolbar$ .

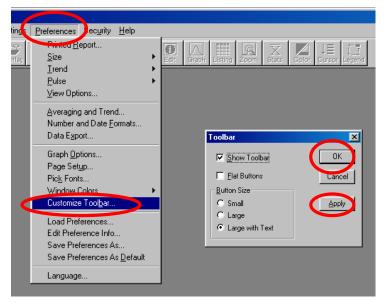

After selecting options,  $\neg \ominus \rightarrow Apply$  to implement changes, then  $\neg \ominus \rightarrow OK$ .

#### 8.3.13 Selecting the Language

To select the language,  $\mathcal{D} \rightarrow \mathbf{Preferences}$  then  $\mathcal{D} \rightarrow \mathbf{Language}$ .

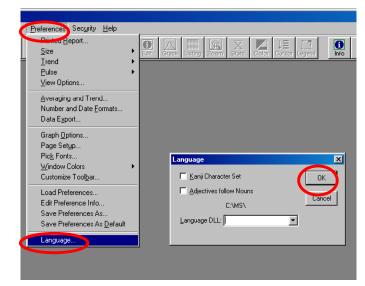


The available languages will appear in the drop down menu of the language dialog box. After selecting the language,  $\mathcal{A} \rightarrow OK$ .


#### 8.3.14 Saving Preferences

At this moment all the options for the Preferences have been selected.

All the options for the **Preferences** can be saved for future use. To save the **Preferences**,  $\mathcal{D}$ ->Settings on the Main Menu bar and then  $\mathcal{D}$ ->Save Preferences As.

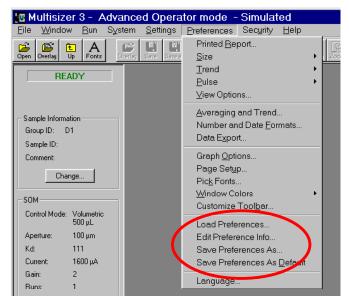

After assigning a name to identify it, the **Preferences** file can be saved in the sub-folder **Preferences** located in C:\MS\SOP. Any number of **Preferences** files can be saved; they can be used at a later time as the options to view the results and print the analysis report or for creating a **Standard Operating Procedure** (see <u>Section 9</u>).

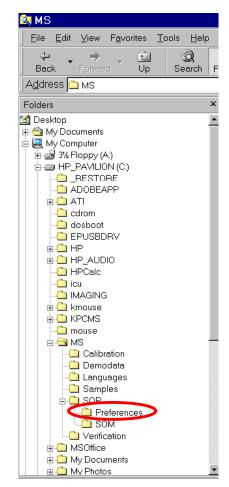

To use a Preferences file previously saved, ↔ Preferences on the Main Menu bar, then ↔ Load Preferences. To change a Preferences file, ↔ Horder Preferences Info.

The **Preferences** options can also be saved as the **Default Preferences**. These default options will appear every time the Multisizer 3 program is started. To save **Preferences** as default,  $\neg \partial \rightarrow Preferences$  on the Main Menu bar, then  $\neg \partial \rightarrow Save$  **Preferences As Default**.

#### Back to TOC

# SECTION 9 CREATING A STANDARD OPERATING PROCEDURE (SOP)

#### 9.1 Introduction

This section describes how to create and use **Standard Operating Procedures (SOP)**. This feature of the software allows you to combine **Preferences** files and **Standard Operating Methods (SOM)** files into one single file. This new file will include all the necessary settings and options described in any user defined **SOP**. Creating and saving SOP's is specially useful when more than one operator is using the Multisizer 3 or when the instrument is used for more than one type of application. The use of SOP's also ensures the consistency of the analyses by preventing an unauthorized operator to change any of the settings or options included in the **SOP**, (see <u>Section 10</u>).

#### 9.2 Creating and Saving an SOP

Before creating an SOP, it is necessary to create and save the **Standard Operating Method (SOM)** file and the **Preferences** file (see <u>Section 6.2 to 6.4</u> and <u>Section 8</u>).

🔟 Multisizer 3 - Simulated	To create the SOP file, on the
<u>F</u> ile <u>W</u> indow <u>R</u> un System S	ettings Preferences Security Help main menu 🕀 -> Settings and in
Open Overlay Up Fonts Overlag	Load an SOP the drop down list select Create
	an SOP.
READY	Create an SOP
	Change SOM
	Load an SOM
Sample Information	Save an SOM
Group ID: D1	Enter Sample Info
Sample ID:	Load Sample Info
Comment:	Convert Pulses to Size Settings
Change	Sizing Threshold
	Set Current and <u>G</u> ain
- SOM	Select <u>Electrolyte and Dispersant</u>
Control Mode: Volumetric	Control Sample Options
500 µL	Eile Name Generation
Aperture: 100 μm	Load Background Run
Kd: 111	Remove Background
Current: 1600 µA	Advanced
Gain: 2 Buns: 1	Diagnostics
Wait Time: 0 s	Status Panel
Save: Yes	
	Create a Standard Operating Procedure
	SOM File
The dialog box to select	SUM File Save
the necessary files for	
creating the SOP will	Cancel
appear.	
••	✓ Do not allow operator to change SOM settings     Clear <u>A</u> II
	Select
	Do not allow operator to change preferences
	iver both of allow operator to change presences
	Comment:

## 9.2.1 Selecting the SOM file.

In the **SOM** file section of the **Create a Standard Operating Procedure** dialog box ↔→**Select**.

Create a Standard Operating Procedure	×
SOM File	Sa <u>v</u> e
✓ Do not allow operator to change SOM settings Preference File Select	Cancel Clear <u>A</u> ll
Do not allow operator to change preferences	
<u>C</u> omment:	

A dialog box will display the folders under C:\MS. Select the f and ∽⊕**→Op** 

ie folder <b>SOP</b> Open.	Look jn: MS Calibration Demodata Languages Samples SOP Verification File name: Files of type: SOM (*.som) Recent Directories: C:MS	
Load a Standard Ope Look jn: 🔄 SOP Preferences SOM	erating Method (SOM) 🛛 ? 🗙	A dialog box will display the folders under C:\MS\SOP. Select the folder <b>SOM</b> and ℃ <b>→Open</b> .
File name: Files of type: SOM (*.som) Recent Directories: C:\MS\SOP	Cancel	

Load a Standard Operating Method (SOM)


? ×

After selecting the folder **SOM**,  $\mathcal{D} \rightarrow \mathbf{Open}$ . All the previously saved **SOM** files will be shown (see <u>Section 6.4</u>). Select the necessary file to be included in the **SOP** and  $\mathcal{D} \rightarrow \mathbf{Open}$ .

Load a Standard Operating Method (SOM) 💦 🛛 🔀	
Look in: 🔁 SOM 💽 🖛 🖻 📸 🎹 🗸	
demo.som	
File name:	
Files of type: SOM (*.som)	
Recent Directories: C:\MS\SOP\SOM	
<u>G</u> et Info	
Create a Standard Operating Procedure	×
SOM File demo.som C:\MS\SOP\SOM\demo.som	Sa <u>v</u> e Cancel
Do not allow operator to change SOM settings	Clear <u>A</u> ll
Preference File Select	
Comment:	

#### 9.2.2 Selecting the Preferences File

In the **Preference** file section of **Create a Standard Operating Procedure** dialog box *√*<sup>0</sup>→**Select**.

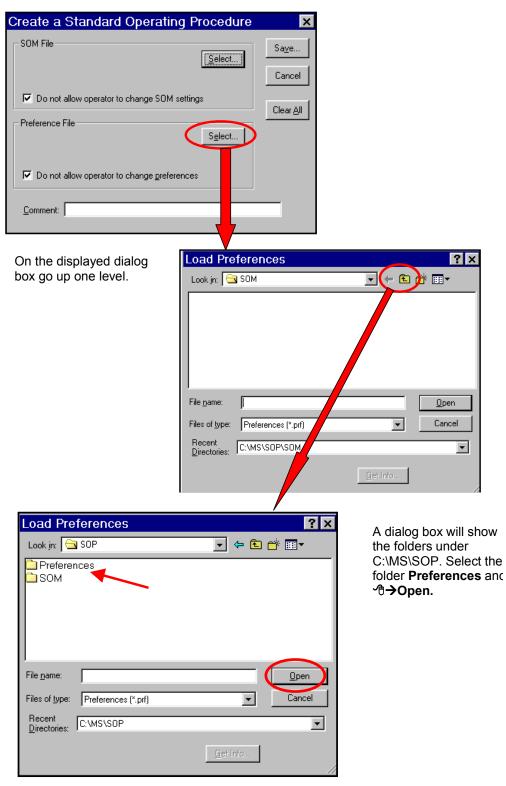

Cr

After selecting the folder **Preferences**, **\*∂→Open**. All the previously saved **Preferences** files will be shown (see <u>Section 8.3.14</u>). Select the necessary file to be included in the **SOP** and **\*∂→Open**.

	Load Prefe	erences			? ×
	Look jn: 🔂 P	references	•	🗢 🗈 🔿	<u>.</u>
	PDEMO				
	File <u>n</u> ame:	DEMO			<u>O</u> pen
	Files of <u>type</u> :	Preferences (*.prf)			Cancel
	Recent Directories:	:\MS\SOP\Preferences	3		•
	<u>D</u> irectories.		<u>G</u> et Ir	nfo	
eate a Standar	d Oneratin	na Procedure	×		
SOM File demo.som C:\MS\SOP\SOM\demo		Select	Sa <u>v</u> e Cancel		
🔽 Do not allow operator	to change SOM	settings			
Preference File DEMO.PRF C:\MS\SOP\Preferences IV Do not allow operator	\DEMO.PRF	S <u>e</u> lect	Clear <u>A</u> ll		
Comment: Demo SOP			_		

## 9.2.3 Saving the SOP

At this moment the two files needed to create an **SOP** have been selected. A comment can be added and it will be shown in the Info box for the **SOP** (see <u>Section 9.2</u>). To save the **SOP** ö→Save.


Create a Standard Operating Procedure	×
SOM File demo.som <u>Select</u> C:\MS\SOP\SOM\demo.som ✓ Do not allow operator to change SOM settings Preference File DEMO.PRF <u>Select</u> C:\MS\SOP\Preferences\DEMO.PRF	Cancel
Comment: Demo SOP	

SOM
 SOM
 SOffication
 MSOffice
 My Documents
 My Photos

F


×

٠

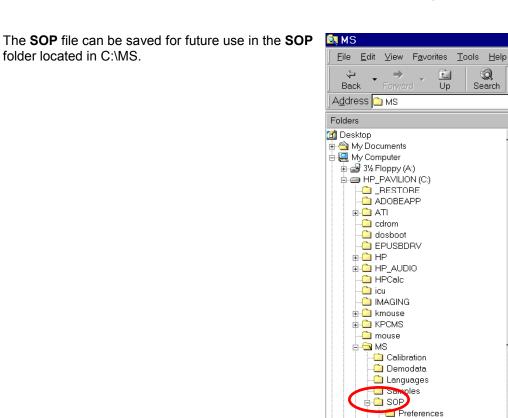

#### 9.3 Using an SOP

Any number of SOP files can be created and saved for future use (see <u>Section 9.1</u>). To use a previously saved SOP file,  $\textcircled{} \rightarrow$ **Settings** on the main menu and select. Load an **SOP.** 

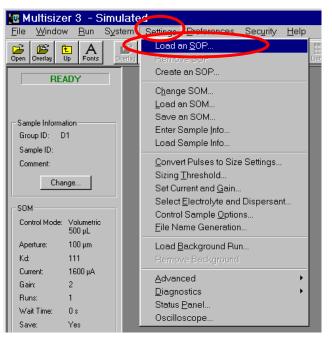

A dialog box will show all the **SOP** files previously saved. Select the required **SOP**. The **Get Info** button shows a display of all the information about the selected SOP file. To use the SOP file for analyzing samples in the Multisizer 3,  $\Im \rightarrow OK$ .

Load a S	tandard Ope	rating Proc	edure (SO)	P) ?×
Look jn: 🧲	SOP		- 🗧 🖶 💣	
Preferent				
, File <u>n</u> ame:	Demo		<b>(</b>	<u>O</u> pen
Files of <u>type</u> :	SOP (*.sop)			Cancel
Recent <u>D</u> irectories:	C:\MS\SOP			•
			<u>G</u> et Info	

#### 9.4 Removing the Active SOP

If during a working session an **SOP** file is in use and a different file is needed, just follow instructions in <u>Section 9.2</u> and load a new **SOP** file. This action will remove the previous file from use. When there is no need for using an SOP file, the current file in use can be removed..  $\textcircled{\rightarrow}$  **Settings** on the main menu and select **Remove SOP**.

🖪 Multisizer 3 - Simulated					
<u>F</u> ile <u>W</u> indow <u>R</u> un System	Settings Preferences Sec <u>u</u> rity <u>H</u> elp				
Open Overlay Up Fonts Overlay	Remove SOP (Demo.sop)				
Open         Overlag         Up         Fonts         Overlag           READY         Sample Information         Group ID:         D1         Sample ID:         Comment:         Comment:         Change         Change         Comment:         Comment: <td>Cleate an SOP Change SOM Load an SOM Save an SOM Enter Sample Info Load Sample Info <u>C</u>onvert Pulses to Size Settings Sizing <u>T</u>hreshold Set Current and <u>G</u>ain</td>	Cleate an SOP Change SOM Load an SOM Save an SOM Enter Sample Info Load Sample Info <u>C</u> onvert Pulses to Size Settings Sizing <u>T</u> hreshold Set Current and <u>G</u> ain				
SOM Control Mode: Time 60 s	Select <u>E</u> lectrolyte and Dispersant Control Sample <u>O</u> ptions <u>F</u> ile Name Generation				
Aperture: 100 μm Kd: 111	Load <u>Background Run</u> Remove Background				
Current: 1600 µA Gain: 2 Runs: 1 Wait Time: 0 s Save: Yes	Advanced  Diagnostics Status Panel Oscilloscope				
Print: No					

#### Back to TOC

# SECTION 10 REGULATORY COMPLIANCE

#### 10.1 <u>21 CFR Part 11</u>

The Electronic Records and Electronic Signatures Rule (21 CFR Part 11) was established by the FDA to define the requirements for submitting documentation in electronic form and the criteria for approved electronic signatures. This rule, which has been in effect since August 20, 1997, does not stand in isolation; it defines the standards by which an organization can use electronic records to meet its record-keeping requirements. Organizations that choose to use electronic records must comply with 21 CFR Part 11. It is intended to improve an organization's quality control while preserving the FDA's charter to protect the public. Since analytical instrument systems such as the Multisizer<sup>™</sup> 3 generate electronic records, these systems must comply with the Electronic Records Rule.

Here are described the relevant portions of the 21 CFR Part 11 regulations and their implementation using the Multisizer<sup>™</sup> 3 control software explained. It is important to realize that Implementation and compliance of the rule remains the responsibility of the organization or entity creating and signing the electronic records in question. Proper procedures and practices, such as GLP and GMP, are as much part of overall compliance with these regulations as are the features of the Multisizer<sup>™</sup> 3 control software.

#### 10.2 <u>Electronic records</u>

Section 11.3 subpart A of 21 CFR Part 11 deals with the definitions assigned to electronic records, 'electronic record means any combination of text, graphics, data, audio, pictorial, or other information representation in digital form that is created, modified, maintained, archived, retrieved or distributed by a computer system'. In reality this refers to any digital computer file submitted to the agency, or any information not submitted but which is necessary to be maintained. Public docket No 92S-0251 of the Federal Register (Vol.62, No 54) identifies the types of documents acceptable for submission in electronic form and where such submissions may be made.

#### 10.3 FDA Requirements

In the general comments section of the ruling the following is stated: 'The agency emphasizes that these regulations do not require, but rather permit, the use of electronic records and signatures'. In the introduction to the final ruling the following statement is made The use of electronic records as well as their submissions to FDA is voluntary'.

If electronic submissions are made, Section 11.2 subpart A comes into play: ' persons may us electronic records in lieu of paper records or electronic signatures in lieu of traditional signatures....provided that: (1) The requirements of this part are met; and (2) The document or parts of a document to be submitted have been identified in public docket No. 92S-0251'.

The Multisizer<sup>™</sup> 3 control software have been designed to allow users to comply to the electronic records and signatures rule. Any organization deciding to employ electronic signatures must declare to the FDA their intention to do so.

#### 10.4 Implementation of electronic records and electronic signatures

Section 11.3 Subpart A describes two classes of systems; "closed systems" and "open systems". A closed system is one "in which system access is controlled by persons who are responsible for the content of electronic records". In other words, the people and organization responsible for creating and maintaining the information on the system are also responsible for operating and administering the system. In contrast, an open system

is one "in which system access is not controlled by persons who are responsible for the content of electronic records".

A typical Multisizer<sup>™</sup> 3 installation will need to have a procedure designed to ensure proper operation, maintenance and administration for system security and data integrity. Anyone who interacts with the system, from administrators to users, must abide by these procedures. Therefore the ultimate responsibility is with the organization generating electronic records and signatures. The Multisizer<sup>™</sup> 3 software is a component, albeit a vital one, of the overall process.

#### 10.5 Controls for electronic records

Subpart B, Section 11.10 describes the controls to be applied to a "closed system". Section 11.30 describes the controls for an "open system", which include "those identified in § 11.10, as appropriate, and additional measures such as document encryption and use of appropriate digital signature standards". Since a typical Mulitsizer™ 3 system can be regarded as a closed system, additional controls for open systems will not be discussed in this document. The primary thrust of these controls is "to ensure the authenticity, integrity, and, when appropriate, the confidentiality of electronic records, and to ensure that the signer cannot readily repudiate the signed record as not genuine". In other words, to protect the data and to make it difficult for someone to say that this is not their "signature". Many of the controls described in Section 11.10 refer to written procedures (SOP's) required of an organization by the agency, for the purpose of data storage and retrieval, access control, training, accountability, documentation, record keeping, and change control. The other controls are addressed either by the Multisizer™ 3 software itself, or in combination with end-user procedures.

Of the other controls, perhaps the foremost is described in Section 11.10 Paragraph (a): "Validation of systems to ensure accuracy, reliability, consistent intended performance, and the ability to discern invalid or altered records." It is the complete and overall validation of the system, as developed by the organization, which ensures the integrity of the system and the data within. It is to this end that the features of the Multisizer™ 3 software comply with the specifications of these regulations.

#### 10.6 Establishing an electronic record control with the Multisizer<sup>™</sup> 3

The Multisizer  $^{\text{TM}}$  3 software employs a system of usernames and passwords, consistent with the specifications of Subpart C, Section 11.300, "to ensure that only authorized individuals can use the system, electronically sign a record, access the operation or computer system input or output device, alter a record, or perform the operation at hand".

#### 10.6.1 Setting up security

If you selected the security option as described in <u>Section 3.1</u>, a dialog box appear requiring you to enter a user name and password. If the security option was not selected in **Section 3.1**, it can be selected at this time. On the main menu  $\bigcirc \bigcirc \bigcirc$  Security Set Up.

ĺ.										
I	U M	ultisizer	3 - Si	mulated	- Opera	tor Mode				
	<u>F</u> ile	$\underline{W}{indow}$	<u>B</u> un	System	<u>S</u> ettings	Preference	s Sec <u>u</u> rity	<u>H</u> elp		
	Ē	1	Ê	A	F			rity Set-up		Listing Z
	Open	Overlay	Up	Fonts	Overlag	Save Savea	<sup>ड</sup> Chan	ge <u>U</u> ser Mode	Graph	Listing Z
		R	EADY	,					_	
		mple Infor								
	G	roup ID:	22							
	Sa	ample ID:								
	C C	omment:								

The security level option dialog box is displayed.


Security Set-up	×
<ul> <li>Select security level</li> <li>No security</li> <li>Low security</li> <li>Medium security</li> <li>High security</li> <li>21 CFR Part 11</li> </ul>	OK Cancel

After selecting the required option  $\mathcal{T} \rightarrow OK$ . The **Add User Account** dialog box will appear.

#### 10.6.2 Adding a user account

Add an Administrator account when prompted. The administrator can then add and configure the level of access for all other users, when finished,  $\mathcal{T} \rightarrow OK$ . If more user accounts are needed, they can be added at this moment or at a later time by  $\mathcal{T} \rightarrow Administrator$  on the main menu and then  $\mathcal{T} \rightarrow Add$  a User Account.


Add a l	Iser Account		X
<u>E</u> ull I	Name: Jason Anderson		ОК
Log-in I	Name: Janderson		
	ID: JA	(optional)	Cancel
<u>P</u> as:	word: ****	(4 to 16 Characters)	Clear
	Type: Administrator	<b>•</b>	
	🗖 Disabled		

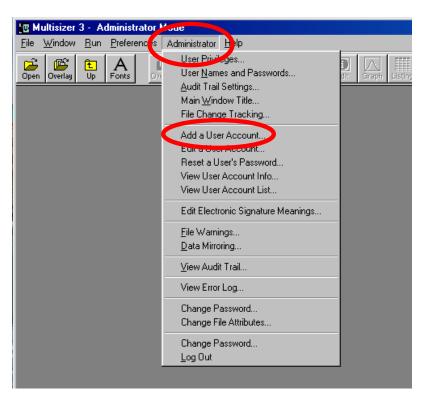

In the Administrator drop down menu select Add a User Account. The Add a User Account dialog box is displayed. Follow the instructions at the beginning of <u>Section</u> 10.7.2

#### 10.6.3 Selecting user names and passwords

In the Administrator drop down menu select User Names and Passwords.

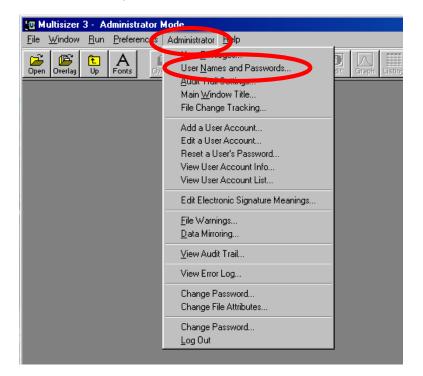

The entire Login process is controlled via the user names and passwords dialog. This ensures that the system login is controlled via time dependant metrics, such as password expiry and dialog inactivity, and has the required password lengths. Select the options in the dialog box and  $\mathcal{O} \rightarrow OK$ .

**√**☺)

Note that the dialog has a button marked 21 CFR 11. When activated this ensures all options are set so that the settings meet with the rules as applicable to 21 CFR part 11.

User Names and Passwords
User names
Log-in name must be at least 4 characters long (>1)
Eull name must be at least 3 characters long (>1)
User passwords
Must be at least 6 characters long (>1)
Will expire every 60 days (>1)
Automatically Log Out after 30 minutes of inactivity (>4)
Close <u>S</u> ignature dialog after 60 seconds of inactivity (>19)
Close Log-in dialog after 60 seconds of inactivity (>19)
Close Log-in dialog after 5 consecutive failed Log-in attempts (>1)
Disable account after 15 consecutive failed Log-in attempts (>1)


#### 10.6.4 File History and Audit Trail

The Multisizer™ 3 software also performs data input and "operational checks", as specified in Subpart B, Section 11.10, "to determine, as appropriate, the validity of the source of data input or operational instruction", and "to enforce permitted sequencing of steps and events". These two features ensure that, as much as possible, valid data are being entered into the system, and all required steps have been completed to perform the task at hand.

The purpose of all such data checking and validation is described in Section 11.10, Paragraph (b): "The ability to generate accurate and complete copies of records in both human readable and electronic form suitable for inspection, review, and copying by the agency". Consequently, strict procedures can be enforced within the Multisizer™ 3 software system to record all changes that are made to data generated from within Multisizer™ 3 software, as defined in Section 11.10, Paragraph (e).

Any file created by the Multisizer<sup>™</sup> 3 software can have 'File history' or auditing enabled. Once auditing has been enabled for a file, it cannot be disabled, nor can it be bypassed. Under these conditions, all changes made to a file are automatically recorded. These changes consist of "computer-generated, time-stamped audit trails to independently record the date and time of operator entries and actions that create, modify, or delete electronic records". When a change to a file is detected, the Multisizer<sup>™</sup> 3 software automatically records the identity of the user making the change, the date and timestamp of the change, the parameter that has changed, the old value and the new value. The user is also required to 're-sign' the record electronically and enter a reason for the change, from a pre-defined list or as free text. This audit trail is then stored as **File History** inside the file itself, such that "record changes shall not obscure previously recorded information", and in a "form suitable for inspection, review, and copying by the agency". This ensures a complete and continuous record of all changes to the file is maintained. Through the file protection and archiving capabilities of the Multisizer<sup>TM</sup> 3 software, it can be ensured, in compliance with Section 11.10, Paragraph (e), that "Such audit trail documentation shall be retained for a period at least as long as that required for the subject electronic records".

To access the **File History** information, open a file as described in <u>Section 7.2</u>, *A*→**Run File** and select **Get File History** 

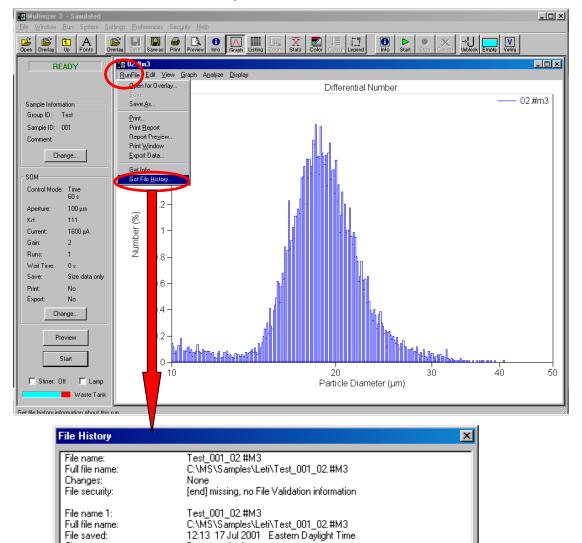

Print.

Changes:

User ID:

User type: Computer user's name:

User log-in name:

Computer's name:

Windows version:

User full name:

Data acquired

Jason Anderson

Janderson

JA Administrator

lobeia 78FRTNL


4.10

Close

#### 10.6.5 Setting Audit Trail


In addition to the auditing associated with the electronic file itself the Multisizer<sup>TM</sup> 3 software offers two additional levels of auditing. First, a general system audits trail that records and stores information at a system level, for example who logged in when, when users were added to the system etc. This is analogous to the security features of a Windows NT<sup>©</sup> operating system; this can be useful for systems using other operating systems. Second, the Multisizer<sup>TM</sup> 3 software generates an error log that stores and records errors associated with the system e.g. missing files, bad communications etc.

In the Administrator drop down menu select Audit Trail Settings.

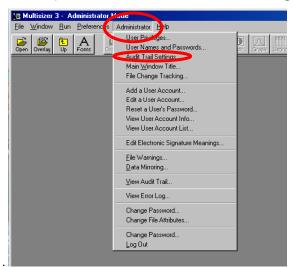

Select the options in the dialog box and  $\mathcal{P} \rightarrow OK$ .

Note that the dialog has a button marked 21 CFR 11. When activated this ensures all options are set so that the settings meet with the rules as applicable to 21 CFR part 11.

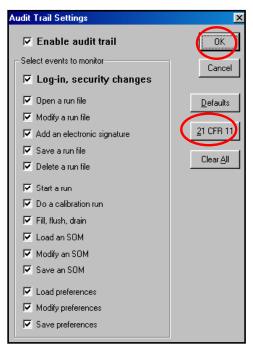

#### **10.6.6 Assigning user privileges**

The level of access to the Multisizer<sup>M</sup> 3 software system is controlled by the rights granted through the **User Privileges** dialog box. The operations granted through these options are linked to the level of security chosen and to the **"User Category"** to which a user is assigned. Users can be classified into five different categories.

- Reviewer
- Operator
- Advanced Operator
- Supervisor
- Administrator

To access the **User Privileges** dialog box, in the **Administrator** drop-down menu select **User Privileges**. At this moment the access level to the software can be assigned to the different categories of users.

If all the user accounts have not been created yet, at this moment they can be created, as well as the **User Names and Passwords** and the **Audit Trial Settings**. These settings can be accessed using the three buttons located in the right lower part of the **User Privileges** dialog box.

User	Privi	lege:	\$						×
Reviewer Operator				Re <sup>.</sup>	viewe Op(	r erator		ΟΚ	
Advanced Operator				T.	Adv	/ance	d Operator		
			Su	pervisor				Sup	Dervisor Cancel
				Administrator I					Administrator
	◄	◄	☑	🔽 Must log in			₽	₽	Can open a file with changes allowed (not read-only)
	◄	▼	Γ	🔲 Name is shown in log-in dialog			◄	▼	🔽 Can edit sample info
	▼	▼	◄	Must enter a password	Γ		Γ	▼	🔽 Can edit size data
	◄	◄	☑	🔽 Can add an Electronic Signature		Γ		◄	🔽 Can change a Kd
	Γ	Γ	☑	🔽 Can edit the aperture list				▼	Can save a file without read-only attribute
	Γ	Γ	☑	🔽 Can calibrate an aperture				◄	🔽 Can overwrite a file
	◄	☑	☑	🔽 Can change the aperture tube				◄	🔽 Can delete a file
		Γ	◄	🔽 Can create an SOP			Γ	Γ	🔽 Can change a file's attributes
	◄	☑	☑	🔽 Can change an SOM		Γ		◄	🔽 Can open a file which has no [end] section
	Γ	☑	☑	🔽 Can save an SOM		Γ	Γ	▼	🔽 Can open a file which has no File Validation section
	⊽	◄	◄	Can change preferences		Γ		⊽	🔽 Can open a file which fails File Validation testing
₹	◄	☑	☑	🔽 Can save preferences	◄	◄	◄	▼	🔽 Can allow Windows to shut down
					▼	◄	◄	◄	🔽 Can exit the Multisizer 3 program
- 5	et defa	aultus	er nri	vileges	- H	ser ac	count	s (ena	abled/total)
	) No						Adm	`	
			2	Set		0/0	) Sup	erviso	r(s) Add a User
C Low security			0/0	) Adv	anced	d Operator(s)			
C Medium security			0/0	) Ope	rator(:	s) <u>U</u> ser Settings			
0	) Hig	h sec	urity (;	21 CFR 11)		0/0	) Rev	riewer	(s) Audit <u>T</u> rail Settings

Select the options in the dialog box  $\mathcal{D}$  **Set** and then  $\mathcal{D}$  **OK**.

Note that the dialog has a button marked High Security 21 CFR 11. When activated this ensures all options are set so that the settings meet with the rules as applicable to 21 CFR part 11.

#### 10.6.7 Electronic Signature

In Subpart A, Section 11.3, an electronic signature is defined as "a computer data compilation of any symbol or series of symbols executed, adopted, or authorized by an individual to be the legally binding equivalent of the individual's handwritten signature". Subpart C, Section 11.100 of the regulation defines the general requirements of such a manifestation. Paragraph (a) states that "each electronic signature shall be unique to one individual and shall not be reused by, or reassigned to, anyone else". These two paragraphs, taken together, mean that an electronic signature is some computer representation of a user's identity, developed to insure the distinct and unique identity of that user. The procedural aspect of Section 11.100 requires that before any such electronic representation is applied, the organization first must "verify" the identity of that individual. The subsequent use of electronic signatures as the "legally binding equivalent of traditional handwritten signatures" then must be "certified" to the agency in writing

Subpart C, Section 11.200, refers to biometric and non-biometric forms of electronic signature. Biometric signatures are defined in Subpart A, Section 11.3 as a "a method of verifying an individual's identity based on measurement of the individual's physical feature(s) or repeatable action(s) where those features and/or actions are both unique to that individual and measurable". Biometrics are generally regarded as techniques such as fingerprints or retinal scans, which are considered to be totally unique to each individual and require specific forms of scanning devices to read and interpret. Non-biometric signatures are those that are computer generated and, as per Section 11.200, "Employ at least two distinct identification components such as an identification code and password". It is this form of electronic signature that is supported by the Multisizer™ 3 software.


#### **10.6.8 Generating electronic signatures**

The Multisizer<sup>™</sup> 3 software employs User ID's and passwords to verify the identification of each user logging into the system. When using this technique, Subpart C, Section 11.300 of the regulation requires "maintaining the uniqueness of each combined identification code and password, such that no two individuals have the same combination of identification code and password". This section also requires that the "identification code and password issuance's are periodically checked, recalled, or revised". Multisizer<sup>™</sup> 3 software supports both of these provisions.

The administration of the system requires that individuals are added to the list of valid Multisizer<sup>™</sup> 3 users via the **Add a User Account** dialog screen. The "identification code" or username of each Multisizer<sup>™</sup> 3 user must be unique. No two users on the same Multisizer<sup>™</sup> 3 system can have the same user name. It is also required that these users supply a password to access the Multisizer<sup>™</sup> 3 software, thus satisfying the requirement to "employ at least two distinct identification components such as an identification code and password". Passwords can be controlled to prohibit the use of duplicates and to force the selection of new passwords after a prescribed period of time.

By the implementation of these features, the Multisizer<sup>™</sup> 3 software can satisfy the requirement that "identification code and password issuance's are periodically checked, recalled, or revised".

Users may be authorized to generate electronic signatures by the administrator. This authorization is granted in the **User Privileges** dialog box. When a user has been authorized, he can electronically sign a file. To generate the electronic signature, open a file as described in <u>Section 7.2</u>, ↔ → Edit, and select Add Electronic Signature.

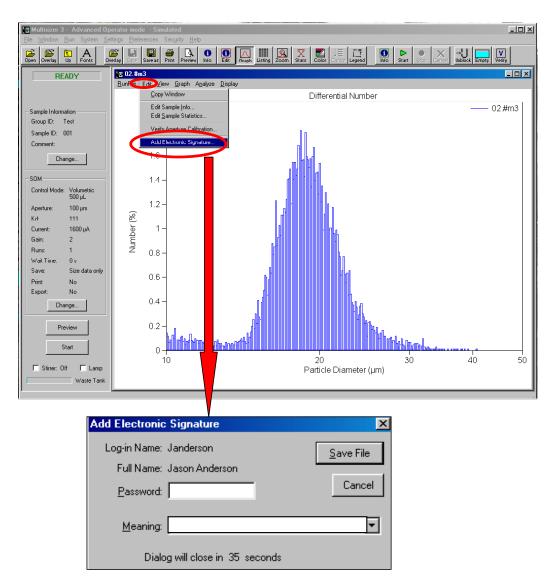

#### **10.6.9** Applying electronic signatures

Subpart C, Section 11.200 stipulates several requirements for the control of electronic signatures. Procedurally, the regulations require that electronic signatures "be used only by their genuine owners" and that they "be administered and executed to ensure that attempted use of an individual's electronic signature by anyone other than its genuine owner requires collaboration of two or more individuals". Through the application of Multisizer<sup>TM</sup> 3 user and password configuration procedures, the system can be configured to "ensure" that inappropriate use of these identifiers can be performed only by the intentional divulgence of security information.

Section 11.200 further specifies the use of electronic signature components during a period "when an individual executes a series of signings during a single, continuous period of controlled system access", and "when an individual executes one or more signings not performed during a single, continuous period of controlled system access". This section of the document represents the "heart" of electronic signature application. To comply with these provisions, the Multisizer<sup>™</sup> 3 software uses the application of the username and password to authenticate the user making and saving the changes, in conjunction with file history and audit trailing, "to independently record the date and time of operator entries and actions that create, modify, or delete electronic records".

#### 10.6.10 Additional security features

The Multisizer  $\mathbb{M}$  3 software offers two important additional levels of security, that whilst not specifically called for by the regulations, make defining and implementing system policies easier.

**Data Mirroring:** The administrator has the option to securely store files in a separate location and in addition to that specified by a user. This is done by enabling the data mirroring option that can only be accessed under administrator mode.

ок
Cancel
ĺ.

File Attributes: Files can be protected to varying degrees using the file attribute options.

Change File Attribu	ıtes			×
D:\Instrument Data\L	_S Data\Ato tech			
File <u>N</u> ame: .\$*		All LS (*.\$*)	]	
Directories: [] [-c-] [-d-] [-e-]	Eiles: NANO G.\$03 NANO G.\$04 Nanog.\$01 Nanog.\$02		Change Change Read-only Hidden Archive System <u>Apply</u>	Files to <u>Show</u> Normal O Only Read-only O Only Hidden O All
Update	Get <u>I</u> nfo			<u>C</u> lose

#### 10.7 <u>Starting the Multisizer 3 software with enabled security.</u>

After all settings for security have been configured, every time the Multisizer 3 software is started, the user is prompted to enter a user name and password. The software will automatically configure according to the privileges granted to the user by the administrator.

After using the Administrator, Supervisor and/or Advance Operator modes, in order to log off you must close down the Multisizer 3 program.

Back to TOC

# **SECTION 11 SPECIFICATIONS**

#### 11.1 **Dimensions and weight**

#### Weight, Dimensions and Power (excluding computer)

Unpacked weight:	45 kg (99 lb)
Width:	45 cm (18 in)
Depth:	60 cm (24 in)
Height:	60 cm (24 in)

#### Input Voltage within Set Ranges:

100 - 120 VAC; 230 - 240 VAC + 10%; single phase

Supply Frequency:	47 to 63 Hz inclusive
Power:	Less than 55 volt-amps (watts)
Fuse types:	240 V, IEC (5x20 mm), Time delay (T), 3.15 A

A The use of other fuse types may result in damage to the instrument.

#### 11.2 **Environmental Conditions**

- This instrument is safe for indoor use only. a)
- Installation category: 11. b)
- Pollution degree: 1. C)

Operating Temperature:	5°C to 40°C
Relative Humidity:	10% to 80% without condensation
Altitude:	Up to 2000 m (5500 ft). Above 7300 m (20,000 ft) unpressurized transport may damage the monitor.
Electrolyte Solutions:	All aqueous and non-aqueous electrolyte solutions devised for use with aperture technology will be suitable for use with the Multisizer 3. Electrolytes should be compatible with glass, fluoropolymers, fluoroelastomers, and polypropylene.

#### Solvents unsuitable for use with silicone rubbers and silicone adhesives are unlikely to be suitable for the Multisizer 3.

#### 11.3 **Operational Parameters**

Particle Size Range:	$0.4\mu m$ to $1200\mu m$ in diameter. $0.0336fL$ to $904.8 \times 10^{6} fL$ or $\mu m^{3}$ in volume. The upper size limit is dictated by solution viscosity and particle density, and the lower limit by the aperture size, electrolyte conductivity and environmental conditions.
Orifice Tube Sizes:	20 $\mu$ m to 2000 $\mu$ m apertures (nominal sizes). Each size allows the measurement of particles in a range of 2% to 60% of aperture diameter.

Aperture	Range	Range
(µm)	(μm)	(fL or μm <sup>3</sup> )
20	0.4 – 12	0.0336 - 905
30	0.6 – 18	0.113 – 3054
50	1.0 – 30	0.524 – 14140
70	1.4 – 42	1.440 – 38800
100	2.0 - 60	4.189 – 113130
140	2.8 – 84	11.49 – 310330
200	4.0 – 120	33.5 – 904830
280	5.6 – 168	92.05 – 2.483x10 <sup>6</sup>
400	8.0 – 240	268 – 7.238 x10 <sup>6</sup>
560	11.2 – 336	735.6 – 19.86 x10 <sup>6</sup>
1000	20 – 600	4189 – 113.1 x10 <sup>6</sup>
2000	40 - 1200	33540 – 904.8 x10 <sup>6</sup>

Number of Channels:	Pulse data is digitized and can be processed to achieve up to 300 size channels for any sleeted range. The number of channels and range can be reprocessed as necessary.	
Digital Pulse Processor:	High-speed digitalization of the signal allows the use of pulse area analysis and other techniques for additional particle characterization. This information may be used at a later time to recalculate the size distribution using different parameters.	
Pulse Data:	Up to 525,000 pulses per analysis.	
Linearity:	± 1 % for diameter	
	± 3 % for volume.	
Aperture Current Range:	30 μA – 6000 μA in 0.2 μA steps	
Aperture Current Accuracy:	± 0.4% of setting	
Polarity error:	Less than 0.5%	
Time mode range:	0.01 to 1000 seconds, selectable in 10 ms increments	
Metering System:	Mercury-free wide range metering pump	
Metered Sample Volumes: (Volumetric mode only)	Continuously selectable from 50µl to 2000µl	
Volumetric pump accuracy:	Better than 99.5%	

## 11.4 Regulatory Compliance

The software is 21 CFR part 11 compliant

Back to TOC

#### 2

21 CFR Part 11, 118

#### A

ACCUVETTE<sup>®</sup>, 21, 31, 37 Administrator Mode, 120, 121, 124, 125 Analysis Report, 48 Analysis Settings, 40 Analytical volume, 52 Aperture, 1, 2, 3, 4, 5, 6, 7, 18, 20, 21, 22, 23, 24, 26, 29, 30, 31, 35, 36, 37, 40, 44, 46, 50, 58, 76, 98, 99, 100, 129 Aperture current Excessive, 7 Aperture List Edit, 35 Aperture Tube, 5, 21, 22, 23, 35 Aperture Tube Changing. See Removing Aperture Tube Removing, 23 Aperture Tube, 18, 19, 21, 22, 26, 35, 46 Fitting, 21 Installing, 24 Setting Up. 22 Aperture Viewer, 18, 30 Apertures Damaged, 7 As Accumulated, 104 Audit Trail, 122, 124 Setting, 124 Average, 68, 69, 70

#### В

Background, 55, 56, 57, 81, 82 Running, 55 Subtracting, 56 Between Cursors, 98, 99, 100 Bin Spacing, 50

#### С

Calibration, 6, 29, 33, 34, 38 Verification, 29 Calibration constant, 6 Calibrator, 31, 33, 39 Choosing, 31 Running, 33 Cautions, 5, 6 Chemical, 6 Electrical, 6 Cells, 28, 44, 104 Change Aperture Wizard, 23, 35 Channels, 2, 4, 40, 50, 86, 98, 99, 130 Cleaning, 58 Compare Statistics, 84 Concentration Index Bar, 32, 38 Connect to Multisizer 3, 14, 17 Connections, 9 Control Mode, 41, 42 Manual, 41 Modal Count, 41 Selecting, 42 Time, 41 Total Count, 41 Volumetric, 41 Control Program, 14 Control Sample, 53, 54, 84 Convert Pulses to Size, 2, 49 Coulter Principle, 1, 2. Current and gain, 28, 43, 44, 46 Current and Gain, 28 Setting, 28 Custom logo, 107

#### D

Data Export, 106 Data Mirroring, 128 Default Preferences, 80, 110 Demodata, 61, 80 Density. 53 Destination Directory, 48 Diameter, 50, 104 Digital Pulse Processing, 2 Dilution Factor, 54 Disconnect from Multisizer 3, 15 Dispersant agent, 45 Distribution, 42, 95 Cumulative <, 97 Differential. 95 Types, 95 Distribution Types, 95

#### Ε

Electrical interference, 7 Electrical Sensing Zone, 1, 2 Electrode, 18, 20 Electrolyte, 5, 6, 7, 9, 11, 21, 25, 26, 31, 35, 37, 40, 44, 45, 46, 52, 55, 58, 59, 129 Electrolyte Selecting, 26 Electrolyte Volume, 52, 81 Electronic Records, 118 Electronic Signature Generating, 126 Applying, 127 Electronic Signatures, 118 Enter Sample Info, 51, 53 Equivalent Spherical Diameter, 2, 53 ESZ. See Electrical Sensing Zone Exporting Data, 47

#### F

FDA, 118 File Opening, 60 File Attributes, 128 File History, 122, 123 Getting, 123 File Name, 47, 53 Generating, 47 Flushing, 46 Fonts, 107 Frontal Panel, 18

#### G

Graph Options, 106

Т

Installation, 8 Instrument Controls, 15 International Standard ISO 13319, 2 Interpolation Points, 86

#### K

Kd, 6, 34, 35, 81

## L

Μ

Language, 109 Linear X axis, 97 Location, 8 Logarithmic X axis, 97 Login, 122

Main Menu Bar, 14, 17, 23, 33, 35, 38 Mass, 3, 52, 53, 95 Metering pump, 58 Metering Pump, 27 Adjusting, 27 Modal value, 31, 33, 39 Multiple runs, 42 Multisizer II Edit, 50 Multi-Tube Overlap, 76, 77

#### Ν

Noise Level, 29, 43 Measuring, 29

#### 0

Operational checks, 122 Optics, 30 Aligning, 30 Overlay, 63, 64, 65, 66, 67 Open, 65, 67 Save, 66 Overlaying Files, 63

#### P

Page Setup, 107 Particle Concentration, 42, 58 Using Large Apertures, 58 Particle Size, 1, 2, 76 Preferences, 80, 94, 101, 102, 103, 105, 106, 107, 108, 109, 110, 111, 114, 115 Costumizing, 94 Saving, 109 Preferences File, 80, 114 Preferences Menu. 80 Preview, 32, 37, 57 Prferences, 80. Menu Printed. See Printed Report, 85, 86 Printer, 93 Printing, 48, 93 Pulse, 1, 2, 46, 89, 90, 94, 102, 103, 104, 118.130 Pulse Graph, 103 Pulse Listing, 103 Pulses, 2, 49, 89, 102 Pulses Graph, 89

#### R

Regulatory Compliance, 118 Report, 3, 9, 48, 80, 82, 83, 85, 86, 87, 88, 89, 90, 91, 92, 93 Report, 80, 111 Printing, 80 Results, 60 Saving, 60 Viewing, 60 Reviewer, 125

## S

Sample, 7, 18, 25, 51, 52, 53, 54, 57, 81, 82, 84, 85, 98. Info Running, 57 Sample Analysis, 40 Sample Compartment, 18, 25 Sample Dispersion, 7 Sample Info Short. 82 Standard, 81 Sample Information, 51 Entering, 51 Sample Platform, 18 Sample Statistics, 54, 84, 85 Sample volume, 52 Saving, 46, 50, 80, 109, 111, 115 Security, 14, 15, 119, 120, 125 Setting Up, 119 Security Level, 14, 120 Settings, 41, 49, 50, 51, 56, 57, 109, 111, 116, 117 Shape Factor, 53 Shutting down the Multisizer 3, 17 Size Bins, 50 Size classes, 2, 4, 86, 100 Size distribution, 2, 3, 4, 42, 50, 53, 76, 78, 82, 85, 88 Size Graph, 82, 94, 95, 98 Size Interpolation, 86, 94, 100 Size Listing, 85, 86, 94, 98 Size Statistics, 83, 88, 94, 99 Size Trend, 72, 73, 74, 105 Creating, 72 SOM. See Standard Operating Method SOP. See Standard Operating Procedure Specific Surface Area, 104 Specifications, 129 Standard Operating Method, 41, 47, 49, 50, 51.111 Changing, 41 Saving, 50 Standard Operating Procedure, 51, 109, 112, 114 Saving, 114 Standard Operating Procedures Creating, 111 Standard Operation Procedure, 50, 80 Statistics, 83, 84, 85, 99

Status Panel, 16, 32, 37, 41, 55, 57 Stirrer, 18, 19, 20 Stirrer Rotation Control, 18 Stirrer Speed Control, 18 System, 27, 51, 58, 59 Cleaning, 58

#### Т

Threshold, 42, 43, 50 Toolbar, 108 Trap, 19 Trend, 72, 73, 74, 75, 91, 92, 94, 101, 102, Trend Graph, 91, 94, 101 Trend Listing, 92, 94, 102

#### U

User Account, 120, 121, 126 Adding, 120 User Names and Passwords, 121, 122 Selcting, 121 User Privileges, 125 Assigning, 125 Granting, 125

#### V

Verfication. See Calibration View Options, 103 Voltage, 1, 4, 6, 11, 12, 13 Volume, 58, 88, 104

#### W

Warnings, 5 Chemical, 5 Electrical, 5 Fire, 5 Mechanical, 5 Windows Colors, 108

#### Х

X Axis, 97 Scale, 97 X-Axis, 104